



**Agilent
Genomic Workbench
Lite Edition 6.5
SureSelect Quality Analyzer**

User Guide



Agilent Technologies

Notices

© Agilent Technologies, Inc. 2010

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

Revision

G3800-90041

Revision A0, December 2010

Agilent Technologies, Inc.
5301 Stevens Creek Blvd.
Santa Clara, CA 95051 USA

Trademarks

Adobe® and Adobe® Reader® are registered trademarks or trademarks of Adobe Systems Incorporated in the United States and/or other countries.

Microsoft® is a registered trademark of Microsoft Corporation in the United States and other countries.

Software Revision

This guide is valid for 6.5 and later revisions of the Agilent Genomic Workbench Lite Edition software, until superseded.

Warranty

The material contained in this document is provided “as is,” and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.

Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

Restricted Rights Legend

U.S. Government Restricted Rights. Software and technical data rights granted to the federal government include only those rights customarily provided to end user customers. Agilent provides this customary commercial license in Software and technical data pursuant to FAR 12.211 (Technical Data) and 12.212 (Computer Software) and, for the Department of Defense, DFARS 252.227-7015 (Technical Data - Commercial Items) and DFARS 227.7202-3 (Rights in Commercial Computer Software or Computer Software Documentation).

Safety Notices

CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

In This Guide...

This guide describes how to use the SureSelect Quality Analyzer program, which is part of the Agilent Genomic Workbench Lite Edition program suite.

1 Getting Started

This chapter describes SureSelect Quality Analyzer and how it fits into the typical SureSelect Target Enrichment research flow. It gives instructions on how to start both the Quality Analyzer and Workflow programs, and includes two example exercises, one that leads you through a typical SureSelect QC analysis experiment and one that shows you how to set up and run a workflow to sequentially and automatically produce QC reports and enrichment analyses.

2 Using SureSelect Quality Analyzer Interactively

This chapter describes how to import and organize sequence read and target interval files into analysis experiments. It describes how to display and analyze the reads and target intervals, and how to create and display QC reports of several types, including summary QC metrics, and reports on read depth and enrichment analyses.

3 Setting Up and Running Workflows

This chapter gives instructions on how to set up and run workflows for producing QC reports sequentially and automatically using read files, chromosome mapping files and target interval files.

4 SureSelect Quality Analyzer Reference

This chapter describes the tabs, panes, commands, menus, shortcut menus, parameter panels and dialog boxes that can appear when you use the SureSelect Target Enrichment application, which includes the Quality Analyzer and Workflow programs. It also contains reference information on the reports that the program creates.


Contents

1	Getting Started	9
	SureSelect Quality Analyzer	10
	Information for Mac Users	12
	Setting Up SureSelect Quality Analyzer	13
	Before you use SureSelect Quality Analyzer	13
	To start the SureSelect Target Enrichment application	14
	To start SureSelect Quality Analyzer	17
	To set the application type	18
	To select where data and results are stored	19
	Example Exercise 1 – Interactive	20
	Before you start the example exercise	20
	Step 1. Import data and create an experiment	21
	Step 2. Analyze data and display results	23
	Setting Up SureSelect Quality Analyzer Workflows	27
	To start the Workflow program	27
	Example Exercise 2 – Workflow	30
	Before you start the example exercise	30
	Step 1. Set up the workflow	31
	Step 2. Run the workflow and review results	34
	Getting Help	35
	To get help within Agilent Genomic Workbench	35
	To contact Agilent Technical Support	36
	To learn about Agilent products and services	36
2	Using SureSelect Quality Analyzer Interactively	37
	Importing Files	38
	To import sequence read files	38
	To import target genomic regions as a track	41

To import a genome build	43
Setting Up QC Experiments	44
To create a new experiment	44
To select an experiment for analysis	46
To deselect an experiment	47
To view the data in an experiment	47
To view the intervals in a target interval track	50
To export tracks	50
To remove a track from the program	51
To customize the appearance of an experiment in the Genomic Viewer	51
To delete an experiment	54
To save an experiment	54
To copy what appears on your screen to the clipboard	54
Creating Target Enrichment QC Reports	55
To analyze read quality	55
To view and save reports	56
3 Setting Up and Running Workflows	59
Creating and Managing Workflows	60
To create a new workflow	60
To edit an existing workflow	60
To save a workflow	61
To save a workflow to a new name	61
To delete a workflow	61
To set a password for an existing workflow	61
Setting Up/Editing Workflows	62
To select target interval input	63
To select sequencing input	65
To select reports	67
To rename the experiment	68
Running Workflows	69

To run a workflow	69
To run a series of workflows	70
To monitor workflow runs	71
To abort a single workflow	72
To abort all workflows	72
To display the results of a workflow	73
4 SureSelect Quality Analyzer Reference	75
SureSelect Target Enrichment Main Window	76
Switch Application menu	78
SureSelect Target Enrichment tab bar	79
Command ribbon – Home tab	80
Command ribbon – Help tab	82
Quality Analyzer Tab Window	83
Command ribbon – Quality Analyzer tab	85
Navigator overview	87
Navigator – Experiment pane	88
Genomic Viewer – overview	91
Genomic Viewer – Genome View	93
Genomic Viewer – Chromosome View	95
Genomic Viewer – Gene View	96
Genomic Viewer – Tab View	98
Workflow Tab Window	100
Command ribbon – Workflow tab	102
Workflow Navigator	103
Parameter Panels – Overview	105
Experiment Parameter Panel	106
Import Track Parameter Panel	107
Select Imported Track Parameter Panel	108
Sequencing Input Parameter Panel	109
Text Summary Report Parameter Panel	112
Workflow Run button	114

Summary Console/Progress tabs – Overview	115
Summary Console tab	116
Named workflow tabs	118
Dialog Boxes	120
Choose Track Color	120
Create Experiment	124
Create Workflow	126
Export	127
Export Tracks	128
Go to Gene/Genomic Location	129
Import Genome Build	130
Import Track	131
Open	132
Provide Workflow Identifier	133
Set Password	134
Select SEQ Files	135
Specify Input Information	136
Track	139
User Preferences – Tracks tab	144
View Preferences	147
Target Enrichment Reports	149
QC metrics	150
Read depth distribution	152
Enrichment analysis results – UCSC	153
Enrichment analysis results – XLS report	154

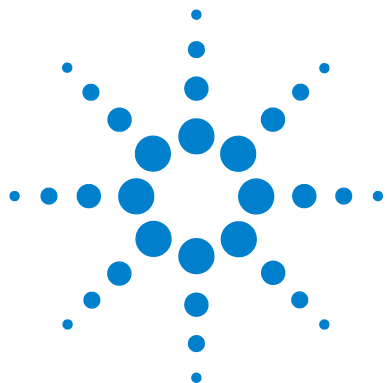

1

Getting Started

SureSelect Quality Analyzer	10
Information for Mac Users	12
Setting Up SureSelect Quality Analyzer	13
Example Exercise 1 – Interactive	20
Setting Up SureSelect Quality Analyzer Workflows	27
Example Exercise 2 – Workflow	30
Getting Help	35

This chapter describes SureSelect Quality Analyzer and how it fits into the typical SureSelect Target Enrichment research flow. It gives instructions on how to start the program, and describes the main program window. It describes the files that you must have before you can use the program, and includes an example exercise that leads you through a typical SureSelect QC analysis experiment.

In addition, this chapter shows how you can set up a workflow to sequentially and automatically produce QC reports and enrichment analyses from the required files. A second exercise helps you understand the Workflow process.

You can also find out how to get additional help on other Agilent Genomic Workbench Lite Edition programs.

For information on how to run a Feature Extraction workflow, see the *Workflow User Guide*.

To find more information on all the Agilent Genomic Workbench Lite Edition programs, see the *Product Overview Guide*. Click the Open Application tab if it is not already open. Access to the guide is in the top right corner of the Open Application window.



SureSelect Quality Analyzer

SureSelect Quality Analyzer is a program that lets you assess the effectiveness of the pull-down of targeted genomic fragments when you use the Agilent SureSelect Target Enrichment system. This system uses libraries of biotinylated RNA oligonucleotide “baits” to harvest genomic DNA fragments of interest for sequencing. This forms the basis of a powerful selection method that lets you focus your sequencing efforts.

The selective nature of the process makes it ideal for targeted resequencing that uses next-generation sequencing technology. After you sequence the harvested fragments, you can use SureSelect Quality Analyzer to calculate statistical metrics and analyze read depth. You can also do an enrichment analysis, and display the results in the UCSC Genome Browser next to the annotation tracks of your choice.

SureSelect Quality Analyzer is part of the Agilent Genomic Workbench suite, a collection of essential programs that supplies a single solution for all of your genomics data processing needs. [Figure 1](#) shows a typical sequencing process that uses the Agilent SureSelect Target Enrichment system.

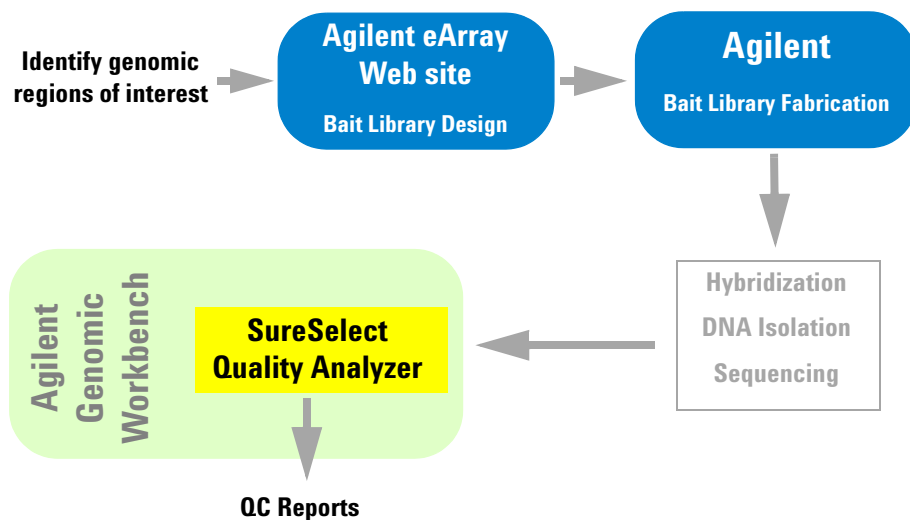


Figure 1 Typical sequencing process with the SureSelect Target Enrichment system

After you identify the genomic regions of interest, you use the Agilent eArray Web site to design a library of oligonucleotide baits to the desired regions. You submit the library to Agilent, who supplies the biotinylated RNA oligonucleotides in the library to you in any desired quantity. You use the bait library to enrich your sample DNA for the desired target genomic regions, and sequence the resulting DNA fragments.

You import the sequence reads into Agilent Genomic Workbench, and use SureSelect Quality Analyzer to do several crucial quality-related analyses:




- **QC metrics** – Includes overall statistics, such as the percentage of reads in targeted regions, the overall fold enrichment, and the average read depth.
- **Read depth analysis** – Shows the overall distribution of reads over the targeted regions and targeted bases. Distributions are available both as tables and as graphs.
- **Enrichment analysis** – Shows the degree of enrichment for target intervals in the genome of interest. The program creates a WIG format file, and uploads it to the UCSC Genome Browser, where it appears graphically in the User Track area. For information on the WIG file format, go to genome.ucsc.edu/FAQ/FAQformat.

You can also create a text version of the enrichment analysis results as an *.xls file.

Like the CGH and ChIP applications, the SureSelect Target Enrichment application lets you set up and run workflows to sequentially and automatically import the required files and use them to produce reports from the quality-related analyses described above.

Information for Mac Users

The content of this User Guide applies to both the Windows and Mac versions of Agilent Genomic Workbench Lite Edition. Both of these versions have the same features. However, when you use the Mac version of the program, please note the following:

Windows command	Equivalent Mac command
Right-click	<ul style="list-style-type: none">• Command-click (-click)• On Macs with trackpads, other options are available. On certain machines, you place two fingers on the trackpad while you press the button below the trackpad. See the user guide for your specific machine.• If you have a third-party mouse that has more than one button, you may be able to use one of the buttons as a right mouse button.
Control-click	Control-click (Same as the Windows command)
Shift-click	Shift-click (Same as the Windows command)
 (Close button)	 (Close button)

Setting Up SureSelect Quality Analyzer

This section describes what you must do before you can use SureSelect Quality Analyzer, and explains how to start the program. It also includes a description of the main program window as it appears for the SureSelect Target Enrichment application type.

Before you use SureSelect Quality Analyzer

- You must have several types of files available before you can use SureSelect Quality Analyzer:
 - Sequence read file** – A gzipped or *.zip archive that contains the nucleotide sequences and locations of the fragments in your DNA sample, along with related header information. The program supports these read file types:

File Type	Comments
Illumina GA Export	Revised version of the ELAND file format that contains additional fields, including quality scores for all bases. You cannot import a multiplexed read file into Agilent Genomic Workbench. You must first separate the file into individual indexed sample files, then import these.
AB SOLiD ma	Color space *.ma file produced by the Applied Biosystems SOLiD system. The AB SOLiD alignment software produces this CSFASTA format file, and places sequence alignment information in the FASTA headers. Color calls are converted to bases when the file is imported.

NOTE

You can import sequence read files in the gzipped or *.zip formats, but not in the .zipx format. In addition, Agilent Genomic Workbench cannot handle a compressed file whose uncompressed file is greater than 4G. In this case, import the uncompressed file directly.

1 Getting Started

To start the SureSelect Target Enrichment application

- **Chromosome mapping file** – A *.txt file that contains the names of the chromosomes that are referenced in the read file, correlated with the names of the chromosomes in the desired genome build of your species of interest. For details, see [“To import sequence read files”](#) on page 38.
- **Target intervals file** – A *.bed format track file that contains a list of the target genomic intervals. Typically, these are the target genomic intervals to which you designed your SureSelect Target Enrichment baits. If you used eArray to create the bait library, you can download this file from eArray when you download the library. You can also create the target intervals file manually.

To start the SureSelect Target Enrichment application

In Agilent Genomic Workbench, content as well as function are partitioned by experimental application type. When you select an application type, only the tools and content that are relevant to the selected type are accessible. The content and functionality for SureSelect Quality Analyzer are only available when you select **SureSelect Target Enrichment** as the application type.

1 Start Agilent Genomic Workbench.

If upon starting the program you see unfamiliar dialog boxes and don't know how to proceed, please refer to “To start the Agilent Genomic Workbench user/client software” in the *Installation Guide*.

2 Click **Switch Application**.

3 Select **SureSelect Target Enrichment**.

The tools and content relevant to this application type become available.

For details on the other options that are available, see [“Switch Application menu”](#) on page 78.

When you first switch the application to the SureSelect Target Enrichment application, you see the window on the next page for the Home tab.

To start the SureSelect Target Enrichment application

The Tab bar and Switch Application button remain the same for every tab. But the Command ribbon changes for each tab, and the Tab panes may or may not stay the same.

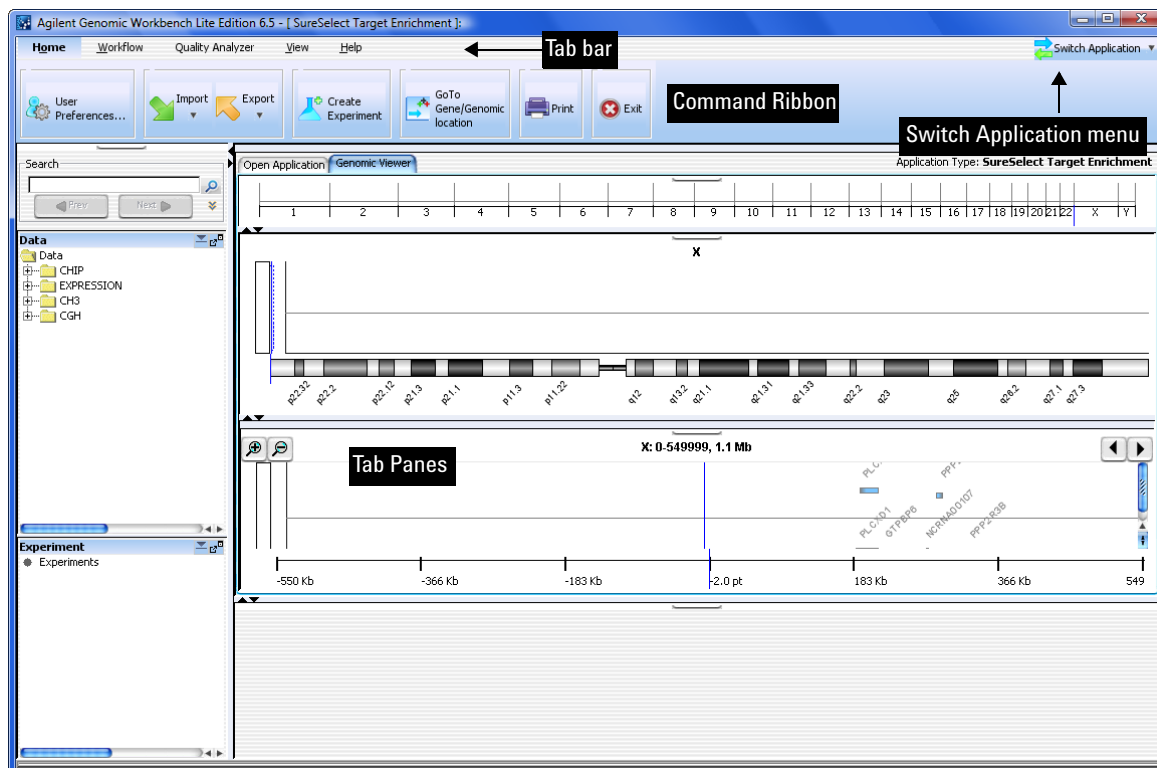


Figure 2 SureSelect Target Enrichment Main Window – Home tab

1 Getting Started

To start the SureSelect Target Enrichment application

The SureSelect Target Enrichment Main Window has these elements:

Element	Description
Title bar	Displays the program name and the currently selected application type.
Switch Application menu	Lets you select the experimental application type. This is one of the main settings for Agilent Genomic Workbench, since functionality as well as design, sample, and sequence data are partitioned by application type. See “Switch Application menu” on page 78.
Tab bar	Contains the tabs that let you use the main functions of the SureSelect Target Enrichment program. See “SureSelect Target Enrichment tab bar” on page 79.
Command ribbon	<p>Contains the commands that you use to access the main functionality of the SureSelect Target Enrichment application. A different command ribbon appears for each tab. The Command ribbons for the Quality Analyzer tab window and the Workflow tab window are described in separate sections.</p> <p>See these topics:</p> <ul style="list-style-type: none">• “Command ribbon – Home tab” on page 80• “Command ribbon – Help tab” on page 82• “Quality Analyzer Tab Window” on page 83• “Workflow Tab Window” on page 100
Tab panes	<p>These change depending on the tab you choose.</p> <p>When you select either the Home, Quality Analyzer, View, Tool or Help tabs, the Navigator and Genomic Viewer panes appear. See “The SureSelect Target Enrichment Main Window has these elements:” on page 16.</p> <p>When you select the Workflow tab, the Workflow Navigator, Parameter Panels and Summary Console/Progress tabs panes appear. See “To start the Workflow program” on page 27.</p>

To start SureSelect Quality Analyzer

SureSelect Quality Analyzer is part of the SureSelect Target Enrichment application. It includes the functions of the Quality Analyzer tab and the Workflow tab.

- 1 Start the SureSelect Target Enrichment application.

See [“To start the SureSelect Target Enrichment application”](#) on page 14.

- 2 Click the **Quality Analyzer** tab.

The Quality Analyzer tab window appears. See [Figure 3](#).

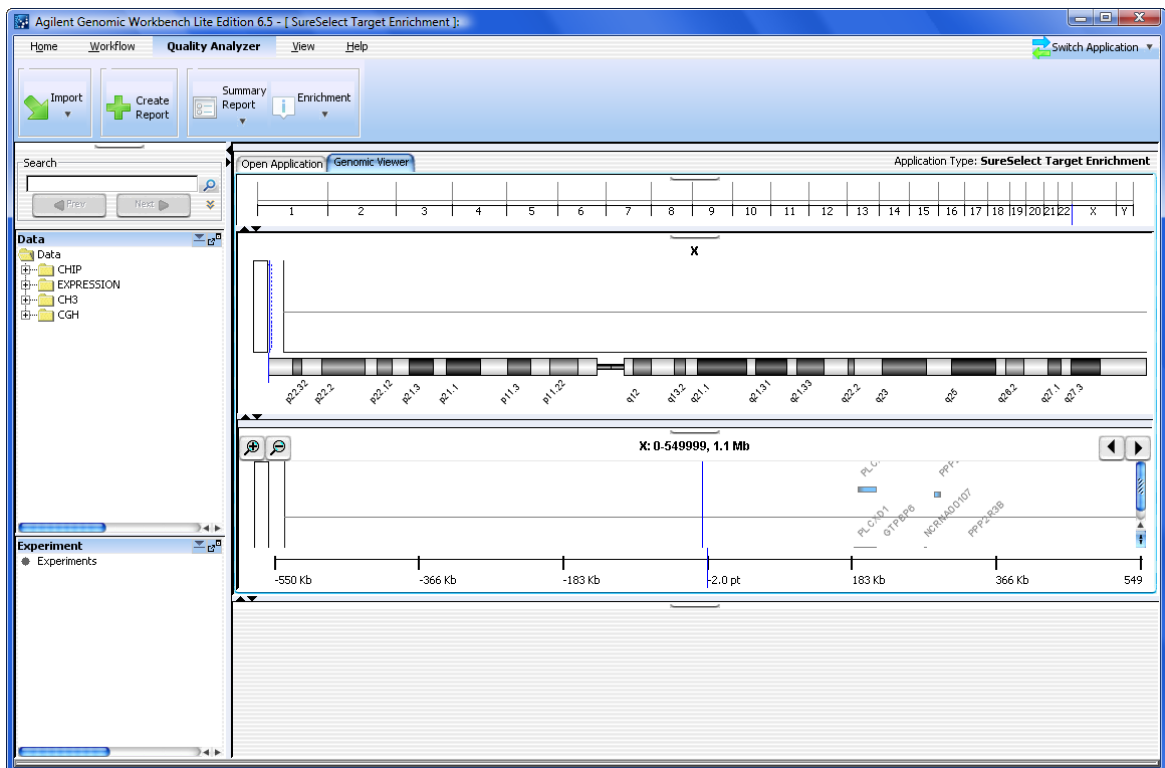


Figure 3 Quality Analyzer tab window

1 Getting Started

To set the application type


The tab window contains the elements described in [Table 1](#).

Table 1 Quality Analyzer tab window elements

Window Element	Description
Command ribbon – Quality Analyzer	Contains the commands for importing sequence read files and target interval tracks, doing QC analyses, and creating QC reports. See “Command ribbon – Quality Analyzer tab” on page 85.
Tab pane – Navigator	Contains lists of the designs, sample data, sequence data, analysis experiments, and other related items that are available to you. A search feature is also available. See these topics: <ul style="list-style-type: none">• “Navigator overview” on page 87• “Navigator – Experiment pane” on page 88
Tab pane – Genomic Viewer	<p>Shows the genome of the selected organism at several levels of detail. Genes, tracks, probes, baits, and data of many types can appear in the Genomic Viewer in the context of the genome.</p> <p>For general information on how to use the Genomic Viewer, see the <i>Data Viewing User Guide</i>.</p> <p>For information on the specific features of the Genomic Viewer that are relevant to the Quality Analyzer tab, see “Genomic Viewer – overview” on page 91.</p>

To set the application type

In Agilent Genomic Workbench, content as well as function are partitioned by experimental application type. When you select an application type, only the tools and content that are relevant to the selected type are accessible. The content and functionality for SureSelect Quality Analyzer are only available when you select **SureSelect Target Enrichment** as the application type.

- 1 Start Agilent Genomic Workbench.
- 2 Click  **Switch Application**.
- 3 Select the desired application type.

The selected experimental application type appears in Application Type. The tools and content relevant to this application type become available.

You can use the tools in Agilent Genomic Workbench Lite Edition to work with experimental data and other related files for many different application types. The program supports these application types:

Application Type	Description
CGH	Analysis of data for comparative genomic hybridization studies
ChIP-on-chip	Analysis of data for chromatin immunoprecipitation studies
CH3	Analysis of data for methylation studies
SureSelect Target Enrichment	QC analysis of target fragment pull-down

To select where data and results are stored

By default, sequence data and analysis results are stored on your computer in the **data** folder, within the folder that contains the Agilent Genomic Workbench program. You can change this location.

- 1 In the **Home** tab, click **User Preferences**.

The User Preferences dialog box appears.

- 2 Click the **Miscellaneous** tab.

The Miscellaneous tab appears. See [“User Preferences – Miscellaneous tab”](#) on page 88.

- 3 In Data Location, click **Browse**.

An Open dialog box appears.

- 4 Select the desired location, then click **Open**.

The selected location appears in Data Location.

- 5 Click **OK**.

Example Exercise 1 – Interactive

In this exercise, you create and view a PDF format QC report that contains summary QC statistics and the results of read depth analysis. You use the UCSC Genome Browser to view the results of enrichment analysis and the Genomic Viewer to view reads and target intervals.

Before you start the example exercise

- Download the example files for this exercise. Follow these steps:
 - a Go to earray.chem.agilent.com.
The login page of the eArray Web site appears. You do not need to log in.
 - b Under Additional Information, click **Download Agilent Genomic Workbench**.
A page of available Agilent Genomic Workbench files appears.
 - c Under Sample Data, next to Sample files for SureSelect Target Enrichment Quality Analyzer, right-click **Download**.
 - d Click **Save Target As** and save the file to your desktop.
The eArray Web site transfers a zip format file to your computer.
 - e Extract the zip file to your desktop.
The folder of example files appears on your desktop. Do not extract the zip format files within this folder.
- Familiarize yourself with the Agilent Genomic Workbench Quality Analyzer interface. See “[The SureSelect Target Enrichment Main Window has these elements:](#)” on page 16 and “[Quality Analyzer Tab Window](#)” on page 83.
- Note that the main steps in the example exercises appear in the first column of the table. More detailed instructions appear in the second column. Additional notes and comments appear in the third column.

Step 1. Import data and create an experiment

In this step you import the read and target interval files (track files) and then create an experiment to contain them.

Example Exercise 1 – Import data and create an experiment

Step	Detailed Instructions	Notes/Comments
1 Start SureSelect Quality Analyzer.	<p>a Start Agilent Genomic Workbench Lite Edition 6.5. The main window appears.</p> <p>b Click Switch Application > SureSelect Target Enrichment. SureSelect Target Enrichment appears in Application Type. The appropriate tools and panes appear in the main window of the program.</p> <p>c Click Quality Analyzer. The Quality Analyzer main window appears.</p>	<ul style="list-style-type: none"> If upon starting the program you see unfamiliar dialog boxes and don't know how to proceed, please refer to "To start the Agilent Genomic Workbench software" in the <i>Installation Guide</i>. You can use the SureSelect Target Enrichment application free of charge. You do not need to purchase or enter a license.
2 Import the read file and chromosome mapping file found in the sample files folder you downloaded. <ul style="list-style-type: none"> Keep File Type as Illumina GA Export. Under Species Info, keep Select Species set to H. sapiens, and Select Genome Build set to hg18. 	<p>a Click Import > Reads. A dialog box appears.</p> <p>b In the folder of sample files that you extracted to your desktop, select the file s_1_1_export_top.zip.</p> <p>c Click Open. The Specify Input Information dialog box appears.</p> <p>d Under Read Info, in File Type, select Illumina GA Export.</p> <p>e Under Map Chromosomes, in Chr Mapping File, click Browse. An Open dialog box appears.</p> <p>f In the folder of sample files that you extracted to your desktop, select the file chrMapping.txt, then click Open. The name of the selected file appears in Chr Mapping File.</p> <p>g Click Import. A progress bar appears.</p>	<ul style="list-style-type: none"> In this step, you import these files: <ul style="list-style-type: none"> Read file – Contains header information, and the sequence data to be analyzed. The program supports read files in several formats. See "To import sequence read files" on page 38. Chromosome mapping file – A *.txt file that associates the names of chromosomes in the read file with the names of chromosomes as they are specified in the selected genome build.

1 Getting Started

Step 1. Import data and create an experiment

Example Exercise 1 – Import data and create an experiment

Step	Detailed Instructions	Notes/Comments
3 Import the target intervals file from your downloaded sample files folder. <ul style="list-style-type: none">Track Name - Example TargetsKeep the Build the same # as for the Reads.	<ul style="list-style-type: none">a Click Import > Track. The Import Track dialog box appears.b Change the Build Name to hg18.c In Track Name, type Example Targets.d In Track File, click Browse. An Open dialog box appears.e In the folder of sample files that you extracted to your desktop, select SureSelect_All_Exon_G3362_with_names.bed, then click Open. The location of the file appears in Track File.f Click OK. A dialog box tells you that the track was successfully imported.g Click OK.	<ul style="list-style-type: none">In this step, you import a BED format track file that contains the desired target genomic intervals.When you select a track file to import, the program only displays *.bed format files.
4 Create a QC analysis experiment with this information: <ul style="list-style-type: none">Name - xxxExampleExpt, where “xxx” are your initials.Read Length - 76Targets - Example TargetsSEQ Sample - s_1_1_export_topSingle End	<ul style="list-style-type: none">a In the Experiment pane of the Navigator, right-click the Experiments node, then click New Experiment. The Create Experiment dialog box appears.b Enter the following information:<ul style="list-style-type: none">Name – Type xxxExampleExpt, where “xxx” are your initials.Read Length – Delete the existing value, then type 76Targets – Select Example TargetsSEQ Sample – Select s_1_1_export_top.c Click OK. A progress bar appears. In the Experiment pane of the Navigator, a new experiment with the name xxxExampleExpt appears.	<ul style="list-style-type: none">In SureSelect Quality Analyzer, an experiment is a required organizational level that relates a specific read file to a specific set of target intervals.

Step 2. Analyze data and display results

Once you have imported the read, chromosome mapping and target interval files and created a new experiment containing them, you analyze the data by selecting the experiment and creating reports.


Example Exercise 1 – Analyze data and display results


Step	Detailed Instructions	Notes/Comments
1 Select the experiment for analysis.	<p>a In the Experiment pane of the Navigator, in the Experiments folder, right-click Example Expt, then click Select Experiment. A dialog box asks if you want to select the experiment.</p> <p>b Click Yes. A progress bar appears. A list of reads appears in Tab View of the Genomic Viewer.</p>	<ul style="list-style-type: none"> Later in this example exercise, you view the reads graphically in the context of the human genome. In the Experiment pane of the Navigator, the name of the selected experiment appears in blue. The reads that appear in Tab View are the ones that are found on the chromosome that appears in Chromosome View.
2 Calculate QC metrics and analyze read-depth and enrichment.	<p>a Click Create Report. A progress bar appears. A dialog box tells you that reports were successfully created.</p> <p>b Click OK.</p>	<ul style="list-style-type: none"> In this step, the program analyzes the sequence data against the intervals in your selected target intervals file. You view the results in subsequent steps.
3 Create and view a PDF format Summary Report. <ul style="list-style-type: none"> Use any name and location you want. 	<p>a Click Summary Report > Save as PDF. A Save dialog box appears</p> <p>b Select a location, and type a name for the report file, then click Save. A dialog box tells you that the file was successfully created, and asks if you would like to view it.</p> <p>c Click Yes. The QC Report opens in Adobe Reader.</p>	<ul style="list-style-type: none"> In this step, you create a PDF format report that contains calculated QC metrics and the results of read depth analysis. For details, see: <ul style="list-style-type: none"> “To view and save reports” on page 56 “QC metrics” on page 150 “Read depth distribution” on page 152 You can also save the report as a text file. See “To view and save reports” on page 56.

1 Getting Started

Step 2. Analyze data and display results

Example Exercise 1 – Analyze data and display results

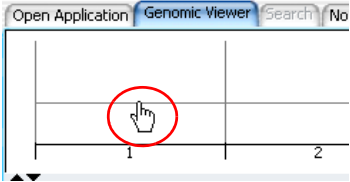Step	Detailed Instructions	Notes/Comments
4	<p>Display enrichment analysis results in the first 20Mb of chromosome 1 in the UCSC Genome Browser.</p> <p>a In the Genomic Viewer, in Genome View, click the center of the Chromosome 1 area.</p>  <p>b Click Enrichment > View in UCSC. Your Web browser opens.</p> <p>c If a security message tells you that the browser has restricted scripts or ActiveX controls, follow the on-screen instructions to allow the blocked content.</p> <p>The UCSC Genome Browser opens in your Web browser.</p> <p>d In position/search, delete the existing chromosomal location, then type <code>chr1:1-20,000,000</code></p> <p>e Click jump.</p> <p>Enriched peaks appear in the User Track area.</p>	<ul style="list-style-type: none">In this step, you transfer enrichment analysis results for chromosome 1 to the UCSC Genome Browser, and view the enriched peaks in the first 20 Mb of that chromosome.You can display the enriched peaks next to your choice of annotation tracks. For instructions, see the online help that is available in the UCSC Genome Browser.You can also save the enrichment analysis results as a *.xls file. See “To view and save reports” on page 56.

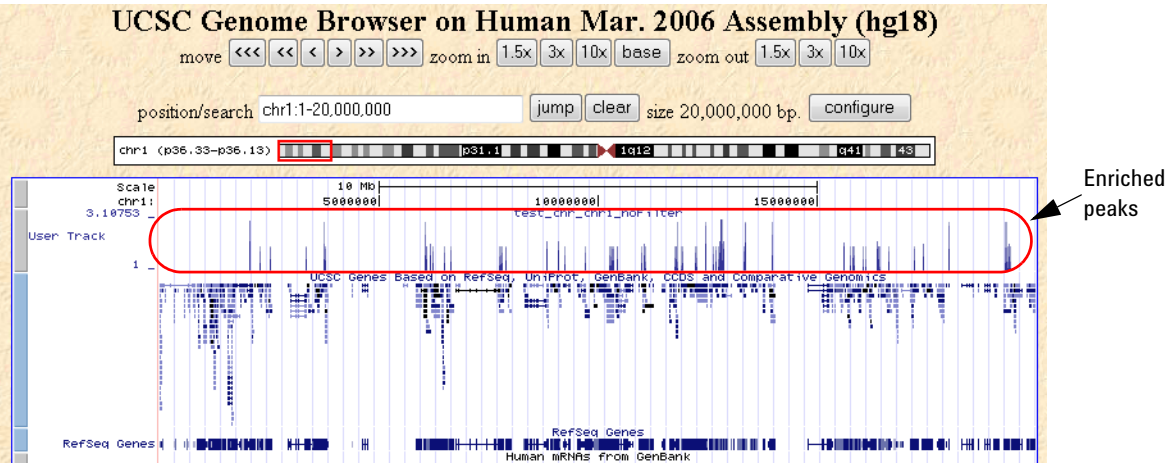

Setting Up SureSelect Quality Analyzer Workflows

This section describes what you must do before you set up SureSelect Quality Analyzer workflows, and explains how to start the program. It also includes a description of the Workflow tab window as it appears for the SureSelect Target Enrichment application type.

If you are endeavoring to set up a workflow without first having created an experiment and analyzed the files interactively with Quality Analyzer, please first read [“Before you use SureSelect Quality Analyzer”](#) on page 13.

To start the Workflow program

The Workflow tab window for the SureSelect Quality Analyzer has a completely distinct user interface from the one for the other tabs for the SureSelect Target Enrichment application.

- 1 Start the SureSelect Target Enrichment application.

See [“To start the SureSelect Target Enrichment application”](#) on page 14.

- 2 Click the **Workflow** tab.

The Workflow tab window appears. See [Figure 4](#).

The Workflow tab window presents a completely different set of panes than the Home, Quality Analyzer, View, Tool and Help tab windows do – Workflow Navigator, Parameter Panels and Summary Console.

1 Getting Started
To start the Workflow program

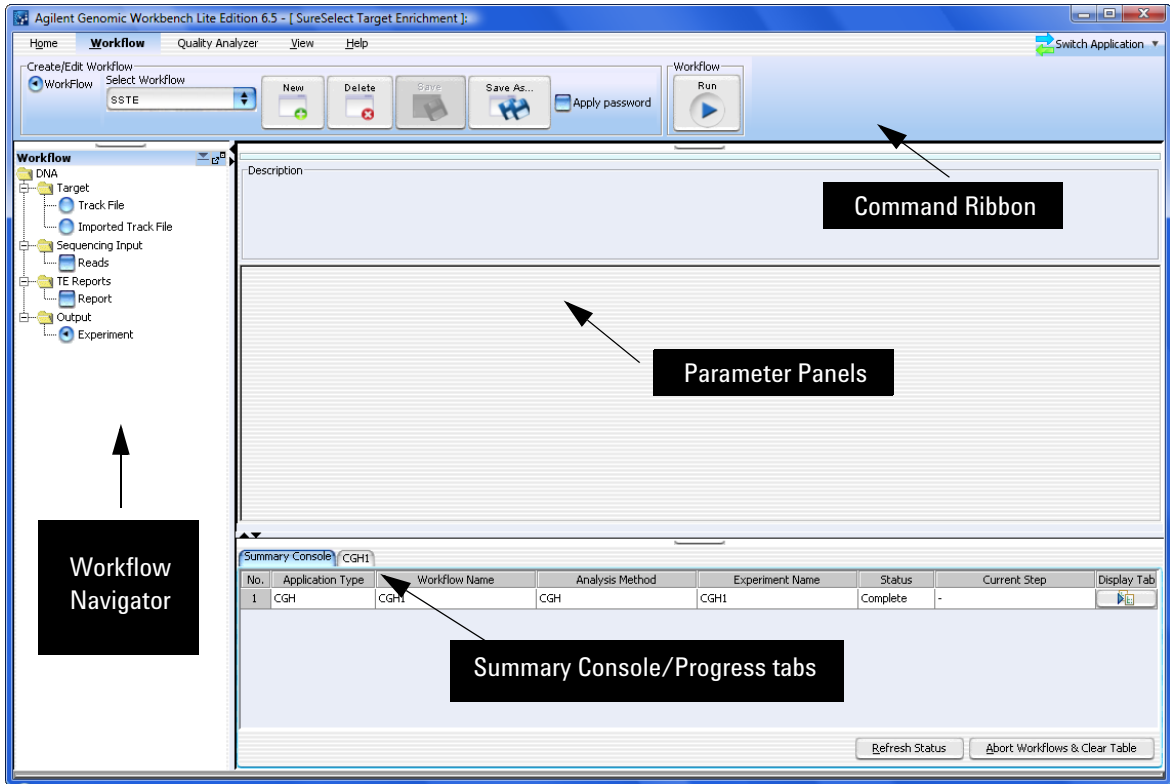


Figure 4 Workflow tab window

Figure 4 shows the tab panes when the Workflow tab is selected. This window contains the elements listed in Table 2 below. More detailed descriptions of each element appear in Chapter 4, “SureSelect Quality Analyzer Reference”.

Table 2 Workflow tab window elements

Element	Purpose
Command ribbon	Contains the commands for setting up, editing, deleting, saving, and running workflows.
Tab pane – Workflow Navigator	Displays current workflow and options. As you select each option, the relevant parameters appear in the parameter panel.
Tab pane – Parameter Panels	Displays parameters for the selection in the Workflow Navigator.
Tab pane – Summary Console/ Progress tabs	For Workflow, this view contains the following tabs: <ul style="list-style-type: none">• Summary Console tab displays the application type, workflow name, experiment name, status, and current step for each workflow completed or in process.• Workflow Progress tab for each named workflow displays the status of the workflow progress, including any errors that might have occurred.

NOTE

In Workflow, the program only lets you set up workflow procedural steps to produce reports, and not explore the data. You must switch to the Quality Analyzer tab to display data and perform further analysis.

Example Exercise 2 – Workflow

In this example exercise, you set up a workflow to automatically perform the three primary steps in the process of producing QC reports on sequencing data. Then you run the workflow.

Before you start the example exercise

- Download the example files for this exercise. Follow these steps:
 - a Go to earray.chem.agilent.com.
The login page of the eArray Web site appears. You do not need to log in.
 - b Under Additional Information, click **Download Agilent Genomic Workbench**.
A page of available Agilent Genomic Workbench files appears.
 - c Under Sample Data, next to Sample files for SureSelect Target Enrichment Quality Analyzer, right-click **Download**, then click **Save Target As**.
 - d Save the file to your desktop.
The eArray Web site transfers a zip format file to your computer.
 - e Extract the zip file to your desktop.
The folder of example files appears on your desktop. Do not extract the zip format files within this folder.
- Familiarize yourself with the parts of the Workflow interface. See “[To start the Workflow program](#)” on page 27 and “[Workflow Tab Window](#)” on page 100.
- Note that the main steps in the example exercises appear in the first column of the table. More detailed instructions appear in the second column. Additional notes and comments appear in the third column.

Step 1. Set up the workflow

You set up a workflow with a similar flow to the one you use to analyze your read files interactively. First, you set up to import the read files, chromosome mapping file and target interval file you need for the QC report and enrichment analyses. Next, you select the report files you intend to generate. Finally, you can choose to rename the experiment that will be automatically created to contain the results of the workflow run.

Example Exercise 2 – Set up the workflow

Step	Detailed Instructions	Notes/Comments
1 Start the workflow for SureSelect Quality Analyzer.	<p>a Start Agilent Genomic Workbench Lite Edition 6.5. The main window of Agilent Genomic Workbench appears.</p> <p>b Click Switch Application > SureSelect Target Enrichment. SureSelect Target Enrichment appears in Application Type. The appropriate tools and panes appear in the main window of the program.</p> <p>c Click Workflow. The Workflow window appears.</p>	<ul style="list-style-type: none"> If upon starting the program you see unfamiliar dialog boxes and don't know how to proceed, please refer to "To start the Agilent Genomic Workbench software" in the <i>Installation Guide</i>. You can use SureSelect Target Enrichment application free of charge. You do not need to purchase or enter a license.
2 Create a new workflow called "xxx Example Workflow," where "xxx" are your initials. <ul style="list-style-type: none"> Do not apply a password. 	<p>a Click New. A dialog box appears.</p> <p>b In Enter Workflow Name, type the name, "xxx Example Workflow".</p> <p>c Since you do not intend to apply a password, click OK.</p> <p>d Set up the workflow in the next steps.</p>	<ul style="list-style-type: none"> If you intend to restrict access to this workflow, follow these steps: <ol style="list-style-type: none"> Click Apply Password. Type in password. Click OK.

1 Getting Started

Step 1. Set up the workflow

Example Exercise 2 – Set up the workflow

Step	Detailed Instructions	Notes/Comments
3 Select target input. <ul style="list-style-type: none"> Select the imported track file (Example Targets from previous exercise) if it has already been imported, OR Select to import the track file in the downloaded samples file folder. <ul style="list-style-type: none"> Build Name: hg18 Track Name: xxx Example Targets, where “xxx” are your initials Track File: *.bed file in downloaded sample files folder 	<ul style="list-style-type: none"> In the Workflow Navigator, open the Target folder. <p>If the file has been imported, follow these steps.</p> <ol style="list-style-type: none"> Click Imported Track File. A parameter panel appears. Next to Imported Track(s), click the arrow and select from the list of imported tracks (Example Targets). <p>If the file SureSelect_All_Exon_G3362_with_names.bed has not been imported, follow the next steps.</p> <ol style="list-style-type: none"> Click Track file. A parameter panel appears. Click the arrows next to Build Name and select hg18. Type Track Name as xxx Example Targets, where “xxx” are your initials. Click Browse, locate the file and click Open. 	<ul style="list-style-type: none"> In this step, you import a BED format track file that contains the desired target genomic intervals. When you select a track file to import, the program only displays *.bed format files. If you worked with the interactive Example Exercise, you have already imported the file. See “Example Exercise 1 – Interactive” on page 20.
4 Enter Read info. <ul style="list-style-type: none"> Use this information: <ul style="list-style-type: none"> File Type: Illumina GA Export Read Length: 76 File Type (2nd one): Single Type Species: H. sapiens Genome Build: hg18 	<ol style="list-style-type: none"> In the Workflow Navigator, open the Sequencing Input folder. Mark Reads. In Read Info next to File Type, select Illumina GA Export, if necessary. Type in a Read Length of 76. For File Type, (2nd one), click Single Type if not selected. 	<ul style="list-style-type: none"> Species and Genome Build should already be filled in with the correct information.

Example Exercise 2 – Set up the workflow

Step	Detailed Instructions	Notes/Comments
5 Select the sequence file. <ul style="list-style-type: none"> Select the imported read file if it has already been imported, or Select to import the read file and chromosome mapping file in the downloaded samples file folder. <ul style="list-style-type: none"> Read Name: xxx s11exporttop Mapping File: chrMapping.txt Sequence File: s_1_1_export_top.zip 	<p>If the file s_1_1_export_top.zip has been imported, follow these steps.</p> <ol style="list-style-type: none"> Click Select already imported Read File. Next to Select Reads, click the arrow and select this file from the list. <p>If the file has not been imported, follow these steps.</p> <ol style="list-style-type: none"> Click Import new Read File. Type the Read Name in the first column for the display. Next to Chr Mapping File, click Browse. In the folder of sample files, select the file chrMapping.txt, then click Open. Next to Sequence File, click Browse. In the folder of sample files, select file s_1_1_export_top.zip and click Open. 	<ul style="list-style-type: none"> In this step, you use these files: <ul style="list-style-type: none"> Sequence file – Contains header information, and the sequence data to be analyzed. The program supports read files in several formats. See “To import sequence read files” on page 38. Chromosome mapping file – A *.txt file that associates the names of chromosomes in the read file with the names of chromosomes as they are specified in the selected genome build.
6 Select these report files and a report location: <ul style="list-style-type: none"> QC Report Processed Intervals Report 	<ol style="list-style-type: none"> Open the TE Reports folder. Mark QC Report. Mark Processed Intervals Report. Click Browse and select a location to store the report. 	<ul style="list-style-type: none"> Do not clear the Overwrite if already exists check box. You can select only two reports at a time.
7 Enter a description for the workflow experiment.	<ol style="list-style-type: none"> Open the Output folder. Click Experiment. Type an Experiment Description if you want. 	<ul style="list-style-type: none"> The name of the workflow automatically populates Experiment Name. You have the option to change it. To view the reports generated during a workflow run, see “To view and save reports” on page 56.
8 Save the workflow.	<ul style="list-style-type: none"> Click Save. <p>Notice the Run button is activated.</p>	

1 Getting Started

Step 2. Run the workflow and review results

Step 2. Run the workflow and review results

Now you are ready to run and monitor the workflow, and review results in the Quality Analyzer window and in the reports generated. During the run, the program analyzes the sequence data against the intervals.

Example Exercise 2 – Run the workflow and review results

Step	Detailed Instructions	Notes/Comments
1 Run the workflow. <ul style="list-style-type: none">• Use “xxx Example Workflow 1” as the name for the run, where “xxx” are your initials.• This name will be the name of the workflow progress tab in the Summary Console pane.	<ul style="list-style-type: none">a On the Workflow ribbon, click Run.b Change the name that is there.c Click OK.	<ul style="list-style-type: none">• After the workflow starts, an experiment is created in the Experiments pane of the Navigator in the Quality Analyzer window.• The Summary Console tab at the bottom of the window displays the status of the current workflow(s).
2 Monitor the workflow run. <ul style="list-style-type: none">• Check the “xxx Example Workflow 1” tab.• Check the Summary Console tab.	<ul style="list-style-type: none">a In the workflow progress tabs, click the “xxx Example Workflow” tab.b Review the ongoing operations in the tab.c Click << Summary Console to return to a summary table of the workflow and experiment.	<ul style="list-style-type: none">• A workflow progress tab is created for the named workflow, which displays the workflow progress and any errors that occur.• Once a workflow has started, you can set up and run additional workflows, which are run in the order you started them. Each workflow will have its own workflow progress tab. See “To run a series of workflows” on page 70.
3 Display results of the workflow both as reports and in the Genomic Viewer.	<ul style="list-style-type: none">a Go to the location you specified for the reports you selected to be generated.b Open the files. Or, if you want to see the results graphically:<ul style="list-style-type: none">a Click Quality Analyzer.b Select the xxx Example Workflow experiment.c Follow the instructions for Example Exercise 1, Step 2, steps 5 and 6.	<ul style="list-style-type: none">• If you want to see enrichment results in the UCSC Genomic Browser, you must mark the WIG file check box before you start the workflow.

Getting Help

To get help within Agilent Genomic Workbench

Help guides are opened in Adobe® Reader® software. Agilent Genomic Workbench has several help resources:

Help Resource	Description/Instructions
SureSelect Quality Analyzer User Guide	<p>This user guide, which you are now reading, supplies comprehensive help on all available Data Viewing tasks. You can access it easily from anywhere within the program.</p> <ol style="list-style-type: none">1 In any tab of Agilent Genomic Workbench, click the Help tab.2 On the Help Ribbon, click Workflow. The SureSelect Quality Analyzer User Guide opens.
Other User Guides	<p>The Help tab in Agilent Genomic Workbench lets you view any of the available user guides that apply to the currently selected application type.</p> <ol style="list-style-type: none">1 Set the desired application type from the Switch Application menu.2 In the Agilent Genomic Workbench tab bar, click Help. The names of the available user guides appear in the command ribbon.3 Click the desired help guide. The selected guide opens.

To contact Agilent Technical Support

Technical support is available by phone and/or e-mail. A variety of useful information is also available on the Agilent Technical Support Web site.

Resource	To find technical support contact information
Agilent Technical Support Web site	<ol style="list-style-type: none">1 Go to http://chem.agilent.com.2 Select a country or area.3 Under Quick Links, select Technical Support.4 Select from the available links to display support information.
Contact Agilent Technical Support by telephone or e-mail (United States and Canada)	Telephone: (800-227-9770) E-mail: informatics_support@agilent.com
Contact Agilent Technical Support by telephone or e-mail (for your country)	<ol style="list-style-type: none">1 Go to http://chem.agilent.com.2 Select Contact Us.3 Under Worldwide Sales and Support Phone Assistance, click to select a country, and then click Go. Complete e-mail and telephone contact information for your country is displayed.

To learn about Agilent products and services

To view information about the Life Sciences and Chemical Analysis products and services that are available from Agilent, go to www.chem.agilent.com.



2 Using SureSelect Quality Analyzer Interactively

Importing Files	38
Setting Up QC Experiments	44
Creating Target Enrichment QC Reports	55

This chapter gives detailed instructions on how to assess the effectiveness of the pull-down of DNA fragments when you use the SureSelect Target Enrichment system. It describes how to import the necessary files into the program, and how to organize the files into *experiments* for analysis. It also describes how to create and export SureSelect QC reports of several different types. In addition, it describes how to use the Genomic Viewer in the Quality Analyzer tab panes to display reads and target intervals in the context of your genome of interest.



Importing Files

You must import a sequence read file, a chromosome mapping file, and a target intervals file before the program can calculate QC metrics and analyze read depth and enrichment. If the applicable genome build for the species of interest is not available in the program, you must also import genome build files. Refer to these topics in this section:

Type of file	See this topic
Sequence read	"To import sequence read files" on page 38
Chromosome mapping	"To import sequence read files" on page 38
Target intervals	"To import target genomic regions as a track" on page 41
Genome build	"To import a genome build" on page 43

To import sequence read files

Sequence read files are gzipped or *.zip archives that contain the base sequences of the fragments in your enriched DNA sample, along with related header information. The program supports these read file formats:

File Type	Comments
Illumina GA Export	Revised version of the ELAND file format that contains additional fields, including quality scores for all bases. You cannot import a multiplexed read file into Agilent Genomic Workbench. You must first separate the file into individual indexed sample files, then import these.
AB SOLiD ma	Color space *.ma file produced by the Applied Biosystems SOLiD system. The AB SOLiD alignment software produces this CSFASTA format file, and places sequence alignment information in the FASTA headers. Color calls are converted to bases when the file is imported.

NOTE

You can import sequence read files in the gzipped or *.zip formats, but not in the .zipx format. In addition, Agilent Genomic Workbench cannot handle a compressed file whose uncompressed file is greater than 4G. In this case, import the uncompressed file directly.

In addition, you must have a *.txt file that maps the names of chromosomes in your sequence read file to the names of the chromosomes in the desired genome build of the species of interest.

- 1 Set the application type to **SureSelect Target Enrichment**. See [“To start the SureSelect Target Enrichment application”](#) on page 14.
- 2 In the Quality Analyzer tab, click **Import > Reads**.
A dialog box appears.
- 3 Select the desired read file, then click **Open**.
The Specify Input Information dialog box appears. See [“Specify Input Information”](#) on page 136.
- 4 Enter the following information:

Group	Instructions/Details
Read Info	<p>The program displays the name of your read file, and the Seq (sequence read file) input type.</p> <ul style="list-style-type: none"> In File Type, select the specific file format that best represents the data in the file.
Species Info	<p>Select the most appropriate species and genome build. The sequence data in your imported read file must be linked to a specific species and genome build.</p> <p>If the desired genome build does not appear, you must import it. See “To import a genome build” on page 43.</p>

2 Using SureSelect Quality Analyzer Interactively

To import sequence read files

Group	Instructions/Details						
Map Chromosomes	<p>You must import a file that maps the names of chromosomes in the read file to the names of chromosomes that are defined in the selected genome build.</p> <p>a Create a tab-delimited text file that contains the desired mapping information. Enter one chromosome per line. Give the name of the chromosome in the read file followed by the name of the matching chromosome in the genome build, separated by a tab character. Do not include a header row.</p> <p>Example file content:</p> <table><tr><td>chrA.fa</td><td>chrA</td></tr><tr><td>chrB.fa</td><td>chrB</td></tr><tr><td>chrC.fa</td><td>chrC</td></tr></table> <p>b In Chr Mapping File, click Browse. An Open dialog box appears.</p> <p>c Select the desired chromosome mapping file, then click Open. The name of the selected file appears in Chr Mapping File.</p>	chrA.fa	chrA	chrB.fa	chrB	chrC.fa	chrC
chrA.fa	chrA						
chrB.fa	chrB						
chrC.fa	chrC						

5 Click **Import**.

The program imports the read file and the chromosome mapping file. A progress bar appears.

Before you can analyze your sequence data, you must also import a BED format track file that defines the target genomic regions. See [“To import target genomic regions as a track”](#) on page 41.

To import target genomic regions as a track

To analyze the sequence data in a read file, the program must correlate reads with target genomic intervals. Typically, these intervals are those to which you have designed your SureSelect Target Enrichment baits. To define the target intervals, you import them as a BED format track file. If you used eArray to create baits, you can download the library as a BED format file, and use this file directly in SureSelect Quality Analyzer. You can also create the file manually.

1 Set the application type to **SureSelect Target Enrichment**. See [“To start the SureSelect Target Enrichment application”](#) on page 14.

2 In the Quality Analyzer tab, click **Import > Track**.

The Import Track dialog box appears. See [“Import Track”](#) on page 131.

3 Enter the following information:

Parameter	Instructions/Details
Species	Select the species that is associated with the genomic intervals in your BED format track file.
Build Name	Select the genome build that is associated with the genomic intervals in your BED format track file.
Color	<p>(Optional) To customize the display color of the intervals in the target intervals track, follow these steps:</p> <ul style="list-style-type: none"> a Click Change. The Choose Track Color Dialog box appears. b In the Swatches tab, select the desired color, then click OK. You can also use the HSB and RGB tabs to define or adjust the color using the HSB and RGB color scales, respectively. See “Choose Track Color” on page 120.
Track Name	Type a name for the track as you want it to appear in reports and lists, and in Gene View of the Genomic Viewer.
Track File	<p>The track file must be a BED format track file that contains the target genomic intervals.</p> <ul style="list-style-type: none"> a Click Browse. An Open dialog box appears. b Select the desired *.bed track file, then click Open. The name of the selected file appears in Track File.

2 Using SureSelect Quality Analyzer Interactively

To import target genomic regions as a track

4 Click OK.

A dialog box tells you that the track was imported successfully.

5 Click OK.

NOTE

- The name, species, genome build, description, and display color of the track can only be entered as you import the track. You cannot edit this information later.
 - You can combine tracks to create a single merged track. Import the individual tracks, then export all of the desired tracks as a single *.zip file. See [“To export tracks”](#) on page 50. You can then import this merged file as the target intervals file.
-

To import a genome build

If you want to import a read file, but the genome build that applies to the reads is not available in the program, you must import the relevant Agilent-supplied genome build files before you import the read file.

1 In the **Home** tab, click **Import > Genome Build**.

The Import Genome Build dialog box appears. See “[Import Genome Build](#)” on page 130.

2 Enter the following information:

Item	Instructions/Details
Species	Type the name of the species to which the imported genome build applies.
Build Name	Type the name of the genome build, as you would like it to appear in lists, dialog boxes, and the like.
Refseq file	This file contains the names and locations of genes in the genome of the selected species. a Click Browse . b Select the desired Agilent-supplied Refseq file, then click Open .
Cytoband file	This file contains the names and locations of cytobands in the genome of the selected species. a Click Browse . b Select the desired Agilent-supplied cytoband file, then click Open .

3 Click **OK**.

A dialog box tells you that the genome build has been successfully imported.

4 Click **OK**.

CAUTION

Use only Agilent-supplied genome build files. You can download these files from the eArray Web site. For more information, see the online help on the eArray Web site.

Setting Up QC Experiments

To create QC reports on your sequence read data, you must set up an experiment. An experiment is a required level of organization that links a specific read file with a specific set of genomic targets.

To create a new experiment

An experiment links a specific sequence read file with the desired target genomic intervals. Target genomic intervals are defined in a BED format track file. Before you can create an experiment, you must import the sequence read file and the track file into the program. See [“To import sequence read files”](#) on page 38 and [“To import target genomic regions as a track”](#) on page 41.

- 1 Set the application type to **SureSelect Target Enrichment**. See [“To start the SureSelect Target Enrichment application”](#) on page 14.
- 2 In the Experiment pane of the Navigator, right-click the **Experiments** node (or folder), then click **New Experiment**.

The Create Experiment dialog box appears. See [“Create Experiment”](#) on page 124.

- 3 Enter the following information:

Parameter	Instructions/Details
Name	Type a name for the experiment. The program uses this name to identify the experiment in the Experiment pane of the Navigator.
Description	(Optional) Type a description for the experiment.
Read Length	Type the number of bases in each read in your sequence read file. The program assumes all reads in the file have the same length.
Targets	Select the track that contains the definitions of the desired genomic targets. If the desired track does not appear in the list, you must import it. Close the dialog box, then see “To import target genomic regions as a track” on page 41.

Parameter	Instructions/Details
SEQ Sample	Select the desired read file. If the desired read file does not appear in the list, you must import it. Close the dialog box, then see “To import sequence read files” on page 38.
Paired	If you have two output files in which the reads are paired, mark this option, then select the paired file. The program imports the two files separately, then merges them before analysis.

4 Click OK.

A progress bar appears. The program adds the experiment to the Experiment pane of the Navigator. To view the experiment, and to create QC reports, you must select the experiment. See [“To select an experiment for analysis”](#) on page 46.

To select an experiment for analysis

To view and analyze the reads in an experiment, and to create QC reports, you must first select the desired experiment to make it active. Only one experiment can be active at a time.

1 Set the application type to **SureSelect Target Enrichment**. See [“To start the SureSelect Target Enrichment application”](#) on page 14.

2 Click **Quality Analyzer**.

SureSelect Quality Analyzer tab appears.

3 In the Experiments pane of the Navigator, in the Experiments folder, right-click the name of the desired experiment, then click **Select Experiment**.

A Confirm dialog box asks if you want to select the experiment.

4 Click **Yes**.

A progress bar appears. The program activates the experiment. In the Experiment pane of the Navigator, the name of the selected experiment appears in blue.

You can now also create and view reports based on the reads in the experiment. See [“Creating Target Enrichment QC Reports”](#) on page 55.

After you select an experiment, the reads in the experiment can appear in all views of the Genomic Viewer. See [“To view the data in an experiment”](#) on page 47.

To deselect an experiment

When you deselect an experiment, its reads no longer appear in the Genomic Viewer, and it is no longer activated for analysis and report creation. However, no files are deleted.

Do one of the following:

- In the Experiment pane of the Navigator, right-click the experiment that you want to deselect, then click **Deselect**.
- Select another experiment. See [“To select an experiment for analysis”](#) on page 46.

To view the data in an experiment

After you select an experiment, the reads in the experiment can appear in all views of the Genomic Viewer.

- 1 Set the application type to **SureSelect Target Enrichment**. See [“To start the SureSelect Target Enrichment application”](#) on page 14.
- 2 In the View tab, under Custom Data, mark **Show**.
- 3 Select the desired experiment. See [“To select an experiment for analysis”](#) on page 46.

The sequence reads from the experiment appear in all views of the Genomic Viewer. [Figure 5](#) shows the Genomic Viewer with reads on human chromosomes 3, 6, and 11.

2 Using SureSelect Quality Analyzer Interactively
To view the data in an experiment

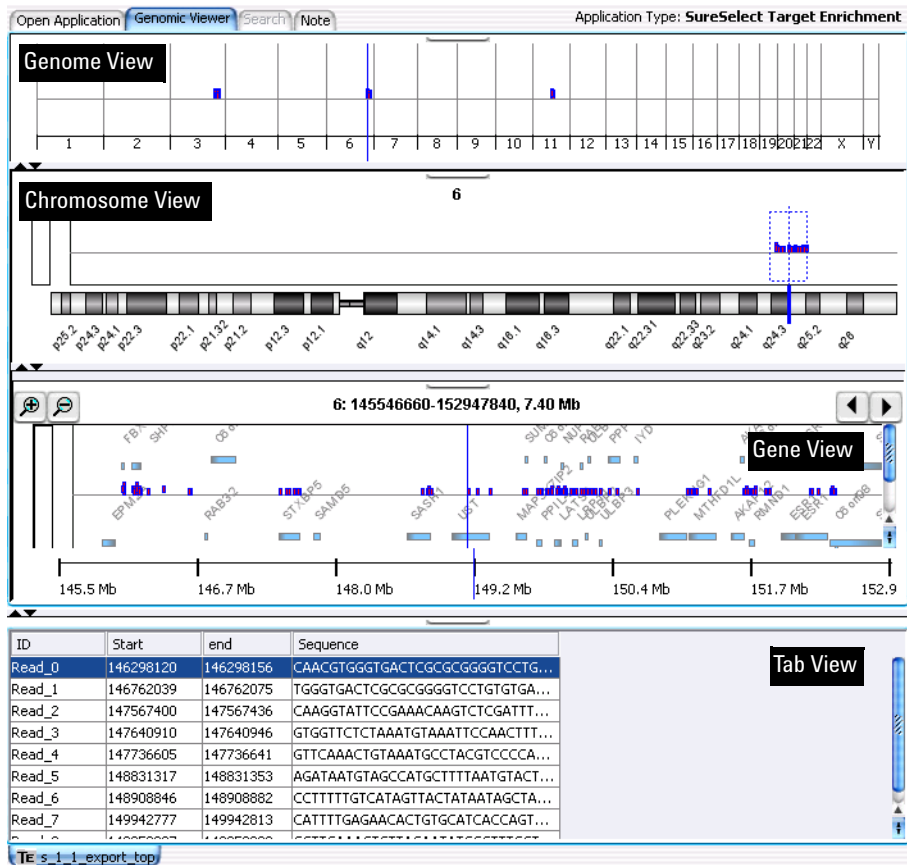






Figure 5 Genomic Viewer. with sequence reads from an experiment

The following read information appears in the Genomic Viewer:

View	Description/Instructions
Genome View	<p>Shows all of the chromosomes of the organism. For chromosomes that have reads associated with them, red/blue bars appear at the approximate loci of reads. When you select a chromosome, its reads appear in all of the other views.</p> <p>Also, to view the results of enrichment analysis in the UCSC Genome Browser, you must first select a chromosome in Genome View. The program subsequently transfers results only for the selected chromosome.</p> <ul style="list-style-type: none"> To select a chromosome, click anywhere within its boundaries. <p>See “Genomic Viewer – Genome View” on page 93.</p>
Chromosome View	<p>Shows the chromosome that was selected in Genome View. A diagram of the selected chromosome appears. Red/Blue bars appear above this diagram that show the approximate locations of reads.</p> <ul style="list-style-type: none"> In the area where the reads appear, drag the pointer through a region, or click a specific location. The region appears in Gene View in greater detail. <p>See “Genomic Viewer – Chromosome View” on page 95.</p>
Gene View	<p>Shows the selected region of the chromosome.</p> <p>You can adjust Gene View in several ways:</p> <ul style="list-style-type: none"> To re-center Gene View on a specific read or gene, click the desired read or gene. To zoom in, click . To zoom out, click . To scroll the view, click  or . <p>See “Genomic Viewer – Gene View” on page 96.</p>
Tab View	<p>Displays a list of the reads that appear in Gene View.</p> <p>See “Genomic Viewer – Tab View” on page 98.</p>

For further instructions on how to use the Genomic Viewer, see the *Data Viewing User Guide*. To open this guide, click **Help > Data Viewing**.

To view the intervals in a target interval track

After you import a set of target intervals as a track, and link the track and a read file in an experiment, you can view the list of intervals in it.

- 1 In the Experiment pane of the Navigator, in the Experiments folder, double-click the name of an experiment that contains the desired target intervals track.
- 2 Right-click the name of the target intervals track, then click **View Details**.

The Track dialog box appears, with header information and a list of intervals. See [“Track”](#) on page 139.

NOTE

You can also use the Genomic Viewer to see the intervals in the genomic region that appears in Gene View. Drag the bottom border of Gene View downward until Gene View is approximately twice its original size. The target intervals track appears at the bottom of gene view.

To export tracks

You can export some or all of the tracks that are available in the program. The program exports all of the intervals in the selected tracks as a single BED format file. This makes it possible to combine tracks. You can then import the combined track and use it in an analysis experiment.

- 1 In the Home tab, click **Export > Tracks**.

The Export Tracks dialog box appears. See [“Export Tracks”](#) on page 128.

- 2 In **Select Tracks**, mark the tracks that you want to export.
- 3 Click **OK**.

An Export dialog box appears.

- 4 Select a location for the exported file, then click **Export**.

The program exports all selected tracks in a single BED format file. You can open the file in a compatible genome browser, and you can also import the BED file into Agilent Genomic Workbench as a new track. See [“To import target genomic regions as a track”](#) on page 41.

To remove a track from the program

If you import a track in error, or if you no longer need one, you can permanently remove it from the program.

- 1 In the **Home** tab, click **User Preferences**.
The Tracks tab of the User Preferences dialog box appears. See “[User Preferences – Tracks tab](#)” on page 144.
- 2 In the list of tracks, in the Delete column, mark the track that you want to delete. You can mark as many tracks as you want.
- 3 Click the **Delete** button.
A dialog box asks if you are sure you want to delete the selected track(s).

CAUTION

When you delete a track, you permanently remove it from the program. To restore a track, you must import it again.

- 4 Click **Yes**.
The program deletes the track.
- 5 Click **OK**.



To customize the appearance of an experiment in the Genomic Viewer



The table below describes options that can be useful when you view sequence reads in the Genomic Viewer:

Customization option	Instructions/Details
To view a different chromosome in Chromosome, Gene, and Tab Views.	<ul style="list-style-type: none">• In Genome View, click the desired chromosome.
To select a specific chromosomal region to view in Gene View	<ul style="list-style-type: none">• In Chromosome View, in the plotting area where reads appear, drag the pointer across the desired region.

2 Using SureSelect Quality Analyzer Interactively

To customize the appearance of an experiment in the Genomic Viewer

Customization option	Instructions/Details
To move the cursor to a specific genomic location	<p>The cursor is a thin blue line that appears in Genome, Chromosome, and Gene Views, and indicates the center of the currently selected region.</p> <p>a In the Home tab, click Go to Gene/Genomic Location. The Go to Gene/Genomic Location dialog box appears. See “Go to Gene/Genomic Location” on page 129.</p> <p>b Do one of the following:</p> <ul style="list-style-type: none"> In RefSeq by Symbol, type a valid RefSeq gene symbol (such as <i>BRCA1</i> or <i>CTSB</i>), then click Go. In Genomic Location, select a chromosome, type a base location, then click Go.
To scroll Gene View	<ul style="list-style-type: none"> Click the scroll buttons at the top of Gene View, as desired. <div style="display: flex; align-items: center; margin-top: 10px;">  – Scrolls Gene View left. </div> <div style="display: flex; align-items: center; margin-top: 10px;">  – Scrolls Gene View right. </div> <p>Note: You can also click anywhere within the gene display area of Gene View to reposition the cursor to that location.</p>
To show only certain chromosomes in Genome View	<ul style="list-style-type: none"> Right-click anywhere within Genome View, then mark or clear the check boxes next to the names of the desired chromosomes. To close the chromosome selection menu, click anywhere outside of the menu.
To show or hide reads in Genome, Chromosome, and Gene Views.	<ul style="list-style-type: none"> In the View tab, under Custom Data, mark or clear Show, as desired.
To show or hide genes in Gene View	<p>a In the Home tab, click User Preferences. The User Preferences dialog box appears. See “User Preferences – Tracks tab” on page 144.</p> <p>b In the Tracks tab, under Visualization Parameters, in Genes, mark or clear Show Gene Symbols in Gene View, as desired.</p>
To show or hide the intervals in the target intervals track	<p>a In the Home tab, click User Preferences. The User Preferences dialog box appears. See “User Preferences – Tracks tab” on page 144.</p> <p>b In the Tracks tab, under Visualization Parameters, in Tracks, mark or clear Show Annotations, as desired.</p> <p>c Click OK.</p>

Customization option	Instructions/Details
To change the width of a column in Tab View.	<p>You can change the width of any column, but it can be especially useful to expand the Sequence column to view the full sequence of reads.</p> <ul style="list-style-type: none"> In the column heading row of Tab View, drag the right edge of the desired column to the left or to the right, as desired.
To increase the size of Gene View	<p>You can expand Gene View to see the target interval track, which can be hidden. Do one of the following:</p> <ul style="list-style-type: none"> If Gene View appears horizontally, drag its bottom border down until you can see the target interval track. If Gene View appears vertically, drag the right border of the main program window to the right, or the left border of Gene View to the left.
To change the zoom level in Gene View	<p>You can increase the zoom level in Gene View to see the boundaries of individual reads, and how they overlap other reads.</p> <ul style="list-style-type: none"> To zoom in, click . To zoom out, click .
To switch between horizontal and vertical orientation	<p>Because of the shape of most monitors, it can be more convenient to use the Genomic Viewer in its default (horizontal) orientation. However, you can use the Genomic Viewer in vertical orientation.</p> <ol style="list-style-type: none"> In the View tab, click View Preferences. The View Preferences dialog box appears. See “View Preferences” on page 147. Under View Alignment, in Orientation, select Horizontal or Vertical, as desired. Click OK.

For additional information about the Genomic Viewer, see the topics referenced in the table above, and the *Data Viewing User Guide*.

To delete an experiment

- 1 In the Experiment pane of the Navigator, right-click the name of the desired experiment, then click **Delete**.

A dialog box asks if you want to delete the experiment.

CAUTION

When you delete an experiment, you permanently remove the link between the particular sequence read file and the target intervals file. To restore the experiment, you must create a new one. However, the program does not delete any data files that you have imported, or any reports that you have saved.

- 2 Click **OK**.

To save an experiment

The program automatically saves experiments as you create them. No additional action is required on your part.

To copy what appears on your screen to the clipboard

You can copy one or all parts of the main program window to the clipboard. You can then paste the copied image into a document in a word processing, graphics, or presentation program.

- 1 In the View tab, click **Copy**, then select the pane that you want to copy. To copy all panes, including the navigator, select **All**.
- 2 In a word processing, graphics, or presentation program, paste the contents of the clipboard into the desired document.

NOTE

This procedure transfers the selected item as a bitmapped image. This image contains only what appears on your screen. It does not contain any hidden panes, or any areas to which you must scroll.

Creating Target Enrichment QC Reports

SureSelect Quality Analyzer can calculate QC metrics and analyze read depth and enrichment, based on the sequence reads in the currently selected experiment.

This section contains the following topics:

- [“To analyze read quality”](#) on page 55
- [“To view and save reports”](#) on page 56

To analyze read quality

After you import read data and target intervals files, and set up and select an experiment, the program can calculate overall QC metrics, and can also analyze read depth and enrichment.

- 1 Set the application type to **SureSelect Target Enrichment**.
- 2 Import files and set up and activate an experiment as described in [“Importing Files”](#) on page 38 and [“Setting Up QC Experiments”](#) on page 44.
- 3 In the Quality Analyzer tab, click **Create Report**.

The program calculates QC metrics and analyzes read depth and enrichment. A dialog box tells you that reports were successfully created.

- 4 Click **OK**.

At this point, the program has created all reports internally. To view a report, you must request a specific report. See [“To view and save reports”](#) on page 56.

To view and save reports

After the program calculates QC metrics and analyzes read depth and enrichment, you can view and save PDF and text format reports. You can also view the results of enrichment analysis in the UCSC Genome Browser.

- 1 Analyze read quality as described in “[To analyze read quality](#)” on page 55.
- 2 View and/or save the desired report as described in the table below:

Report	Description/Instructions
PDF summary report	<p>*.pdf file that contains overall QC metrics. Also contains results of read depth analysis both as tables and as a graph.</p> <p>a In the Quality Analyzer tab, click Summary Report > Save as PDF. A Save dialog box appears.</p> <p>b Select a location and type a name for the report file, then click Save. A dialog box tells you that the report was successfully saved, and asks if you would like to view it.</p> <p>c Click Yes. The report opens in Adobe Reader.</p>
Text summary report	<p>*.txt file that contains overall QC metrics. This report also contains a table of results of read depth analysis.</p> <p>a In the Quality Analyzer tab, click Summary Report > Save as Text. A Save As dialog box appears.</p> <p>b Select a location and type a name for the report file, then click Save As. A dialog box tells you that the report was successfully saved, and asks if you would like to view it.</p> <p>c Click Yes. The report opens.</p>

Report	Description/Instructions
Enrichment analysis results – UCSC	<p>Opens the UCSC Genome Browser and lets you view the results of enrichment analysis for a selected chromosome. See “Enrichment analysis results – UCSC” on page 153.</p> <ol style="list-style-type: none"> In the Genomic Viewer, in Genome View, click anywhere within the boundaries of the desired chromosome. In the command ribbon, click Enrichment > View in UCSC. Your Web browser opens. A message can appear that tells you the browser has blocked scripts and ActiveX controls. If this occurs, allow the blocked content. The UCSC Genome Browser opens. In position/search, type the genomic coordinates of a region that contains reads, then click jump. The enrichment analysis results appear as a bar graph in the User Track area. <p>For additional information on how to use the UCSC Genome Browser, see the online help on the site.</p>
Enrichment analysis results – XLS report	<p>*.xls format file that contains enrichment analysis results for a selected chromosome. See “Enrichment analysis results – XLS report” on page 154.</p> <ol style="list-style-type: none"> In the Genomic Viewer, in Genome View, click anywhere within the boundaries of the desired chromosome. In the command ribbon, click Enrichment > Save as XLS. A Save As dialog box opens. Select a location and type an name for the report file, then click Save As. A dialog box tells you that the report was saved, and asks if you want to view it. Click Yes. The report opens in Microsoft Excel. If Microsoft Excel is not installed on your computer, a dialog box asks you to select a program to open the file. If this dialog box appears, select a spreadsheet program that can open *.xls files.

- For information about QC metrics, see [“QC metrics”](#) on page 150.
- For information about the graph of read depth distribution, see [“Read depth distribution”](#) on page 152.

2 Using SureSelect Quality Analyzer Interactively

To view and save reports



3 Setting Up and Running Workflows

Creating and Managing Workflows [60](#)

Setting Up/Editing Workflows [62](#)

Running Workflows [69](#)

This chapter gives instructions on how to set up and run workflows for producing QC reports sequentially and automatically, using read files, chromosome mapping files and target interval files. These workflows resemble those for the CGH and ChIP applications but do not require analytical methods.



Creating and Managing Workflows

In this section you learn to create, edit, save and delete workflows. For a detailed description of the window of the Workflow tab, see [“Workflow Tab Window”](#) on page 100, and for the command ribbon of the Workflow tab, see [“Command ribbon – Workflow tab”](#) on page 102.

To create a new workflow

- 1** Start the Workflow program.
See [“To start the Workflow program”](#) on page 27.
- 2** Click **New**.
The Create Workflow dialog box appears. See [“Create Workflow”](#) on page 126.
- 3** In Enter Workflow Name, type a name.
If you intend to restrict access to this workflow, mark **Apply Password**.
- 4** Click **OK**.
If you marked the Apply Password check box, the Set Password dialog box appears. See [“Set Password”](#) on page 134.
 - Type a password and click **OK**.
- 5** Set up the workflow.
See [“Setting Up/Editing Workflows”](#) on page 62.
- 6** Click **Save**.

To edit an existing workflow

- 1** Next to the **Select Workflow** list, click the right arrow.
- 2** Select a workflow name from the list.
- 3** Edit the workflow.
See [“Setting Up/Editing Workflows”](#) on page 62.

You can also create a new workflow from a selected one by saving it to another name. See “[To save a workflow to a new name](#)” on page 61.

To save a workflow

- In the command ribbon click **Save**.

To save a workflow to a new name

- 1 Next to the **Select Workflow** list, click the right arrow.
- 2 Select a workflow name from the list.
- 3 In the command ribbon click **Save As**.
- 4 Type the name of the workflow, and click **OK**.

To delete a workflow

- 1 Next to the **Select Workflow** list, click the right arrow.
- 2 Select a workflow name from the list.
- 3 Click **Delete**.

To set a password for an existing workflow

- 1 Next to the **Select Workflow** list, click the right arrow.
- 2 Select a workflow name from the list.
- 3 If you intend to restrict access to this workflow, mark **Apply Password**.
The Set Password dialog box appears. See “[Set Password](#)” on page 134.
- 4 Type a password and click **OK**.

Setting Up/Editing Workflows

This section describes how to set up/edit a workflow for automatic analysis of sequence read files. You can set up or edit a workflow in the Workflow Navigator. Note that the SSQA workflow consists of four parts: Target, Sequencing Input, TE Reports and Output.

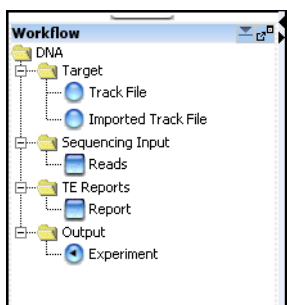


Figure 6 Workflow Navigator

To select target interval input

To analyze the sequence data in a read file, the program must correlate reads with target genomic intervals. Typically, these intervals are those to which you have designed your SureSelect Target Enrichment baits. To define the target intervals for the workflow, you import them as a BED format track file. If you used eArray to create baits, you can download the library as a BED format file.

When you set up a workflow, you must select its source of target interval data.

- In the Workflow Navigator, under Target, click the option button next to the source of target interval data.

See the tables below for a description of the available options.

Table 3 Target Interval Input for SureSelect Quality Analyzer

Option	Description
Track file	Lets you import a track file and use it as a target
Imported track	Lets you select from already imported track files

To import a track file as target input

- 1 Open the Target folder, and click **Track File**.

The Import Track parameter panel appears. See [“Import Track Parameter Panel”](#) on page 107.

- 2 Enter the information in the table below.

Parameter	Instructions/Details
Species	Select the species that is associated with the genomic intervals in your BED format track file.
Build Name	Select the genome build that is associated with the genomic intervals in your BED format track file.

3 Setting Up and Running Workflows

To select target interval input

Parameter	Instructions/Details
Color	(Optional) To customize the display color of the intervals in the target intervals track, follow these steps: a Click Change . The Choose Track Color Dialog box appears. b In the Swatches tab, select the desired color, then click OK . You can also use the HSB and RGB tabs to define or adjust the color using the HSB and RGB color scales, respectively. See “Choose Track Color” on page 120.
Track Name	Type a name for the track as you want it to appear in reports and lists, and in Gene View of the Genomic Viewer.
Track File	The track file must be a BED format track file that contains the target genomic intervals. a Click Browse . An Open dialog box appears. b Select the desired *.bed track file, then click Open . The name of the selected file appears in Track File.

To select an imported track file as target input

- 1 Open the Target folder, and click **Imported Track File**.
The Select Imported Track parameter panel appears. See [“Select Imported Track Parameter Panel”](#) on page 108.
- 2 From the **Imported Track(s)** list, select the imported track whose target intervals you intend to display.

To select sequencing input

Sequence read files are gzipped or *.zip archives that contain the base sequences of the fragments in your enriched DNA sample, along with related header information. The program supports these read file formats:

File Type	Comments
Illumina GA Export	Revised version of the ELAND file format that contains additional fields, including quality scores for all bases. You cannot import a multiplexed read file into Agilent Genomic Workbench. You must first separate the file into individual indexed sample files, then import these.
AB SOLiD ma	Color space *.ma file produced by the Applied Biosystems SOLiD system. The AB SOLiD alignment software produces this CSFASTA format file, and places sequence alignment information in the FASTA headers. Color calls are converted to bases when the file is imported.

In addition, you must have a *.txt file that maps the names of chromosomes in your sequence read file to the names of the chromosomes in the desired genome build of the species of interest.

In Workflow, you have the option of importing a new sequence read file or selecting from the already existing read files.

1 Open the Sequencing Input folder, and mark **Reads**.

The Sequencing Input Parameter Panel appears. See [“Sequencing Input Parameter Panel”](#) on page 109.

2 Enter the following information:

Group	Instructions/Details
Read Info	<ul style="list-style-type: none"> In Species, select the species that is represented in the imported sequence read file. In Build, the available genome builds for the selected species appear in the list. Select the build that applies to the reads in the imported file. If the desired genome build does not appear, you must import it. See “To import a genome build” on page 43. In File Type, select the specific file format that best represents the data in the file. In Read Length, type the number of base pairs in the typical fragment.

3 Setting Up and Running Workflows

To select sequencing input

Group	Instructions/Details
File Type	<ul style="list-style-type: none">Click Simple Type or Pair Type. If you have two output files in which the reads are paired, mark Pair Type, then select the paired file. The program imports the two files separately, then merges them before analysis.
Seq File N	<ul style="list-style-type: none">Click Select already imported Read File, then select the file from the Selected Reads list, ORClick Import new Read File,Click Browse to select the Sequence File.Type a Read Name (display name) for the Sequence File.You must import a file that maps the names of chromosomes in the read file to the names of chromosomes that are defined in the selected genome build.<ul style="list-style-type: none">Create a tab-delimited text file that contains the desired mapping information. Enter one chromosome per line. Give the name of the chromosome in the read file followed by the name of the matching chromosome in the genome build, separated by a tab character. Do not include a header row. <p>Example file content:</p> <pre>chrA.fa chrA chrB.fa chrB chrC.fa chrC</pre> <ul style="list-style-type: none">In Chr Mapping File, click Browse. An Open dialog box appears.Select the desired chromosome mapping file, then click Open. The name of the selected file appears in Chr Mapping File.

NOTE

You can import sequence read files in the gzipped or *.zip formats, but not in the .zipx format. In addition, Agilent Genomic Workbench cannot handle a compressed file whose uncompressed file is greater than 4G. In this case, import the uncompressed file directly.

To select reports

After the workflow program calculates QC metrics and analyzes read depth and enrichment, it automatically saves the information in the formats you have chosen. You can also view the results of enrichment analysis in the UCSC Genome Browser using the workflow-generated WIG file.

- 1 In the Workflow Navigator, under **TE Reports**, mark **Report**.

The Text Summary Report Parameter Panel appears. See [“Text Summary Report Parameter Panel”](#) on page 112.

- 2 Mark the reports you intend to generate.

NOTE

You can select any two reports, but no more than two for any workflow.

- 3 In the parameter panel, click **Browse** and set the location for each report.
- 4 If you do not want to overwrite an existing report, clear **Overwrite if already exists**.

See [Table 4](#) for a description of the available reports.

Table 4 Sure Select Quality Analyzer analysis and enrichment reports

Report	Description/Instructions
Text Summary Report	*.txt file that contains overall QC metrics. This report also contains a table of results of read depth analysis.
QC Report	*.pdf file that contains overall QC metrics. Also contains results of read depth analysis both as tables and as a graph.
Processed Intervals Report	*.xls format file that contains enrichment analysis results for a selected chromosome. See “Enrichment analysis results – XLS report” on page 154. If Microsoft Excel is not installed on your computer, a dialog box asks you to select a program to open the file. If this dialog box appears, select a spreadsheet program that can open *.xls files.

Table 4 Sure Select Quality Analyzer analysis and enrichment reports (continued)

Report	Description/Instructions
WIG file	After the file is generated by the workflow, open it in the UCSC Genome Browser to view the results of enrichment analysis for a selected chromosome. See “Enrichment analysis results – UCSC” on page 153. For additional information on how to use the UCSC Genome Browser, see the online help on the site.

To rename the experiment

When you create and name a workflow, the experiment that will be created when you run the workflow automatically takes on the name of the workflow. You can change its name and description.

- 1 In the Workflow Navigator, click **Experiment**.
The Experiment Parameter Panel appears. See [“Experiment Parameter Panel”](#) on page 106.
- 2 Type a new **Experiment Name**.
- 3 In **Experiment Description** type any new descriptive material you choose.

Running Workflows

When you run a workflow, all of the selected reports are generated automatically, but to view the reads and target intervals graphically, you must use the Genomic View and UCSC Genomic Browser.

To run a workflow

- 1 Create a new workflow or select one from the list.
See [“To create a new workflow”](#) on page 60 or [“To edit an existing workflow”](#) on page 60.
- 2 On the Workflow ribbon, click **Run**.
- 3 At the prompt, type a name into the **Provide Workflow Identifier** dialog box, and then click **OK**.

This name becomes the name of the progress tab in the Summary Console/Progress tabs pane. The default name is the name of the workflow. Entering a new name does not change the name of the experiment or the workflow. See [“Provide Workflow Identifier”](#) on page 133.

An experiment is created in the Experiments pane of the Home tab Navigator. The experiment folder lists the Experiment Name you specified when you set up the workflow.

The Summary Console tab at the bottom of the window in the Summary Console/Progress tabs pane displays the status of the workflow run. See [“Summary Console tab”](#) on page 116.

A workflow progress tab is created in the Summary Console/Progress tabs pane for the named workflow and displays the progress of the run and any errors that occur. A not-yet-completed run has “running” displayed in its progress tab. A completed run has “completed” displayed in its progress tab. See [“Named workflow tabs”](#) on page 118.

NOTE

The Summary Console and progress tabs include workflows for CGH, ChIP and SureSelect Target Enrichment. For information on the CGH and ChIP workflows, see the *Workflow User Guide*.

To run a series of workflows

Once a workflow run has started, you can set up and run additional workflows, which are run in the order in which you started them. Each workflow will have its own progress tab.

On 64-bit computers, you can run up to three workflows simultaneously. On 32-bit computers, workflows are processed sequentially, not simultaneously.

To run a series with different workflows

- 1 Run the first workflow.

Follow the instructions in [“To run a workflow”](#) on page 69.

- 2 Create a new workflow or select an existing one from the list.

See [“To create a new workflow”](#) on page 60 or [“To edit an existing workflow”](#) on page 60.

- 3 In Output, click **Experiment**.

- 4 Change the name of the experiment if the named experiment exists in the Experiment pane.

- 5 Save the workflow.

- 6 Click **Run**.

- 7 Repeat steps 2-5 until you are complete.

See [“To monitor workflow runs”](#) on page 71.


To run a series with the same workflow

- 1 Run the first workflow.

Follow the instructions in [“To run a workflow”](#) on page 69.

- 2 Click **Run** again.

An Input dialog box appears with the name of the experiment in the workflow plus an increment of 1. This will be the name of the experiment created for the second run of the same workflow. You can change the name if you want.

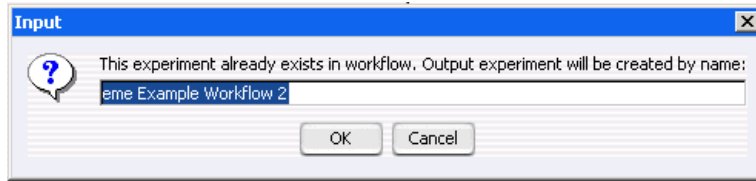

3 Click OK.

The Provide Workflow Identifier dialog box appears. See [“Provide Workflow Identifier”](#) on page 133.

4 Enter the name for the second workflow progress tab, and click OK.

5 Repeat steps 2 through 4 until you are complete.

See [“To monitor workflow runs”](#) on page 71.

The Summary Console tab at the bottom of the window in the Summary Console/Progress tabs pane displays the status of all the in-progress or completed workflow(s). See [“Summary Console tab”](#) on page 116.

To monitor workflow runs

You can monitor workflow runs in the Summary Console/Progress tabs pane by viewing the Summary Console tab and/or the progress tabs.

- Click **Summary Console** to monitor the status of all the workflows.
See [“Summary Console tab”](#) on page 116.
- Click the button under **Display Tab** in the Summary Console row to go to the progress tab for the selected run, or in the progress tabs, click a named workflow tab.

A progress tab displays the workflow progress and any errors that occur. A not-yet-completed workflow run has “running” displayed in its progress tab. A completed workflow run has “completed” displayed in its progress tab. See [“Named workflow tabs”](#) on page 118.

- Click **Close Tab** in a progress tab to stop that run and remove the workflow from the list.
- Click **< Summary Console** to return to a summary table of each workflow and experiment.

3 Setting Up and Running Workflows

To abort a single workflow

- Click **Abort Workflows & Clear Table** in the Summary Console tab to stop all workflow runs, and remove workflows and workflow experiments from the run list.

To abort a single workflow

- 1 In the Summary Console/Progress tabs pane, click the progress tab to abort that workflow.
- 2 Click **Abort Workflow**.

You are asked to confirm that you want to abort the workflow. Note that the Abort Workflow button will not appear unless the workflow is currently running.

The workflow is aborted, and the workflow progress tab and workflow status in the Summary Console tab are removed.

To abort all workflows

- 1 In the Summary Console/Progress tabs pane, click the **Summary Console** tab.
- 2 Click **Abort Workflows & Clear Table**.

You are asked to confirm that you want to abort all workflows.

- 3 Click **Yes**.

A warning that workflow information will be lost is displayed for each workflow.

- 4 Click **Yes** to confirm each.

All workflows are aborted and the Summary Console/Progress tabs pane is cleared of all workflow information.

To display the results of a workflow

After you configure a workflow and run it, the visual results are displayed in the Genomic Viewer pane of the Quality Analyzer window, and the reports are found in the location(s) you specified. You can use other programs to open the reports created by the workflow.

To view the results in Genomic Viewer:

1 Click **Quality Analyzer.**

The workflow results appear as a node under the experiment created during the workflow run.

2 Select the experiment that contains the results.

3 Click **View, and mark **Show** in Custom Data.**

The reads now appear in Genomic Viewer, and the target intervals appear in Gene View below the reads.

To view the reports created during the workflow run:


- Open the report(s) you specified when you set up the workflow.

Result/Report	Do this to open the report.
Text Summary Report	<p>For this report, the program creates a *.txt file that contains a table of QC metrics and the results of read depth analysis. You can open this file with a text editor or word processor.</p> <ol style="list-style-type: none"> 1 Go to the location you specified to contain the Text Summary Report. 2 Open the report with Microsoft Word.
QC Report	<p>For this report, the program creates a PDF format report that contains tables of QC metrics. It also contains the results of read depth analysis shown both as tables and as a graph.</p> <ol style="list-style-type: none"> 1 Go to the location you specified to contain the PDF QC Report. 2 Open the report with Adobe Acrobat Reader.

3 Setting Up and Running Workflows

To display the results of a workflow

Result/Report	Do this to open the report.
Processed Intervals Report	<p>For this report, the program creates an *.xls file that contains enrichment analysis results for all the chromosomes. You can view this file in Microsoft Excel.</p> <ol style="list-style-type: none">1 Go to the location you specified to contain the XLSReport.2 Open the report with Microsoft Excel.
WIG File	<p>For this report, the program creates a track file that contains enrichment analysis results for the entire genome.</p> <ol style="list-style-type: none">1 Right-click Genomic Viewer, and select UCSC Genomic Browser.2 Go to the location you specified to contain the WIG file, and select the file.3 Click Open. <p>The enrichment results appear in the UCSC Genomics Browser.</p>

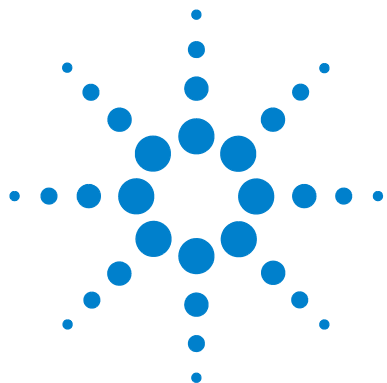

4

SureSelect Quality Analyzer Reference

SureSelect Target Enrichment Main Window 76

Quality Analyzer Tab Window 83

Workflow Tab Window 100

Dialog Boxes 120

Target Enrichment Reports 149

This chapter describes the parts of the Agilent Genomic Workbench Lite Edition main window that are relevant to the SureSelect Target Enrichment application, including descriptions of both the Quality Analyzer tab and Workflow tab windows. It also describes the menus, shortcut menus, panes, and dialog boxes that can appear. In addition, it contains details about the reports the program can produce.

You can find more information on the Agilent Genomic Workbench Lite Edition programs in the *Product Overview Guide*. Click the Open Application tab if it is not already open. Access to the guide is in the top right corner of the Open Application window.

For general help with the Genomic Viewer and Navigator panes of the user interface, see the *Data Viewing User Guide*. In addition, separate user guides cover each of the main programs in Agilent Genomic Workbench Lite Edition. To view these other guides, select the relevant application type, then click the **Help** tab. You can view some of the applications only if you have purchased a license.



SureSelect Target Enrichment Main Window

When you first switch the application to the SureSelect Target Enrichment application, you see this window for the Home tab. The Tab bar and Switch Application button remain the same for every click of a tab. But the Command ribbon changes for each tab, and the Tab panes may or may not stay the same.

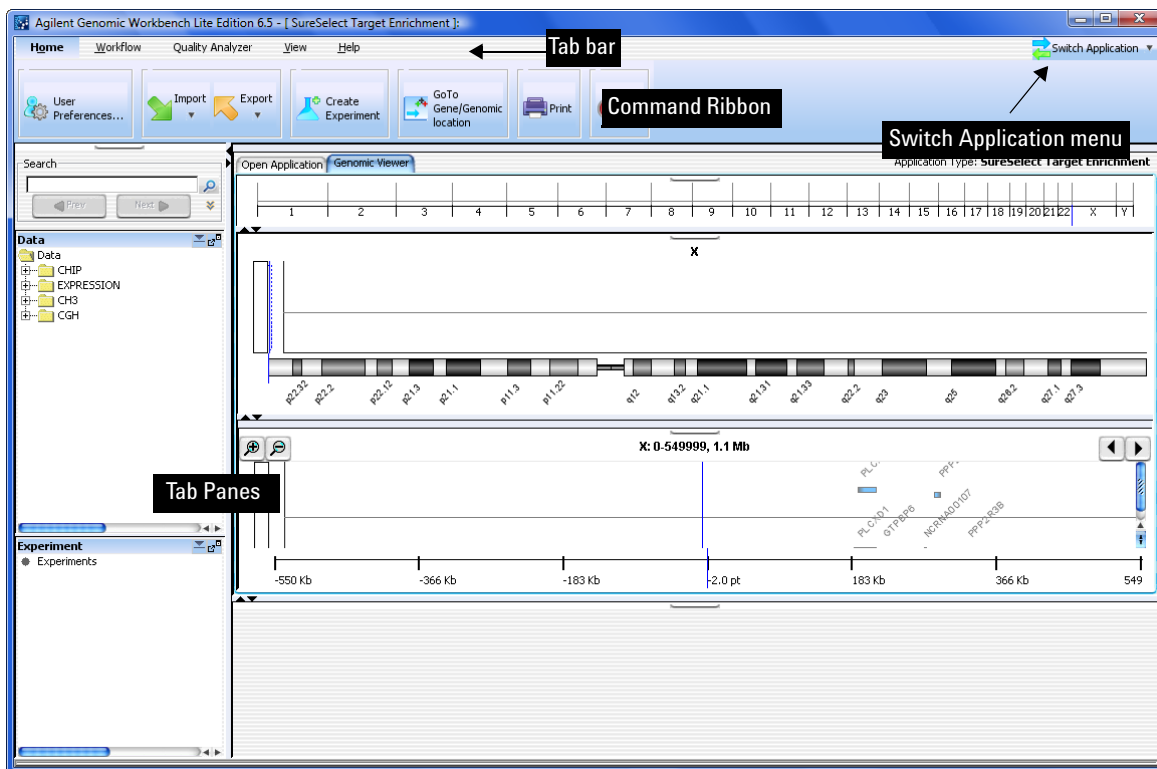


Figure 7 SureSelect Target Enrichment main window – Home tab

The SureSelect Target Enrichment main window has these main elements:

Element	Description
Title bar	Displays the program name and the currently selected application type.
Switch Application menu	Lets you select the experimental application type. This is one of the main settings for Agilent Genomic Workbench, since functionality as well as design, sample, and sequence data are partitioned by application type. See “Switch Application menu” on page 78 and “To start the SureSelect Target Enrichment application” on page 14.
Tab bar	Contains the tabs that let you use the main functions of the SureSelect Target Enrichment program. See “SureSelect Target Enrichment tab bar” on page 79.
Command ribbon	<p>Contains the commands that you use to access the main functionality of the SureSelect Target Enrichment program. A different command ribbon appears for each tab. The Command ribbons for the Quality Analyzer tab window and the Workflow tab window are described in separate sections.</p> <p>See these topics:</p> <ul style="list-style-type: none"> • “Command ribbon – Home tab” on page 80 • “Command ribbon – Help tab” on page 82 • “Quality Analyzer Tab Window” on page 83 • “Workflow Tab Window” on page 100
Tab panes	<p>These change depending on the tab you choose.</p> <p>When you select either the Home, Quality Analyzer, View, Tool or Help tabs, the Navigator and Genomic Viewer panes appear. See “Quality Analyzer Tab Window” on page 83.</p> <p>When you select the Workflow tab, the Workflow Navigator, Parameter Panels and Summary Console panes appear. See “Workflow Tab Window” on page 100.</p>

Switch Application menu

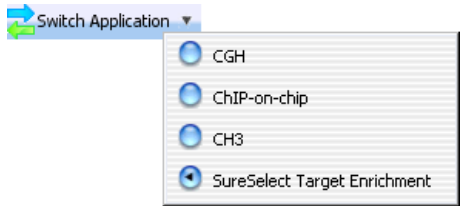


Figure 8 Switch Application menu

Purpose: Lets you select the main experimental application type for the program. Design, sample, and sequence data, as well as program functionality, are partitioned by application type. See [“To start the SureSelect Target Enrichment application”](#) on page 14. To use the Target Enrichment Quality Analyzer, you select **SureSelect Target Enrichment** as the application type.

To open: Click **Switch Application**.

These application types are available:

Application Type	Description
CGH	Data analysis for comparative genomic hybridization studies.
ChIP-on-chip	Data analysis for chromatin immunoprecipitation studies.
CH3	Data analysis for methylation studies.
SureSelect Target Enrichment	Lets you do a QC analysis of target fragment pull-down.

SureSelect Target Enrichment tab bar

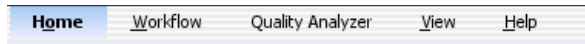


Figure 9 Tab bar for SureSelect Target Enrichment

The tabs that are available in the tab bar vary by application type. To create and view Target Enrichment QC reports, you set the application type to SureSelect Target Enrichment. For this application type, these tabs appear in the tab bar:

Tab	Description
Home	Lets you set user preferences, import and export files, create experiments, view data, genes, and tracks, print certain panes, and exit the program. See “Command ribbon – Home tab” on page 80. Note: To calculate Target Enrichment QC metrics, and to produce reports, you do not need to use this tab. All necessary commands are available in the Quality Analyzer and Workflow tabs.
Workflow	Lets you set up importing files, creating experiments and producing QC reports for automatic processing. See “Command ribbon – Workflow tab” on page 102.
Quality Analyzer	Lets you calculate QC metrics based on selected sequence read data, and analyze read depth and enrichment. You can also create PDF and text QC reports. See “Command ribbon – Quality Analyzer tab” on page 85.
View	Lets you set view preferences, and customize the appearance of the main program window. For details, see the <i>Data Viewing User Guide</i> .
Help	Lets you view the user guides that are relevant to the SureSelect Target Enrichment application type. See “Command ribbon – Help tab” on page 82.

Command ribbon – Home tab



Figure 10 Command ribbon – Home tab for SureSelect Target Enrichment

Purpose: Lets you set user preferences, import and export files, create experiments, view data, genes, and tracks, print certain program panes, and exit the program.

To open: Set the application type to **SureSelect Target Enrichment**, then click **Home**.

User Preferences Opens the User Preferences dialog box, where you can customize the appearance of genes and tracks in the Genomic Viewer, select a default data location, and select other program options. See [“User Preferences – Miscellaneous tab”](#) on page 142 and [“User Preferences – Tracks tab”](#) on page 144.

Import Opens a menu with these options:

Option	Description
Reads	Opens a dialog box where you can import a sequence read file and a chromosome mapping file. See “To import sequence read files” on page 38.
Track	Opens the Import Track dialog box, where you can set up and import a target intervals track file. See “To import target genomic regions as a track” on page 41 and “Import Track” on page 131.
Genome Build	Opens the Import Genome Build dialog box, where you can set up and import a new Agilent-supplied genome build. See “To import a genome build” on page 43 and “Import Genome Build” on page 130.

Export Opens a menu with a Tracks option. This option opens the Export Tracks dialog box, where you can select one or more tracks to export as a single BED format file. See [“To export tracks”](#) on page 50 and [“Export Tracks”](#) on page 128.

Create Experiment	Opens the Create Experiment dialog box, where you can begin the creation of a Target Enrichment QC analysis experiment. Experiments are required organizational units that link a read file with related files and information. See “Create Experiment” on page 124 and “To create a new experiment” on page 44.
Go to Gene/Genomic Location	Opens the Go to Gene/Genomic Location dialog box, where you can enter a gene symbol or genomic location. This lets you move the cursor of the Genomic Viewer to a specific locus. See “To customize the appearance of an experiment in the Genomic Viewer” on page 51 and “Go to Gene/Genomic Location” on page 129.
Print	Opens a Print dialog box, where you can print the contents of several program panes. See the <i>Data Viewing User Guide</i> .
Exit	Exits the program.

Command ribbon – Help tab

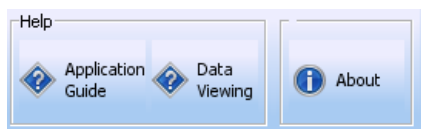


Figure 11 Help tab for SureSelect Target Enrichment

Purpose: Lets you view the user guides for Agilent Genomic Workbench that are relevant to the selected application type.

To open: Set the desired application type, then click **Help**.

Help

All guides open in Adobe Reader. These buttons can appear.

Application Guide

Opens the following User Guide:

- **SureSelect Target Enrichment** – Opens the *SureSelect Quality Analyzer User Guide*. This guide describes how to use the SureSelect Quality Analyzer application (including Workflow) of Agilent Genomic Workbench to assess the effectiveness of fragment pull-down for target enrichment experiments. This application does not require a license.

Data Viewing

Opens the *Data Viewing User Guide*. This guide describes how to import, organize, manage, export, and display data and other content (experiments, gene lists, tracks) within Agilent Genomic Workbench. It is targeted for users who have no DNA Analytics application license(s).

An additional guide is available in the Open Application tab of the program. The *Product Overview Guide* gives an overview of the capabilities within Agilent Genomic Workbench. It also describes how to start each of the component programs and find help, and how to enter your license information. To open this guide, click the **Open Application** tab, then click **Product Overview**.

About

Opens a dialog box that displays title, copyright, and version information for the program.

Quality Analyzer Tab Window

You will find descriptions here of the Quality Analyzer command ribbon, and the Navigator and Genomic Viewer panes, even though the latter appear when you click any of the SureSelect Target Enrichment tabs, except for the Workflow tab.

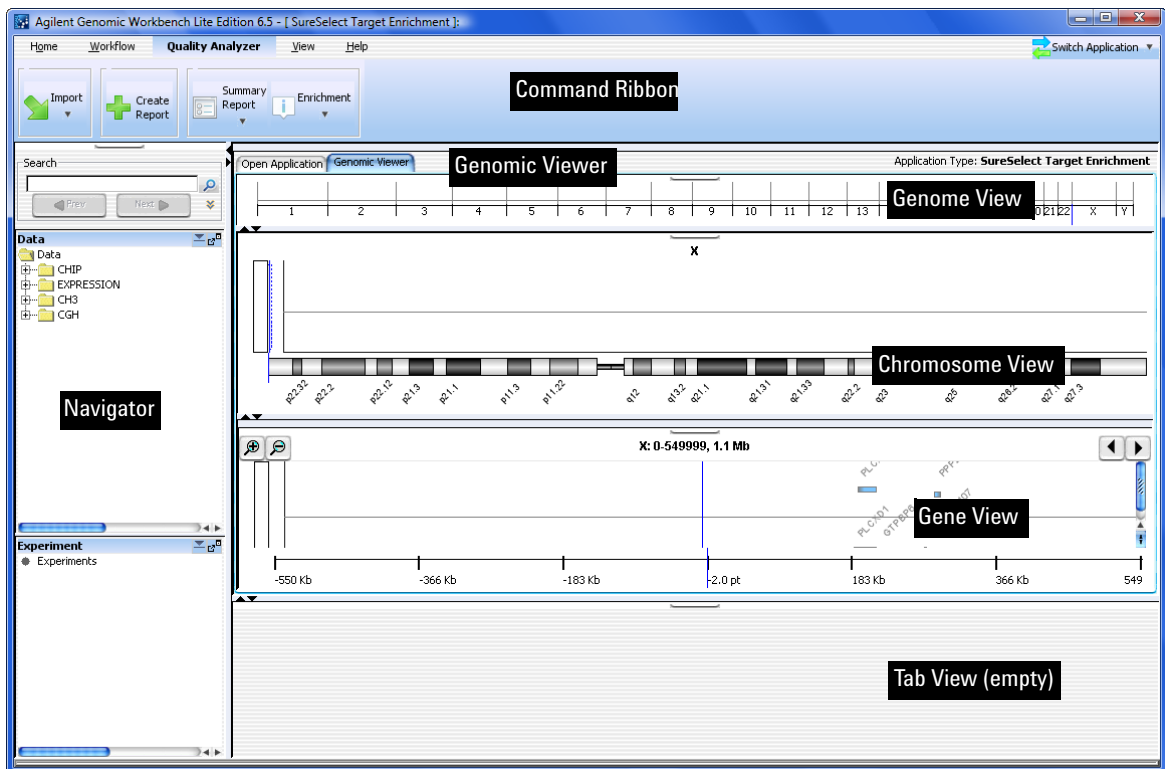


Figure 12 Quality Analyzer tab window

Table 5 Quality Analyzer tab window elements

Window Element	Description
Command ribbon	Contains the commands for importing sequence read files and target interval tracks, doing QC analyses, and creating QC reports. See Appendix 4 , “Command ribbon – Quality Analyzer tab,” starting on page 85.
Tab pane – Navigator	Contains lists of the designs, sample data, sequence data, analysis experiments, and other related items that are available to you. A search feature is also available. See these topics: <ul style="list-style-type: none">• “Navigator overview” on page 87• “Navigator – Experiment pane” on page 88
Tab pane – Genomic Viewer	<p>Shows the genome of the selected organism at several levels of detail. Genes, tracks, probes, baits, and data of many types can appear in the Genomic Viewer in the context of the genome.</p> <p>For general information on how to use the Genomic Viewer, see the <i>Data Viewing User Guide</i>.</p> <p>For information on the specific features of the Genomic Viewer that are relevant to the Quality Analyzer tab, see “Genomic Viewer – overview” on page 91.</p>

Command ribbon – Quality Analyzer tab

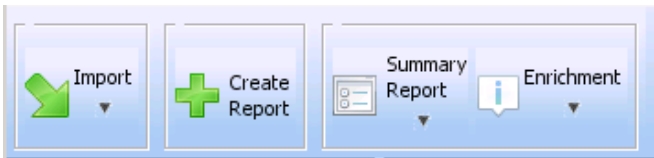


Figure 13 Command ribbon – Quality Analyzer tab

Purpose: Lets you import sequence read files and target interval tracks, do QC analyses, and create QC reports.

To open: Set the application type to **SureSelect Target Enrichment**, then click **Quality Analyzer**.

Import Opens a menu with these options:

Option	Description
Reads	<p>Opens a dialog box where you can import a sequence read file. You must import a gzipped or *.zip file that contains a read file in one of these formats:</p> <ul style="list-style-type: none">• Illumina GA Export – Revised version of the ELAND file format that contains additional fields, including quality scores for all bases. You cannot import a multiplexed read file into Agilent Genomic Workbench. You must first separate the file into individual indexed sample files, then import these.• AB SOLiD ma – Color space *.ma file produced by the Applied Biosystems SOLiD system. The AB SOLiD alignment software produces this CSFASTA format file, and places sequence alignment information in the FASTA headers. Color calls are converted to bases when the file is imported. <p>Note: You cannot import a read file in the .zipx format. See “To import sequence read files” on page 38.</p>
Track	<p>Opens the Import Track dialog box, where you can set up and import a target intervals track. The track must be a *.txt file in BED format. See “To import target genomic regions as a track” on page 41.</p>

Create Report Analyzes the data in the currently selected experiment. To view the results, you use the **Summary Report** and **Enrichment** commands (see below).

Summary Report Opens a menu with these options:

Option	Description
Save as PDF	Opens a Save dialog box, where you can select a location for a PDF QC report file. This report contains calculated QC metrics. It also contains the results of read depth analysis both as tables and as a graph. See these topics: <ul style="list-style-type: none">• “To view and save reports” on page 56• “QC metrics” on page 150• “Read depth distribution” on page 152.
Save as Text	Opens a Save As dialog box where you can select a location for a *.txt format report. This report contains calculated QC metrics. It also contains tables of the results of read depth analysis. See “To view and save reports” on page 56 and “QC metrics” on page 150.

Enrichment Opens a menu with these options:

Option	Description
View in UCSC	Opens the UCSC Genome Browser in your Web browser, and displays the results of enrichment analysis in the User Track area. This command lets you view these results one chromosome at a time. You must select the desired chromosome in Genome View of the Genomic Viewer before you click this command. See “To view and save reports” on page 56 and “Enrichment analysis results – UCSC” on page 153.
Save as XLS	Opens a Save As dialog box, where you can select a location for a *.xls format report. This report contains the results of enrichment analysis for each target interval in the selected chromosome. You select the desired chromosome in Genome View of the Genomic View before you click this command. You can use a spreadsheet program to view this report.

NOTE

To view or save summary or enrichment reports, you must first import a read file and a target intervals file, create and select an analysis experiment, and analyze the data in the selected experiment.

Navigator overview

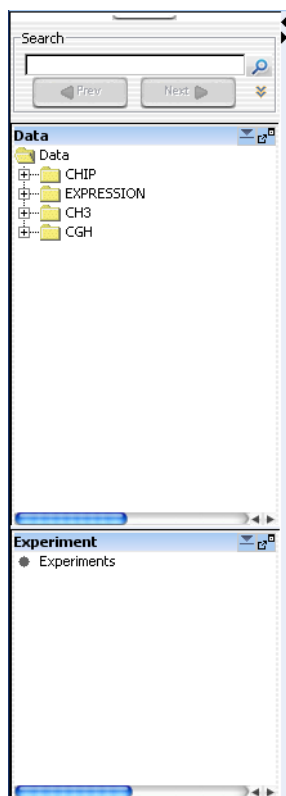


Figure 14 SureSelect Target Enrichment Navigator

For the SureSelect Target Enrichment application type, the Navigator contains several panes that show the content that is available to you:

Pane of Navigator	Description
Search	Lets you search all available panes of the Navigator for items that match a search term. For information about this pane, see the <i>Data Viewing User Guide</i> .

Pane of Navigator	Description
Data	You do not use this pane of the Navigator when you use the Quality Analyzer tab.
Experiment	Lets you create, select, and view experiments for Target Enrichment Quality Analyzer. See “Navigator – Experiment pane” on page 88.

Navigator – Experiment pane











Figure 15 Navigator – Experiment pane

Purpose: In the SureSelect Target Enrichment application type, the Experiment pane of the Navigator lets you create, view, and analyze Target Enrichment QC analysis experiments, which are organizational units that link sequence read data with target intervals and other information. You create an experiment as a prerequisite to QC analysis.

Experiment pane – Icons and special text

These icons and special text items can appear in the Experiment pane of the Navigator:

Item	Description
	(Available if the Experiment pane is not minimized) Minimizes the experiment pane to the bottom of the Navigator.
	(Available if the Experiment pane is minimized) Restores the Experiment pane to its position in the middle of the Navigator.

Item	Description
	If the Experiment pane is docked in the Navigator, this button detaches the Experiment pane from the Navigator and opens it in a new, separate window. If the experiment pane is in its own separate window, this button re-attaches the Experiment pane to the Navigator.
	A collapsed folder whose contents are hidden. Folders can contain data items and/or other folders.
	An expanded folder. Its contents appear below the name of the folder in a hierarchical “tree” format.
	Expands a folder to show its contents.
	Collapses a folder to hide its contents.
	A data node that represents a specific imported read file or target interval track. It also represents the Experiments node (folder) when no experiments are present.
blue text	The name of an experiment that is selected for analysis.
red text	An item that matches the search term in a search of the Navigator.
highlighted text	The currently selected search result from a search of the Navigator.

Experiment pane – Actions and shortcut menus

- Double-click the **Experiments** folder to expand it. Double-click it again to collapse it.
- Right-click the **Experiments** folder to open a shortcut menu with a New Experiment option. This option opens the Create Experiment dialog box, where you can create a new QC analysis experiment. See “[Create Experiment](#)” on page 124 and “[To create a new experiment](#)” on page 44.

- Right-click the name of an experiment to open a shortcut menu with these options:

Option	Description
Select Experiment	(Available for experiments other than the currently active one) Lets you view and analyze the data in the experiment. The reads and target intervals that are defined for the selected experiment can appear in the Genomic Viewer, and are used for QC analysis. When you select an experiment, the program displays the name of the experiment in blue. You can only select one experiment at a time. See “To select an experiment for analysis” on page 46.
Deselect Experiment	(Available for the currently selected experiment, if any) Removes the reads and target intervals that are defined in the experiment from the Genomic Viewer, and makes them unavailable for QC analysis. See “To deselect an experiment” on page 47. If you deselect an experiment, you can select it again as needed.
Delete Experiment	Removes the experiment from the Experiment pane of the Navigator. To restore a deleted experiment, you must create a new experiment. See “To delete an experiment” on page 54.

- Double-click the name of an experiment to view its contents. Double-click the name of the experiment again to hide its contents. Each experiment contains a genome build folder that contains a link to the uploaded read file that has been selected for the experiment. It also contains a link to the target intervals track that has been selected for the experiment.
- Right-click the name of a target intervals file to open a shortcut menu with a View Details option. This option opens the Track dialog box, where you can view a list of the target intervals in the track, and header information. See [“To view the intervals in a target interval track”](#) on page 50 and [“Track”](#) on page 139.

Genomic Viewer – overview

This section describes the panes that appear in the Genomic Viewer, and the commands, options, and shortcut menus that are available. These descriptions are specific to the SureSelect Target Enrichment application type. Other applications, such as the CGH, ChIP-on-chip, and CH3 data analysis applications, display data and results differently.

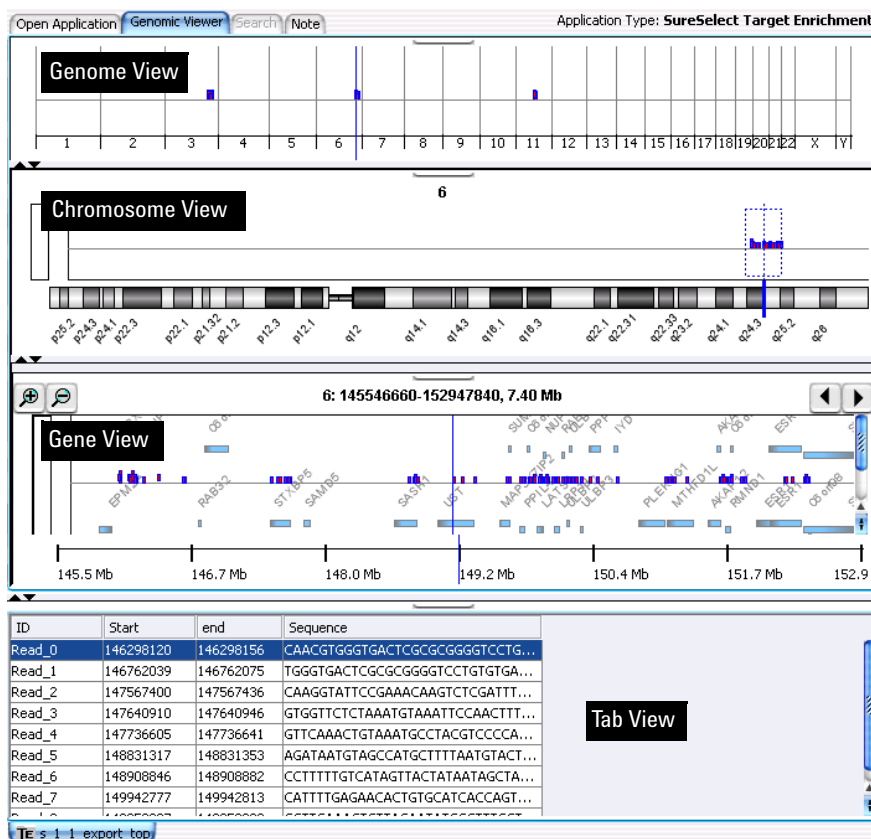


Figure 16 Genomic Viewer – SureSelect Target Enrichment application type

By default, for the SureSelect Target Enrichment application type, the Genomic Viewer appears in horizontal orientation. However, if desired, you can change the orientation to vertical. See [“To customize the appearance of an experiment in the Genomic Viewer”](#) on page 51.

The Genomic Viewer contains four Views:

Pane	Description
Genome View	Shows reads in the context of the genome of the species of interest. See “Genomic Viewer – Genome View” on page 93.
Chromosome View	Shows the reads that are mapped to a selected chromosome. See “Genomic Viewer – Chromosome View” on page 95.
Gene View	Shows an expanded view of a selected chromosomal region. The reads, target intervals, and genes in the selected region appear. See “Genomic Viewer – Gene View” on page 96.
Tab View	Shows a list of the reads on the selected chromosome. See “Genomic Viewer – Tab View” on page 98.

The four views are linked. When you select a location in one view, the other views change to display that location. The selected location appears as a blue line (“the cursor”) in Genome, Chromosome, and Gene Views.

Buttons available in all panes



Detach button. Detaches the view and opens it in a separate window. To re-attach the detached view to the main window, click its close box



Resize buttons. Increases or decreases the size of the given pane. You can also drag the border between panes to resize them.

Genomic Viewer – Genome View

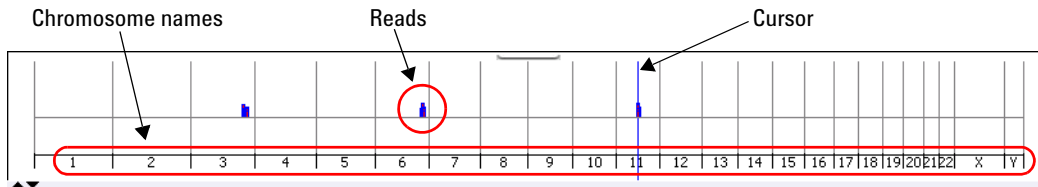


Figure 17 Genome View, with reads on chromosomes 3, 6, and 11. Chromosome 11 is selected.

When an experiment is selected, this view shows reads as blue and red bars in the context of the genome of the organism of interest. The chromosome that you select in this view sets the chromosome for the other views. Also, before you can view the results of enrichment analysis, you must first select the desired chromosome in Genome View. The program lets you view these results for one chromosome at a time. See [“To view and save reports”](#) on page 56.

- Chromosome Names** The names of the chromosomes of the organism, in order. The width of the space that each chromosome occupies is proportional to its size.
- Reads** The reads from the read file appear as blue and red bars. Their placement within the boundary of a chromosome reflects their approximate location within that chromosome.
- Cursor** The cursor appears as a thin blue line at a specific genomic location. To move the cursor, click the desired new location within Genome View. The other views adjust to reflect the new position of the cursor.

Shortcut menu

- Right-click anywhere within Genome View to open a menu with these options:

Option	Description
User Preferences	Opens the User Preference dialog box, where you can customize the appearance of genes and tracks in Gene View and enter the default data storage location. See “User Preferences – Tracks tab” on page 144 and “User Preferences – Miscellaneous tab” on page 142. You can also enter your login credentials for the eArray Web site.
View Preferences	Opens the View Preferences dialog box, where you can customize the orientation of the main program window. See “To customize the appearance of an experiment in the Genomic Viewer” on page 51 and “View Preferences” on page 147.
Select All	Marks the check boxes next to the names of all chromosomes. All chromosomes of the organism appear in Genome View.
Deselect All	Clears the check boxes next to the names of all chromosomes, except for the one in which the cursor currently appears.
Chromosomes	The names of all of the chromosomes of the current species appear. If the check box next to the name of a chromosome is marked, that chromosome appears in Genome View.

Genomic Viewer – Chromosome View

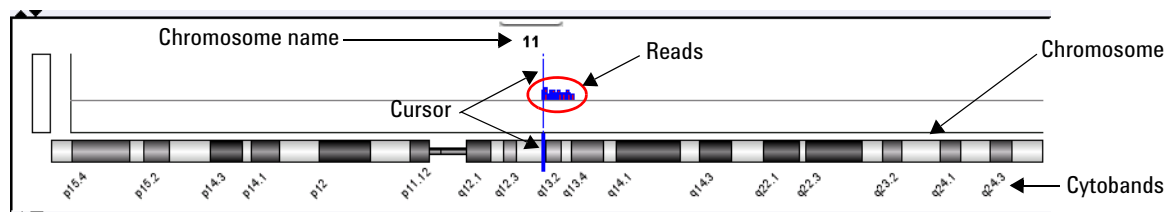


Figure 18 Chromosome View

This view shows an expanded view of the chromosome that is selected in Genome View. When an experiment is selected, reads appear as blue and red bars above the chromosomal locations to which they map. A region that you select in this view becomes the region that you can view in greater detail in Gene View.

Actions and shortcut menus

- Click a location within Chromosome View to move the cursor to the new location. The cursor in all views adjusts to the new location.
- Drag the pointer through the part of Chromosome View where reads appear, to select a region to view in greater detail. The selected region appears in Gene View.
- Right-click anywhere within Chromosome View to open a shortcut menu with these options:

Option	Description
User Preferences	Opens the User Preferences dialog box, where you can customize the appearance of genes and tracks in Gene View and enter the default data storage location. See “User Preferences – Tracks tab” on page 144 and “User Preferences – Miscellaneous tab” on page 142. You can also enter your login credentials for the eArray Web site.
View Preferences	Opens the View Preferences dialog box, where you can customize the orientation of the main program window. See “To customize the appearance of an experiment in the Genomic Viewer” on page 51 and “View Preferences” on page 147.

Genomic Viewer – Gene View

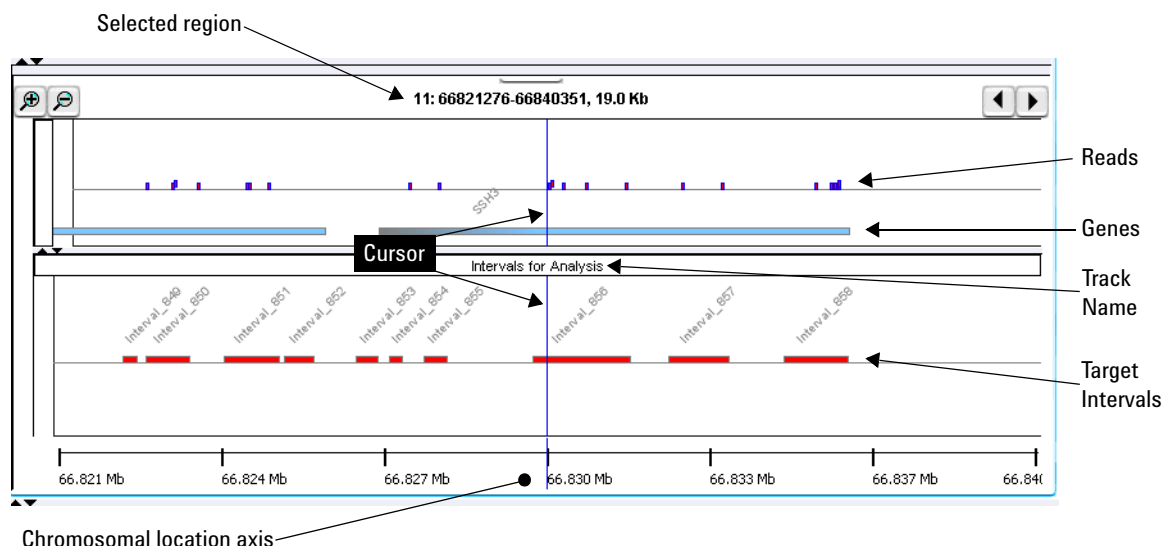


Figure 19 Gene View

Gene View displays the reads, target intervals, and genes that are associated with a specific region of a chromosome. Many zoom levels are available. See [“To view the data in an experiment”](#) on page 47.

Selected region The genomic coordinates of the chromosomal region that appears in Gene View.

Example: In [Figure 19](#), the notation **11:66821276-66840351, 19.0 Kb** tells you that Gene View currently shows chromosome 11, base pairs 66,821,276 to 66,840,351, which is a region that is approximately 19.0 Kb in size.



Zooms in to see a smaller region in greater detail.



Zooms out to see a larger region.



Scrolls left to lower-numbered base pairs on the chromosome.



Scrolls right to higher-numbered base pairs on the chromosome.

Reads	(Appear if an experiment is selected) The reads in the read file from the selected experiment appear as blue and red bars that are aligned with the genomic locations to which they map.
Genes	The genes that map to the currently displayed chromosomal region. A light blue bar indicates the genomic region that is covered by each gene. By default, the name of each gene also appears.
Cursor	The cursor is a thin blue line that indicates the center of the currently displayed chromosomal region. To change the location of the cursor, click one of the scroll buttons, or click anywhere within the gene display area of Gene View.
Track name	(Appears if an experiment is selected) The name of the target intervals track from the currently selected experiment. To see the target intervals track, you must expand Gene View.
Target intervals	(Appear if an experiment is selected) The target intervals from the selected experiment. By default, each target interval appears as a red bar that extends over the applicable genomic region. The name of each interval also appears.
Chromosomal location axis	This axis applies to the reads, genes, and tracks that appear in Gene View. The values on the axis change as you zoom in or out, or go to different chromosomal locations.

Actions and shortcut menus

- Click anywhere within the gene display area of Gene View to move the cursor to a new location.
- Right-click anywhere within the gene display area of the Gene View to open a shortcut menu with these options:

Option	Description
Create Gene List	This option does not apply to SureSelect Quality Analyzer.
Show in UCSC	Opens the View Coordinates in UCSC Browser dialog box, where you can create a user track based on the region that appears in Gene View. From this dialog box, the program can open the UCSC Genome Browser in your Web browser to view the region(s) that you defined. See the <i>Data Viewing User Guide</i> .

Option	Description
Search probes in eArray	Lets you search the eArray Web site for probes that are designed to regions that appear in Gene View, or that you define. For more information see the <i>Data Viewing User Guide</i> .
User Preferences	Opens the User Preference dialog box, where you can customize the appearance of genes and tracks in Gene View and set the default data storage location. See “ User Preferences – Tracks tab ” on page 144 and “ User Preferences – Miscellaneous tab ” on page 142. You can also set your login credentials for the eArray Web site.
View Preferences	Opens the View Preferences dialog box, where you can customize the orientation of the main program window. See “ To customize the appearance of an experiment in the Genomic Viewer ” on page 51 and “ View Preferences ” on page 147.

Genomic Viewer – Tab View



ID	Start	end	Sequence
Read_0	146298120	146298156	CAACGTGGGTGACTCGCGGGGTCCTG...
Read_1	146762039	146762075	TGGGTGACTCGCGGGGTCCTGTGTGA...
Read_2	147567400	147567436	CAAGGTATCCGAAACAAGTCTCGATT...
Read_3	147640910	147640946	GTGGTTCTCTAAATGTAAATCCAACTT...
Read_4	147736605	147736641	GTTCAAACGTAAATGCCTACGTCCCA...
Read_5	148831317	148831353	AGATAATGTAGCCATGCTTTAATGTACT...
Read_6	148908846	148908882	CCTTTTGTATAGTTACTATAATAGCTA...
Read_7	149942777	149942813	CATTTTGAGAACTGTGCATCACCAGT...

TE s_1-1-export_top

Figure 20 Tab View

When an experiment is selected, Tab View shows a list of the reads that are found on the currently selected chromosome.

- ID** The name of each read, as defined in the read file.
- Start** The first base pair on the chromosome to which each read maps.
- End** The last base pair on the chromosome to which each read maps.
- Sequence** The base sequence of each read.

Selected read To select a read, and move the cursor to its location, click anywhere in the row of the desired read. If you select a read, the program highlights it in blue.

Actions and shortcut menus

- Click anywhere within the row of a specific read to select it. When you select a read, the cursor location changes to the location of that read, and the other views adjust accordingly.
- Right-click anywhere in the column heading row to open a shortcut menu with a Scroll to Column option. This option opens the Scroll to Column dialog box, which lets you select a column, and scrolls Tab View so that the selected column is visible.
- In the column heading row, drag the right border of a column to increase or decrease the width of the column. This can be especially useful when you want to inspect the full base sequence of reads.
- Right-click any data item in Tab View to open a shortcut menu with these options:

Option	Description
Find in Column	Opens the Find in Column dialog box, where you can start a search for a text string in the column. See the <i>Data Viewing User Guide</i> .
Search Links	Many links to databases such as LocusLink, PubMed, and the UCSC Genome Browser appear. Each link passes the data item that you right-clicked to the selected database as a search string.
Customize Link	Opens the Customize Search Link dialog box, where you can create a new search link to include in the shortcut menu. See the <i>Data Viewing User Guide</i> .

Workflow Tab Window

The Workflow tab window presents a completely different set of panes than the Home, Quality Analyzer, View, Tool and Help tabs. This section presents a description of the panes for the Workflow Navigator, Parameter Panels and Summary Console/Progress tabs, as well as for the Workflow command ribbon.

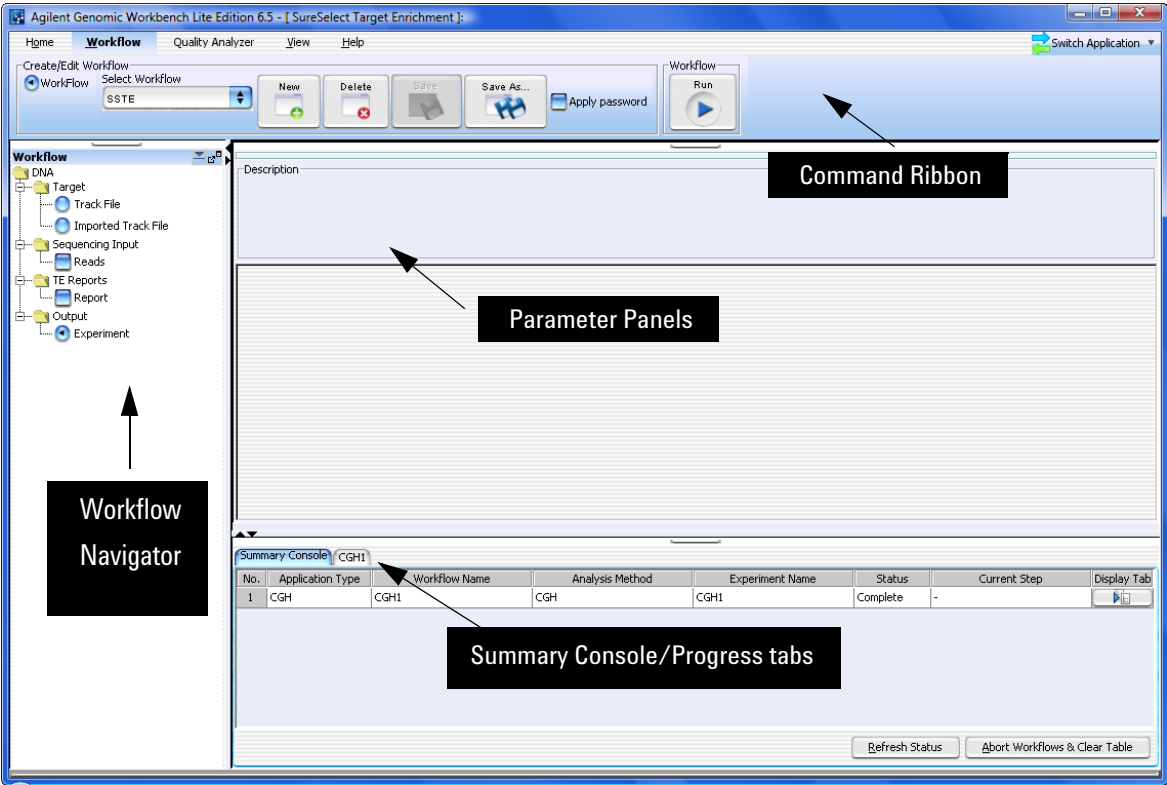


Figure 21 Workflow tab window

Figure 21 shows the tab panes when the Workflow tab is selected. This window contains the elements listed in Table 6 below. More detailed descriptions of each element appear later in this section.

Table 6 Workflow tab window elements

Element	Purpose
Command ribbon	Contains the commands for setting up, editing, deleting, saving, and running workflows. See “Command ribbon – Workflow tab” on page 102.
Tab pane – Workflow Navigator	Displays current workflow and options. As you select each option, the relevant parameters appear in the parameter panel. See “Workflow Navigator” on page 103.
Tab pane – Parameter Panels	Displays parameters for the selection in the Workflow Navigator. See “Parameter Panels – Overview” on page 105.
Tab pane – Summary Console/ Progress tabs	For Workflow, this view contains the following tabs: <ul style="list-style-type: none"> • Summary Console tab that displays the application type, workflow name, experiment name, status, and current step for each workflow completed or in process. See “Summary Console tab” on page 116. • Workflow Progress tab for each named workflow, that displays the status of the workflow progress, including any errors that might have occurred. See “Named workflow tabs” on page 118.

NOTE

In Workflow, the program only lets you set up and run workflow procedural steps to produce reports, and not explore the data. You must switch to the Quality Analyzer tab to display data and perform further analysis.

Command ribbon – Workflow tab

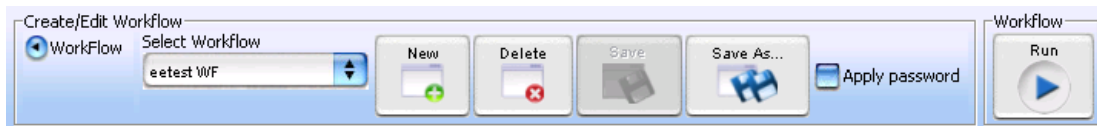


Figure 22 Command ribbon – Workflow tab

Purpose: Lets you set up, edit, delete, save, and run workflows.

To open: Set the application type to **SureSelect Target Enrichment**, then click **Workflow**.

Create/Edit Workflow

Select Workflow Shows the existing workflows in the program. To use an existing workflow, select one from this list.

New Opens the Create Workflow dialog box. To create a new workflow, type a name for the workflow in Enter Workflow Name, then click **OK**.

Delete (Available only if a workflow is selected.) Opens a Confirm dialog box that asks if you want to delete the workflow. To delete the current workflow from the program, click **Yes**.

Save (Available only if there are unsaved changes in the selected workflow.) Saves the changes in the selected workflow. (Note: This does not save the workflow in a file.)

Save As (Available only if workflow is selected.) Opens a Save As dialog box. To save the selected workflow with a new name, type the new name, then click **OK**. You must type a name that does not match an existing workflow or experiment.

Apply password Lets you type your password for a password-protected workflow.

Workflow

Run Starts the workflow. See “[Workflow Run button](#)” on page 114.

Workflow Navigator

The Workflow Navigator lets you set up QC reports and enrichment analyses so you can run them automatically.

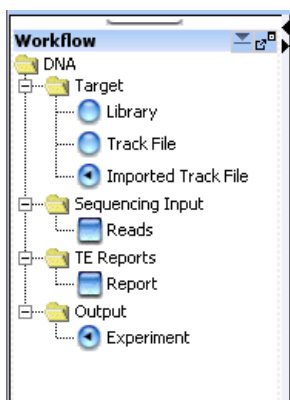


Figure 23 Workflow Navigator for SureSelect Quality Analyzer

To select any workflow option, click the option button or mark the check box next to its name. To display the parameters for a workflow option without changing its selection status, click the name of the option.

Target In Target, you select the target interval file for the workflow. To analyze the sequence data in a read file, the program must correlate reads with target genomic intervals. To define the target intervals, you import them as a BED format track file. You can also create the file manually. See [“To import target genomic regions as a track”](#) on page 41. Select one of these options:

Option	Description
Track File	Configures the workflow to import a track file in BED format directly. When you select this option, the Import Track Parameter Panel appears. Use this panel to select a track file for import. See “Import Track Parameter Panel” on page 107.
Select Imported Track	Configures the workflow to a track file that has already been imported with the Quality Analyzer. When you select this option, the Select Imported Data Parameter Panel appears. Use this panel to select a previously imported track for the workflow. See “Select Imported Track Parameter Panel” on page 108.

Sequencing Input

When you mark the Reads check box, the Sequencing Input Parameter panel appears. Use this panel to select and name a read file, as well as select a chromosome mapping file. See [“Sequencing Input Parameter Panel”](#) on page 109.

TE Reports

When you mark the Report check box the Text Summary Report Parameter panel appears. Use this panel to select the reports to be generated and the location to which they will be saved. See [“Text Summary Report Parameter Panel”](#) on page 112.

NOTE

You can select any two reports, but no more than two for any workflow.

Output

When you double-click Experiment, the Experiment Parameter Panel appears. Use this panel to name and describe the experiment that is created when you run the workflow. See [“Experiment Parameter Panel”](#) on page 106.

Parameter Panels – Overview

As you select options for a workflow in the Workflow Navigator, the parameters available for each option appear in specific parameter panels.

This section describes the parameter panels that appear in Workflow mode, organized alphabetically by name.

Below is a list of Parameter Panels for Workflow.

Table 7 Parameter Panels (in alphabetical order)

Option	Description
Experiment	Use this panel to name and describe the experiment that is created when you run the workflow. See “Experiment Parameter Panel” on page 106.
Import Track	Use this panel to select a track file for import. See “Import Track Parameter Panel” on page 107.
Select Imported Track	Use this panel to select a previously imported track for the workflow. See “Select Imported Track Parameter Panel” on page 108.
Sequencing Input	Use this panel to select and name a read file, as well as select a chromosome mapping file. See “Sequencing Input Parameter Panel” on page 109.
Text Summary Report	Use this panel to select the reports to be generated and the location to which they will be saved. See “Text Summary Report Parameter Panel” on page 112.

Experiment Parameter Panel

The screenshot shows a software window titled "Experiment Parameter Panel". It has a light blue header bar with the title. Below the header is a "Description" section with a text area containing the text: "When workflow run is complete, an experiment with specified name will be created and can be further analyzed from interactive mode." Below this is a large white area with a light gray grid background. In this area, there are two labels: "Experiment Name" and "Experiment Description". To the right of "Experiment Name" is a text input field containing "etest WF". To the right of "Experiment Description" is a larger text area containing "Workflow generated experiment".

Figure 24 Experiment Parameter Panel

Purpose: This parameter panel lets you customize the name of the output experiment for the workflow, and edit the description that the program saves with the experiment.

To open: The Experiment Parameter Panel appears when you click **Experiment** in Output in the Workflow Navigator.

Experiment Name The default name is the name of the workflow, but you can change it. The program creates a new experiment when it runs a workflow, and gives it the name that you enter here. After you run the workflow, the program makes the experiment available in the Quality Analyzer tab.

Experiment Description Type or edit an optional description for the experiment.

Import Track Parameter Panel

Import Track Parameter Panel

Description
Imports the track file which will be used as target.

Species: H. sapiens
Build Name: hg18
Track Name:
Track File:

Color: ■

Figure 25 Workflow – Import Track Parameter Panel

Purpose: This panel lets you import and change the name of a track file that provides the target intervals against which the sequencing reads will be compared. See [“To import target genomic regions as a track”](#) on page 41.

To open: This parameter panel is shown when you click **Track File** under Target in the Workflow Navigator.

- Species** Select the species that is associated with the genomic intervals in your BED format track file.
- Build Name** Select the genome build that is associated with the genomic intervals in your BED format track file.
- Color** (Optional) Customizes the display color of the intervals in the target intervals track. Click **Change** to switch colors. See [“Choose Track Color”](#) on page 120.

- Track Name** Type a name for the track as you want it to appear in reports and lists, and in Gene View of the Genomic Viewer.
- Track File** The track file must be a BED format track file that contains the target genomic intervals. Click **Browse** to select the desired *.bed track file.

Select Imported Track Parameter Panel

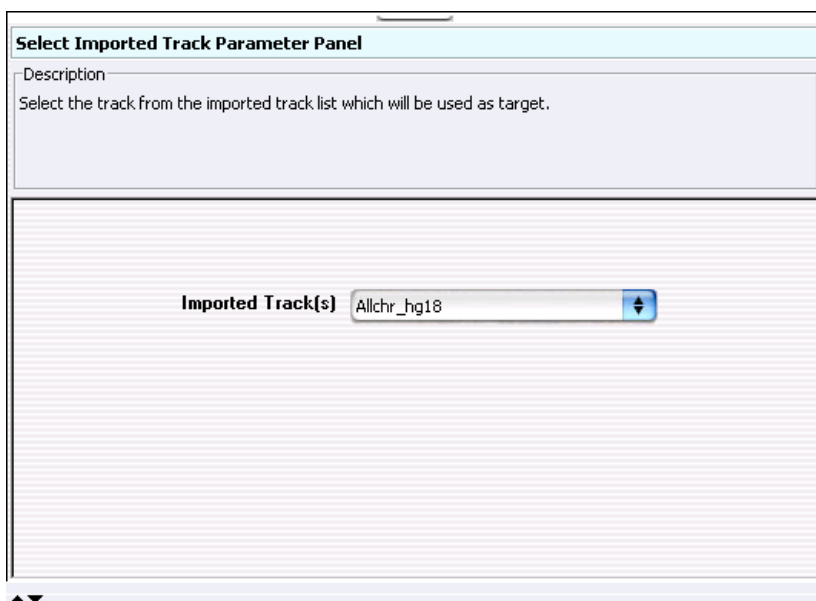


Figure 26 Workflow – Select Imported Track Parameter Panel

Purpose: This parameter panel lets you select previously imported tracks as Target input for the workflow.

To open: The Select Imported Track Parameter Panel appears when you click **Select Imported Track** under Target in the Workflow Navigator.

Imported Track(s) Uses an already imported track for the target intervals. Select the track from the list.

Sequencing Input Parameter Panel

Sequencing Input Parameter Panel

Description
Imports the sequence file.

Read Info.

Species: H. sapiens File Type: Illumina GA Export

Build Name: hg18 Read length: 36

File Type: ☒ Single Type ☐ Paired Type

Seq File 1 Seq File 2

☒ Select already imported Read File ☐ Import new Read File

Read Name:

Chr Mapping File: Browse...

*Provide mapping (tab delimited) between chromosome names in Read file to genome build chromosome names

Sequence File: Browse...

Figure 27 Workflow – Sequencing Input Parameter Panel

Purpose: This panel is used to select sequencing read files and their accompanying information and chromosome mapping files.

To open: This panel is displayed when you mark **Reads** under Sequencing Input in the Workflow Navigator.

Read Info

Species Select the species that is represented in the imported sequence read file.

Build Name The available genome builds for the selected species appear in the list. Select the build that applies to the reads in the imported file.

File Type Select the specific data format that applies to the imported file. The program can import *.zip archives that contain these types of files:

File Type	Comments
Illumina GA Export	Revised version of the ELAND file format that contains additional fields, including quality scores for all bases. You cannot import a multiplexed read file into Agilent Genomic Workbench. You must first separate the file into individual indexed sample files, then import these.
AB SOLiD ma	Color space *.ma file produced by the Applied Biosystems SOLiD system. The AB SOLiD alignment software produces this CSFASTA format file, and places sequence alignment information in the FASTA headers. Color calls are converted to bases when the file is imported.

NOTE

You can import sequence read files in the gzipped or *.zip formats, but not in the .zipx format. In addition, Agilent Genomic Workbench cannot handle a compressed file whose uncompressed file is greater than 4G. In this case, import the uncompressed file directly.

Read Length Average length of the DNA fragments in the sequencing read file

File Type

Single Type Click if you have only one file of reads.

Paired Type If you have two output files in which the reads are paired, click this option, then select the paired file. The program imports the two files separately, then merges them before analysis.

Seq File N (where N = 1, 2, etc.):

Select already imported Read File

Selected Reads Select a Read File from the list of already imported ones.

Import a new Read File

Read Name You can change the name of the selected sequence read file.

Chr Mapping File The name and location of the applicable chromosome mapping file. Because the names of chromosomes in the imported read file can differ from the names of chromosomes in the selected genome build, you must import a chromosome mapping file that relates the chromosome names in these two locations.

Browse – Opens an Open dialog box, where you can select the desired chromosome mapping file.

File format information: Create a tab-delimited text file (*.txt) that contains the desired mapping information. Enter one chromosome per line. Give the name of the chromosome in the read file followed by the name of the matching chromosome in the genome build, separated by a tab character. Do not include a header row.

Example file content:

```
chrA.fa      chrA
chrB.fa      chrB
chrC.fa      chrC
```

Sequence File **Browse** – Opens an **Open** dialog box, where you can select the desired sequence read file.

Text Summary Report Parameter Panel

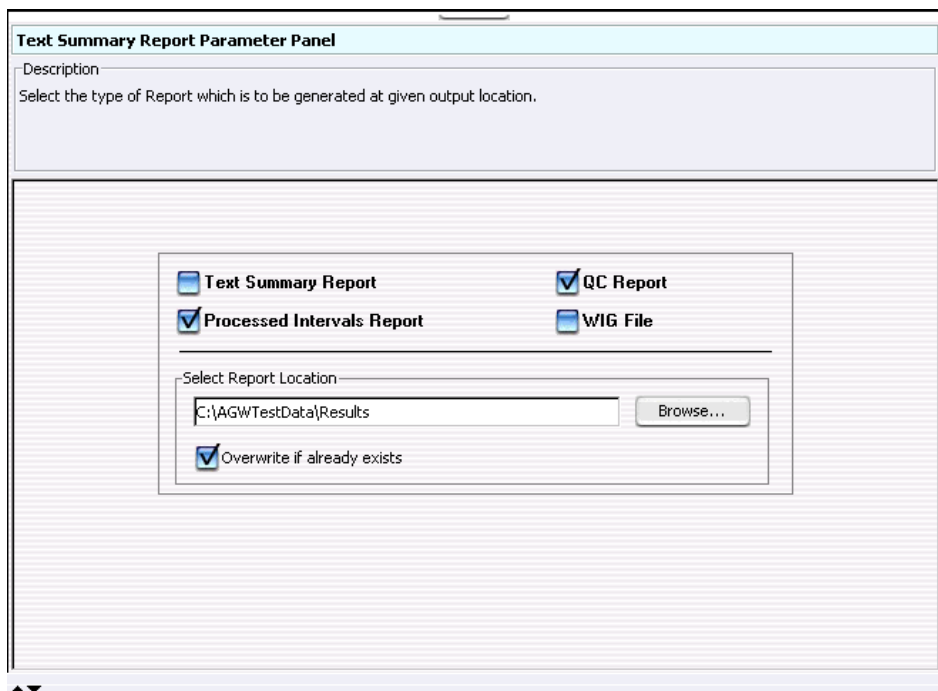


Figure 28 Workflow – Text Summary Report Parameter Panel

Purpose: This parameter panel lets you select the type of report you want to generate, and a location for it.

To open: The Text Summary Report Parameter Panel appears when you mark **Report** under TE Reports in the Workflow Navigator.

NOTE

You can select any two reports, but no more than two for any workflow.

These reports are also described in the Target Enrichment Reports section in this chapter. See [“Target Enrichment Reports”](#) on page 149.

Text Summary Report	For this report, the program creates a *.txt file that contains a table of QC metrics and the results of read depth analysis. You can open this file with a text editor or word processor.
QC Report	For this report, the program creates a PDF format report that contains tables of QC metrics. It also contains the results of read depth analysis shown both as tables and as a graph.
Processed Intervals Report	For this report, the program creates a *.xls file that contains enrichment analysis results for all the chromosomes. You can view this file in Microsoft Excel.
WIG File	The program creates a track file that contains enrichment analysis results for all the chromosomes. You must select one chromosome and then upload the file to the UCSC Genome Browser to see results.
Select Report Location	Click Browse to select a folder to contain the workflow-generated reports.
Overwrite if already existing	Clear if you intend to have new report files overwrite existing ones. (Marked check box is the default.)

Workflow Run button



Figure 29 Workflow Run button

Purpose: This button is used to initiate the current workflow.

To open: This button is located on the Workflow command ribbon.

When you click this button, you are prompted to type a temporary identifier for the workflow. See [“Provide Workflow Identifier”](#) on page 133 and [“To run a workflow”](#) on page 69.

The workflow run then starts, using the Target, Sequencing Input and TE Reports parameters you have set.

A new row is created in the Summary Console for this workflow. See [“Summary Console tab”](#) on page 116.

A new tab is created in the Summary Console/Progress pane. See [“Named workflow tabs”](#) on page 118.

An experiment is created for the workflow in the Experiment pane for interactive review of the results. See [“Navigator – Experiment pane”](#) on page 88.

To run a series of workflows, see [“To run a series of workflows”](#) on page 70.

NOTE

On 64-bit computers, you can run up to three workflows simultaneously. On 32-bit computers, workflows are processed sequentially, not simultaneously.

Summary Console/Progress tabs – Overview

In Workflow mode, workflow progress tabs provide information on the status of each workflow you run. The Summary Console tab displays the workflows you run, and the status of each. Named workflow tabs show information about each single workflow run and a running log of progress through the workflow.


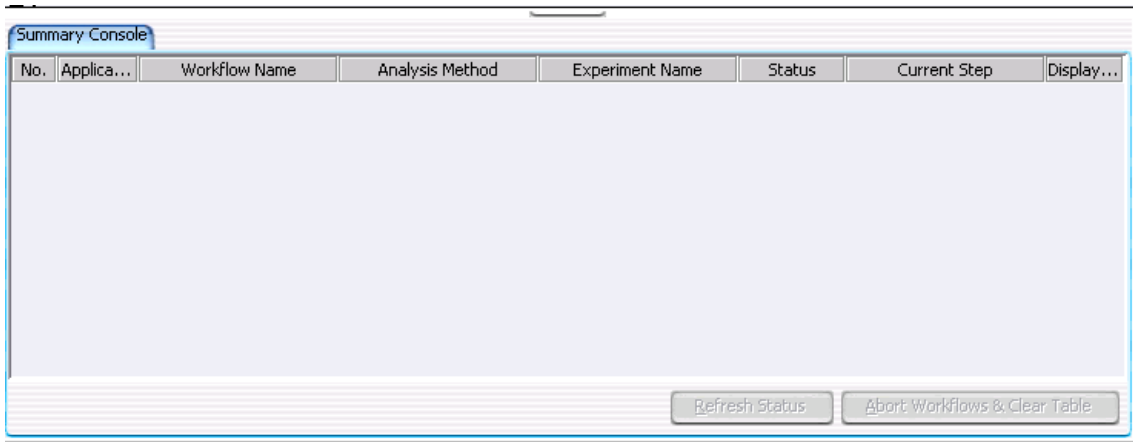


Figure 30 Summary Console before run

For information on the Summary Console and Named workflow tabs during a run or runs, see “[Summary Console tab](#)” on page 116 and “[Named workflow tabs](#)” on page 118.

Summary Console tab

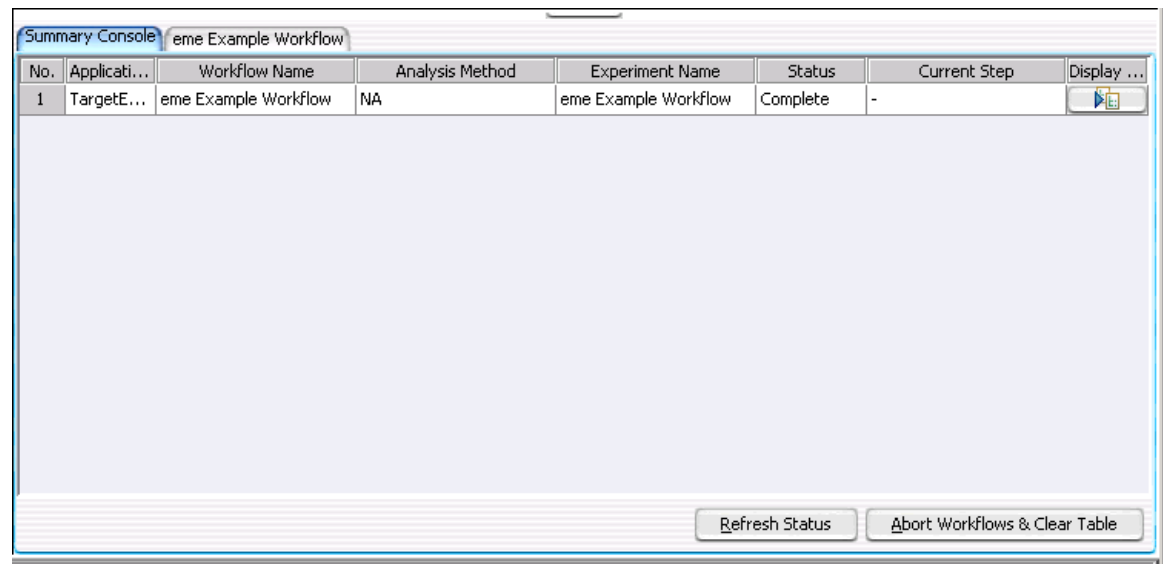


No.	Applicati...	Workflow Name	Analysis Method	Experiment Name	Status	Current Step	Display ...
1	TargetE...	eme Example Workflow	NA	eme Example Workflow	Complete	-	




Figure 31 Summary Console tab after run

The Summary Console tab displays the workflows you run, and gives basic identifying information and the workflow status. It also lets you manage the workflows.

NOTE

The Summary Console and progress tabs include workflows for CGH, ChIP and SureSelect Target Enrichment. For information on the CGH and ChIP workflows, see the *Workflow User Guide*.

No.	(Read-only) The order in which the workflows were started, from first to most recent.
Application Type	(Read-only) The basic application type (CGH or ChIP, for example) for the workflow.
Workflow Name	(Read Only) The name of the workflow you are using for this run.

Analysis Method	(Not applicable to Sure Select Quality Analyzer workflows. Only applicable to CGH or ChIP workflows.) If you are working in the SureSelect Quality Analyzer Workflow tab window and if you have started a CGH or ChIP workflow, you will see the information for the CGH or ChIP workflow in this console.
Experiment Name	(Read-only) The name of the output experiment that the workflow generates. You look at the experiment, including its results, in the interactive tabs.
Status	(Read-only) Indicates whether the workflow is Running or Complete .
Current Step	(Read-only) For a workflow that is running, shows the progress through the steps of the workflow.
Display Tab	In this column, in the row of the desired workflow, click  to open the tab for the selected workflow. This lets you review information about the workflow.
Refresh Status	Updates the status of all workflows. The program also updates the status of the workflows in the workflow list in real time.
Abort Workflows & Clear Table	Opens a Confirm dialog box that asks if you are sure you want to abort all workflows. If you click Yes , the program stops the execution of all workflows, and removes all workflows from the list. It also removes all named workflow tabs.
	(Available if tabs are hidden past the left edge of the pane) Shifts the display of tabs to the left to reveal hidden tabs.
	(Available if tabs are hidden past the right edge of the pane) Shifts the display of tabs to the right to reveal hidden tabs.

Named workflow tabs

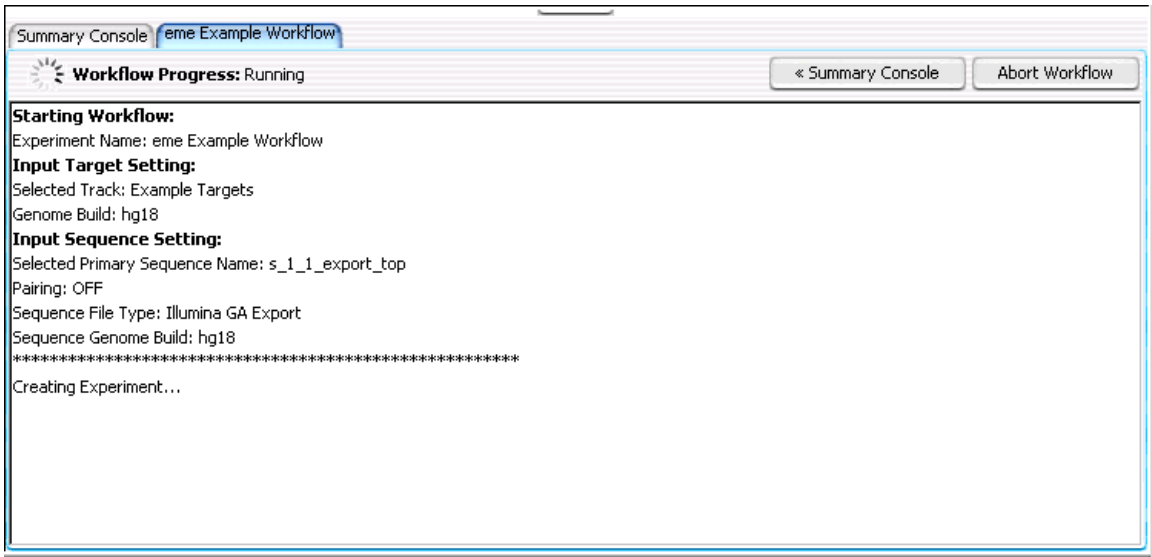


Figure 32 Named Workflow tab

The program creates a separate tab for each workflow that you run. The name displayed is the name you specified in the Workflow Identifier. Each workflow tab shows a log of workflow activity.

Workflow Progress	Shows the status of the workflow, either Running or Complete .
Summary Console	Opens the Summary Console tab, where you see a list of all workflows and check the status of each. See “Summary Console tab” on page 116.
Abort Workflow	Appears only when a workflow is running and that workflow’s progress tab is selected. Click to abort the running workflow and remove its tab from the Summary Console pane.
Close Tab	Removes the current tab, and removes its row in the Summary Console Tab.
Text Box	Displays a running log of workflow activity.



(Available if tabs are hidden past the left edge of the pane) Shifts the display of tabs to the left to reveal hidden tabs.



(Available if tabs are hidden past the right edge of the pane.) Shifts the display of tabs to the right to reveal hidden tabs.

Dialog Boxes

Choose Track Color

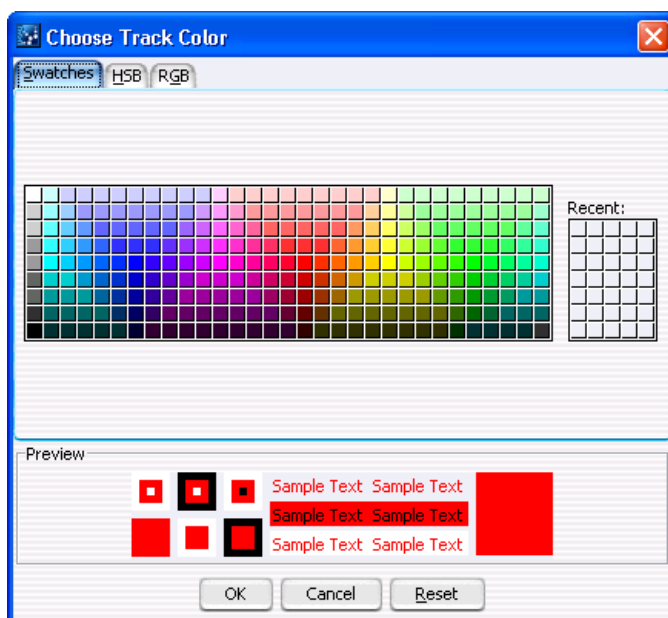


Figure 33 Choose Track Color dialog box – Swatches tab

Purpose: For the SureSelect Target Enrichment application type, lets you select a display color for the bars that represent genomic regions in the target intervals track in Gene View. See [“To import target genomic regions as a track”](#) on page 41.

To open: In the Import Track dialog box (see [“Import Track”](#) on page 131), under Color, click **Change**.

Swatches tab

- Swatches** To select a color, click a swatch. The new color appears in the Preview pane and also under Recent. The color that you select also becomes the selected color in the HSB and RGB tabs, where you can further refine the color.
- Recent** Displays the colors that you have recently selected in the Swatches tab. To select one of these recent colors, click it.

HSB tab

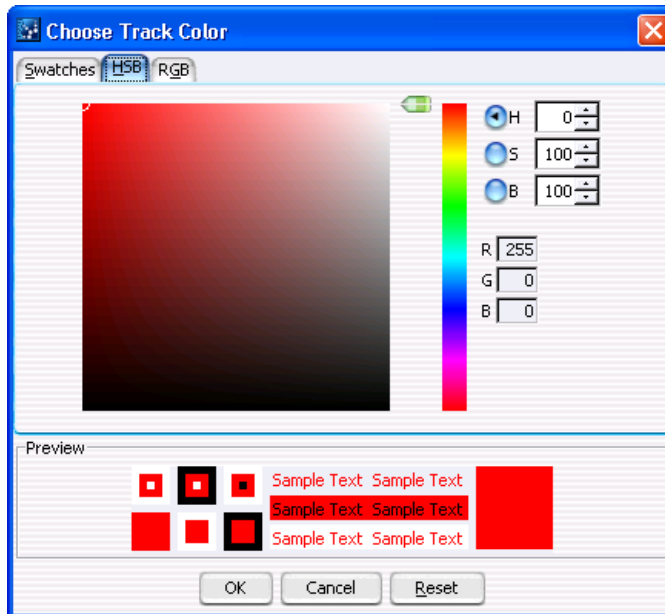


Figure 34 Choose Track Color dialog box – HSB tab

This tab lets you define the hue (H), saturation (S), and brightness (B) levels of for a color. These three values uniquely define a color. The initial color settings in this tab reflect any changes that you have made to the currently selected color in the other tabs. You can set these values in several ways:

- Directly edit the numbers in **H**, **S**, and **B**. You can also click the up or down button to the right of each value to increase or decrease it. In addition, to use the up and down arrow keys on your keyboard to change values, click the number that you want to change, then press the up or down arrow key, as desired.
- Use the green slider to change values. Select **H**, **S**, or **B**, as desired, then drag the slider up or down to change the selected value.
- Use the green slider and the large, square color selection box. Select **H**, **S**, or **B**, then drag the slider up or down to set the desired value. The color selection box shows all of the available colors given the particular setting of the slider. To select a color, click anywhere within the color selection box. This sets the remaining two HSB values.

H Hue – A number from 0 to 359 that represents the basic color. The color spectrum is a 360 degree color circle.

S Saturation – A number from 0 to 100 that represents the intensity of the color. A setting of 100 gives maximum color intensity. A setting of 0 gives no color, and reduces the available color spectrum to grayscale, only.

B Brightness – A number from 0 to 100 that represents the amount of black that is mixed in with the color. A setting of 100 gives maximum brightness, with no black added, and a setting of 0 results in the color black, without regard to the other settings.

R, G, B Each of these values shows the color settings of the selected color using the RGB (Red, Green, Blue) color model. You cannot directly edit the values from HSB tab, but they change when you select a different color. You can set specific RGB values in the RGB tab.

RGB tab

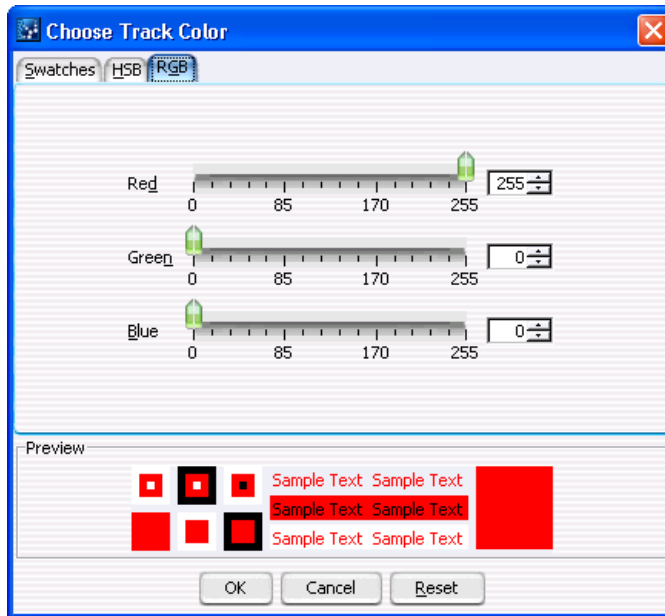


Figure 35 Choose Track Color dialog box – RGB tab

This tab lets you use the RGB (red-green-blue) color model to define a color. In this model, you select the amounts of red, green, and blue to combine to form the desired color. All colors can be defined.

You can change the R, G, or B value of the selected color in two ways:

- In **Red**, **Green**, or **Blue**, drag the green slider to the desired value.
- Directly edit the number in **Red**, **Green**, or **Blue**. You can also click the up and down buttons to the right of the value to increase or decrease it. In addition, to use the up and down arrow keys on your keyboard to change a value, click the number that you want to change, then press the up or down arrow key, as desired.

Red A number from 0 to 255 that represents the amount of red. 0 is the minimum and 255 is the maximum.


Green A number from 0 to 255 that represents the amount of green. 0 is the minimum and 255 is the maximum.

Blue A number from 0 to 255 that represents the amount of blue. 0 is the minimum and 255 is the maximum.

Items that appear in all tabs

- Preview** Shows the selected color in a number of contexts. The right-most diagram in this pane can show two colors:
- **Top color** – The color that is currently selected for the track.
 - **Bottom color** – The color that is currently selected to replace it.
- OK** Accepts any changes that you made to the color, and closes the dialog box.
- Cancel** Closes the dialog box, and discards any changes that you made to the color.
- Reset** Restores the color settings in the dialog box to what they were before you made any changes. The dialog box remains open.

Create Experiment

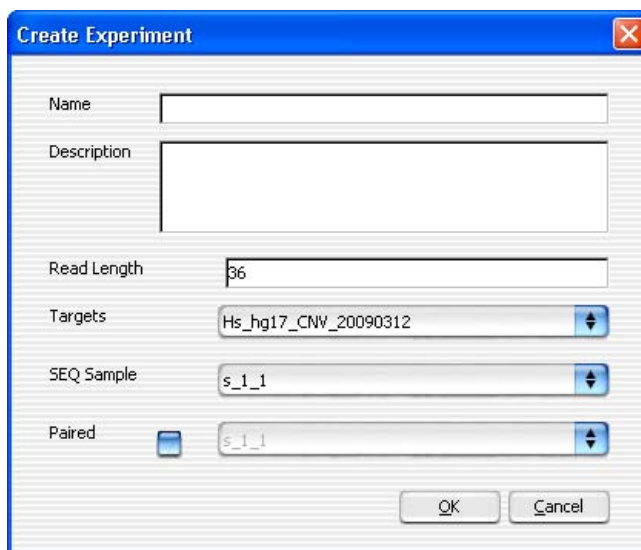

The 'Create Experiment' dialog box is shown with the following fields and values:

- Name: [Empty text box]
- Description: [Empty text box]
- Read Length: 36
- Targets: Hs_hg17_CNV_20090312
- SEQ Sample: s_1_1
- Paired: s_1_1

Buttons: OK, Cancel

Figure 36 Create Experiment dialog box

Purpose: Lets you set up a QC analysis experiment. You use this dialog box to name the experiment, and to select a sequence read file and a set of target intervals for the analysis. See “[To create a new experiment](#)” on page 44.

To open: In the Home tab, click **Create Experiment**. Alternatively, in the Experiment pane of the Navigator, right-click the **Experiments** folder, then click **New Experiment**.

Name	Type a name for the experiment. This name identifies the experiment in the Experiment pane of the Navigator.
Description	(Optional) Type a brief description of the experiment.
Read Length	Type the number of bases in each read in the selected read file. The program assumes all reads in the file have the same length.
Targets	Shows a list of all of the annotation tracks that are available in the program. Select the track that contains the desired target intervals for analysis.
SEQ Sample	Shows a list of the names of all imported sequence read files that are available for analysis. Select the desired sequence read file.
Paired	If you have two output files in which the reads are paired, mark this option, then select the paired file. The program imports the two files separately, then merges them before analysis.
OK	Creates a new experiment with the given name that contains links to the selected sequence read file and target interval track. Creates a new folder in the Experiment pane of the Navigator that bears the name of the experiment.
Cancel	Closes the dialog box, but does not create a new experiment.

NOTE

Before you create an experiment, you must import the desired sequence read file, target file, and chromosome mapping file into the program. See “[Importing Files](#)” on page 38.

Create Workflow

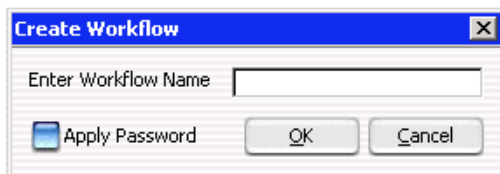


Figure 37 Create Workflow dialog box

Purpose: Lets you set up a new workflow. You use this dialog box to name the workflow, and to apply a password if you want to restrict access to this workflow. See [“To create a new workflow”](#) on page 60.

To open: In the Workflow command ribbon, click **New** or click **Save As**.

**Enter Workflow
Name**

Enter the name of the workflow you want to create. It can be the same as the Experiment name or different.

Apply Password

Mark the check box to restrict access to the workflow. The Set Password dialog box appears. See [“Set Password”](#) on page 134.

Export

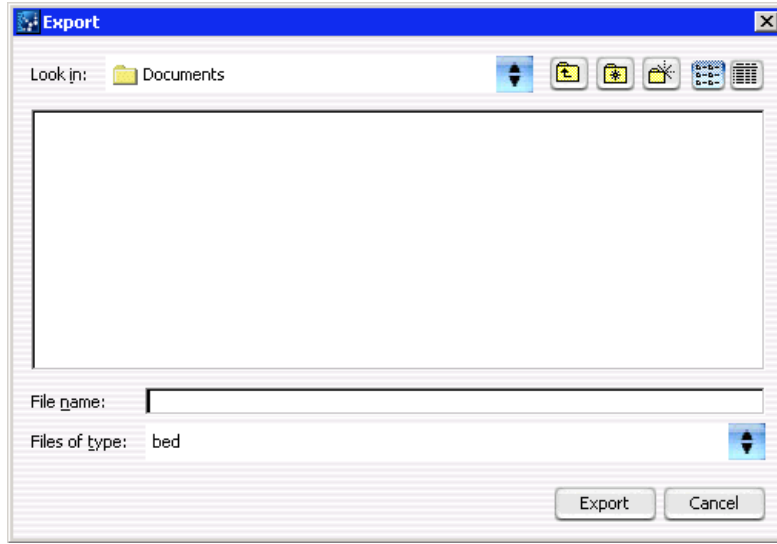


Figure 38 Export dialog box

Purpose: Used to designate a location and file name to export tracks.

To open: This dialog box opens when you click **OK** from the Export Tracks dialog box. See “[Export Tracks](#)” on page 128.

Use the buttons at the top of the dialog box to change the display.

- Look in** Click the arrow and select the folder where you wish to export the file.
- File name** Type the name you wish to use for the exported file.
- Files of type** Shows the type of files displayed in the window. Click the arrow to change the type of files displayed.
- Export** Click to export the currently-selected tracks to the file.
- Cancel** Click to cancel the operation.

Export Tracks

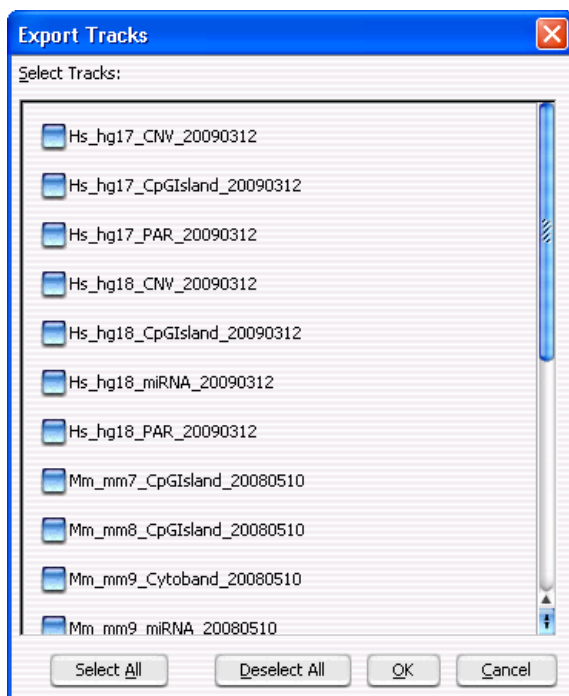


Figure 39 Export Tracks dialog box

Purpose: Lets you select one or more tracks for export as a single BED format track file. You can use this functionality to combine multiple tracks, and also to transfer a track to another computer. In addition, you can view the track in any compatible genome browser. See [“To export tracks”](#) on page 50.

To open: In the Home tab, click **Export > Tracks**.

Select Tracks Shows a list of all of the tracks that are available in the program. Mark the check box next to a track to select it for export.

Select All Marks all tracks for export.

Deselect All Clears the check boxes next to all tracks.

- OK** Opens an Export dialog box, where you can select a location for the exported BED format track file.
- Cancel** Closes the dialog box, but does not export any tracks.

Go to Gene/Genomic Location

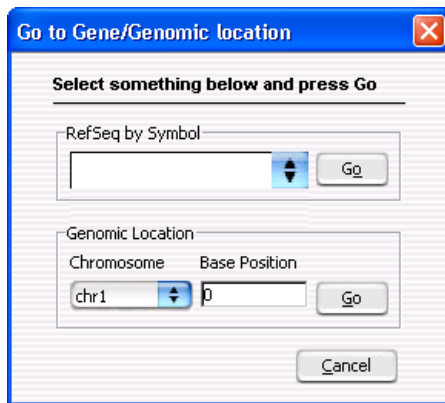


Figure 40 Go to Gene/Genomic Location dialog box

Purpose: Moves the cursor in the Genomic Viewer to a specific location. You can enter either a gene symbol or a genomic location. See [“To customize the appearance of an experiment in the Genomic Viewer”](#) on page 51.

To open: In the **Home** tab, click **Go to Gene/Genomic Location**.

- RefSeq by Symbol** Type a valid RefSeq gene symbol.
- Go** – Moves the cursor to the location of the gene that you entered.
- Genomic Location** **Chromosome** – Select the desired chromosome.
- Base Position** – Type the desired base position on the selected chromosome.
- Go** – Moves the cursor to the specified genomic location.
- Cancel** Closes the dialog box, but does not move the cursor.

Import Genome Build

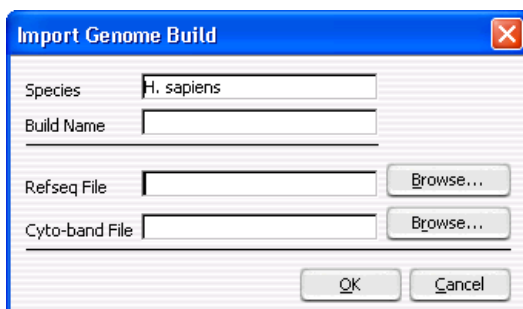


Figure 41 Import Genome Build dialog box

Purpose: Lets you import a new genome build into the program. See [“To import a genome build”](#) on page 43. Genome builds contain the names and locations of genes and cytobands. However, they do not contain any nucleotide sequence data.

To open: In the **Home** tab, click **Import > Genome Build**.

Species The species represented by the new genome build.

Build Name A name for the genome build, as you want it to appear in the program.

RefSeq file The location of the file that contains gene symbol information for the genome.

Browse – Opens an Open dialog box, where you can select the desired file.

Cytoband file The location of the file that contains cytoband information for the genome.

Browse – Opens an Open dialog box, where you can select the desired file.

OK Imports the genome build.

Cancel Closes the dialog box, and does not import a genome build.

CAUTION

Use only Agilent-supplied genome build files. For more information, see the online help on the eArray Web site.

Import Track

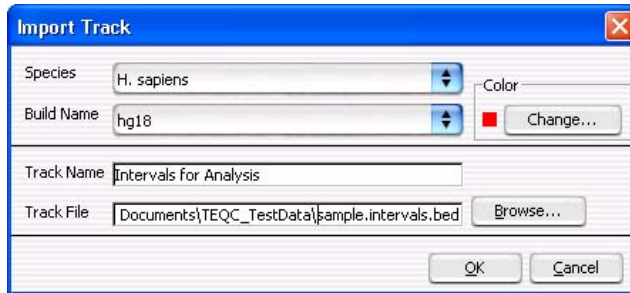


Figure 42 Import Track dialog box

Purpose: Lets you import a BED format track file into the program. For SureSelect Quality Analyzer, this is how you import target genomic intervals for analysis experiments. See [“To import target genomic regions as a track”](#) on page 41.

To open: In the **Quality Analyzer** tab, click **Import > Track**.

- Species** Shows a list of the species for which genome builds are available in the program. Select the desired species.
- Build Name** Shows a list of the genome build(s) that are available for the selected species. Select the desired genome build. If the desired genome build does not appear in the list, you must import it. See [“To import a genome build”](#) on page 43.
- Color** The color swatch shows the currently selected display color for the bars that represent the target intervals in Gene View.
Change – Opens the Choose Track Color dialog box, where you can select a new color for the track. See [“Choose Track Color”](#) on page 120.
- Track Name** The program uses this name to identify the track in Gene View, and also in lists. Type a name for the new track.
- Track File** The location of the BED format track file to be imported.
Browse – Opens a dialog box, where you can import the file.
- OK** Imports the track and closes the dialog box.

Cancel Closes the dialog box, and does not import a track.

Open

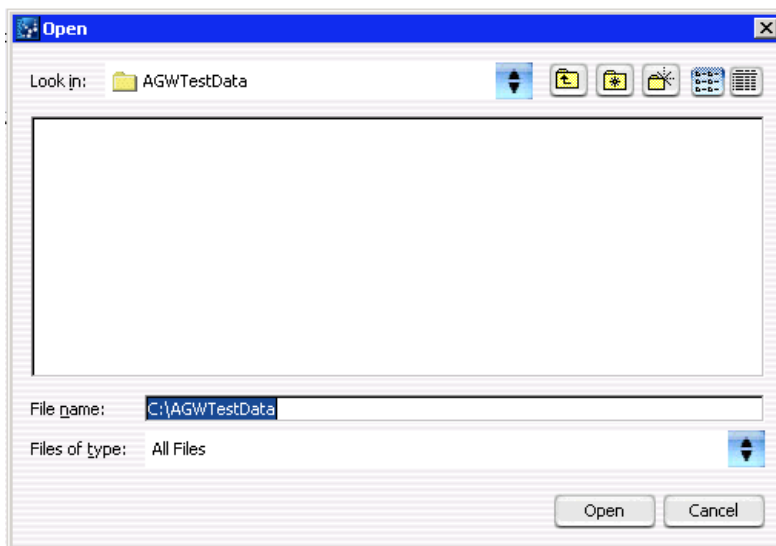


Figure 43 Open dialog box

Purpose: To select the files you want to open, or to identify the location where you want files to be stored.

To open: This dialog box appears when you click **Browse** in a parameter panel or dialog box.

Look in Select folder where the files are located. To browse for a location, click the arrow and browse to the desired folder.

File name Type the name for the file you want to open, or click the file to select it. To select multiple files to open, hold down the **Ctrl** key and click the files to open.

Files of type Expected file type is displayed. Click to change displayed file types.

Open Click **Open** to select the folder or file of interest.

Cancel Click this to cancel the operation.

Provide Workflow Identifier

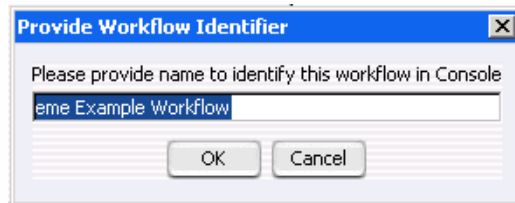


Figure 44 Provide Workflow Identifier dialog box

Purpose: To designate a name used in the named workflow tabs to identify a workflow to be run. See “[Named workflow tabs](#)” on page 118.

To open: This dialog box appears when you click **Run** to run a workflow. See “[Workflow Run button](#)” on page 114, “[To run a workflow](#)” on page 69 and “[To run a series of workflows](#)” on page 70.

**Please provide
name to identify
this workflow in
Console**

Type a name for the workflow progress tab.

OK Click to accept the workflow identifier and start the workflow.

Cancel Click to cancel the operation without running the workflow.

Set Password

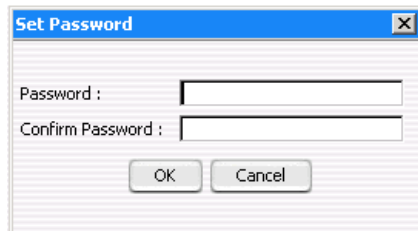


Figure 45 Set Password dialog box

Purpose: To restrict access to a workflow.

To open: Mark **Apply Password** in the Workflow command ribbon or click **New** or **Save As** to bring up the Create Workflow dialog box, mark **Apply Password**, then click **OK**.

Password Type a password.

Confirm Password Type the password again.

OK Click to set password.

Select SEQ Files

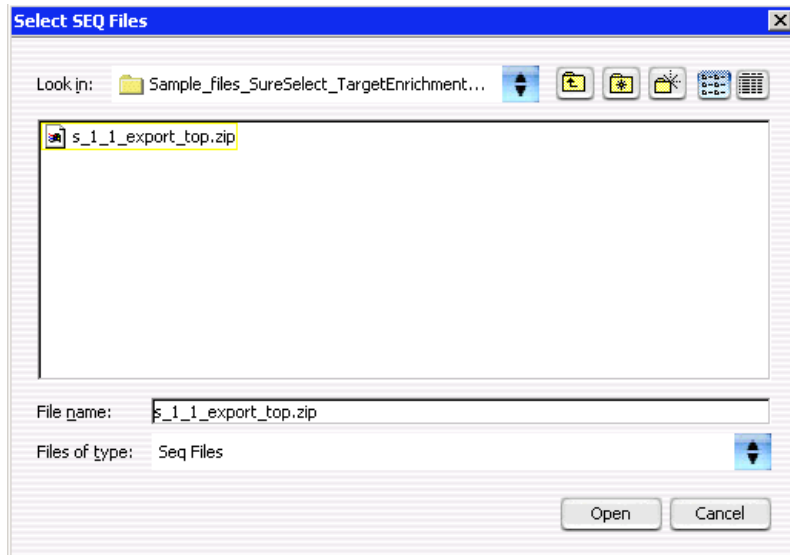


Figure 46 Select SEQ Files dialog box

Purpose: Lets you locate and import a sequence read zip file.

To open: In the Quality Analyzer command ribbon you click **Import** and select **Reads**. See [“Specify Input Information”](#) on page 136 to complete importing the read file.

Look in Select folder where the files are located. To browse for a location, click the arrow and browse to the desired folder.

File name Type the name for the file you want to open, or click the file to select it. To select multiple files to open, hold down the **Ctrl** key and click the files to open.

Files of type Expected file type is displayed. Click to change displayed file types.

Open Click **Open** to select the folder or file of interest.

Cancel Click this to cancel the operation.

Specify Input Information

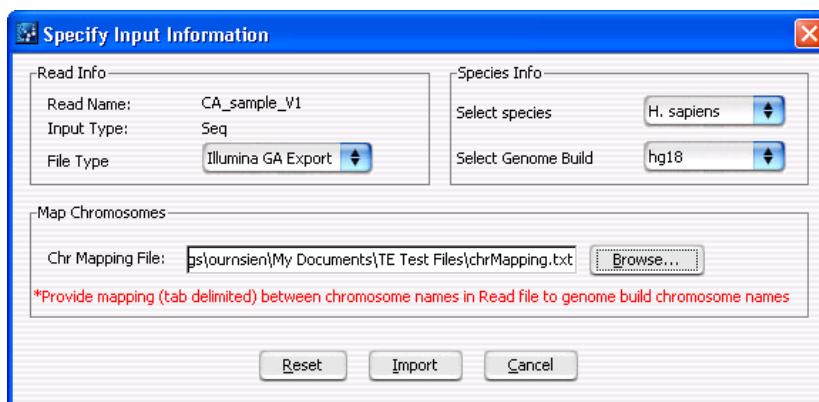


Figure 47 Specify Input Information dialog box

Purpose: Lets you define the attributes of a sequence read file as you import it. See [“To import sequence read files”](#) on page 38 and [“Select SEQ Files”](#) on page 135.

To open: In the **Quality Analyzer** tab, click **Import > Reads**. In the dialog box that appears, select a *.zip file that contains the desired sequence read file, then click **Open**.

Read Info

Read Name (Read-only) The name of the selected sequence read file.

Input Type (Read-only) **Seq** appears, which denotes a file that contains sequence reads.

File Type Select the specific data format that applies to the imported file. The program can import *.zip archives that contain these types of files:

File Type	Comments
Illumina GA Export	Revised version of the ELAND file format that contains additional fields, including quality scores for all bases.
AB SOLiD ma	Color space *.ma file produced by the Applied Biosystems SOLiD system. The AB SOLiD alignment software produces this CSFASTA format file, and places sequence alignment information in the FASTA headers. Color calls are converted to bases when the file is imported.

NOTE

Import sequence read files as *.zip archives. Do not use the gzip format.

Species Info

Select Species Select the species that is represented in the imported sequence read file.

Select Genome Build The available genome builds for the selected species appear in the list. Select the build that applies to the reads in the imported file.

Map Chromosomes

Because the names of chromosomes in the imported read file can differ from the names of chromosomes in the selected genome build, you must import a chromosome mapping file that relates the chromosome names in these two locations.

Chr Mapping File The name and location of the applicable chromosome mapping file.

Browse – Opens an Open dialog box, where you can select the desired chromosome mapping file.

File format information: Create a tab-delimited text file (*.txt) that contains the desired mapping information. Enter one chromosome per line. Give the name of the chromosome in the read file followed by the name of the matching chromosome in the genome build, separated by a tab character. Do not include a header row.

Example file content:

```
chrA.fa      chrA
chrB.fa      chrB
chrC.fa      chrC
```

Other commands

- Reset** Opens a dialog box that asks if you are sure you want to reset the mapping. If you click Yes, the program restores the settings in the Specify Input Information dialog box to what they were before you made any changes.
- Import** Imports the sequence read file into the program, and makes it available for analysis.
- Cancel** Closes the dialog box, but does not import the read file.

Track

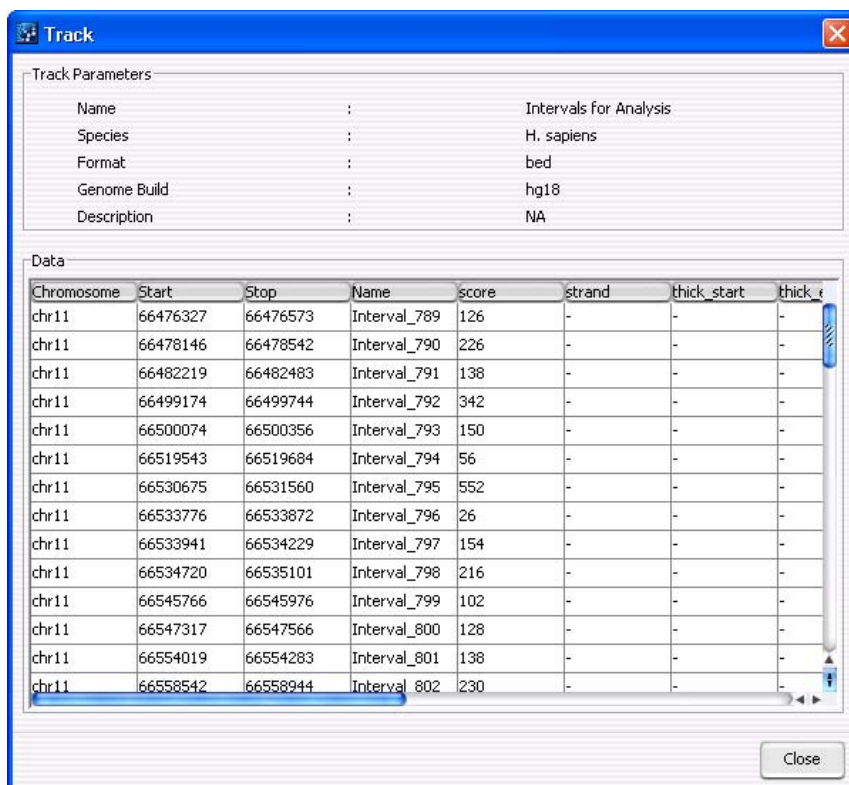


Figure 48 Track dialog box

Purpose: Lets you view the attributes of a specific track that is available in the program. You can view header information, and a list of all of the intervals in the track. All information is read-only. See [“To view the intervals in a target interval track”](#) on page 50.

To open: In the Experiment pane of the Navigator, right-click the name of the target intervals track within a specific experiment, then click **View Details**.

Alternatively, in the **Home** tab, click **User Preferences**. In the list of available tracks, click **Details** in the row of the desired track.

Track Parameters

This header information appears:

Item	Description
Name	The name of the track
Species	The name of the species to which the intervals in the track apply
Format	The format of the imported track file. The program supports BED format track files
Genome Build	The genome build of the applicable species upon which the track is defined
Description	A description of the track, if one is available.

Data

The Data pane of the dialog box shows a list of the intervals in the track, and all of the available BED file columns that are available for them. These columns can appear:

Column	Description
Chromosome	The chromosome on which the given interval is located.
Start	The location of the first base pair in the interval.
Stop	The location of the last base pair in the interval.
Name	The name of the interval that appears when the program displays the track.
Score	A number that represents how darkly the feature appears (higher numbers give darker results). The score can be a number from 0 to 1000.
Strand	The strand to which the annotation applies, which can be + (sense) or – (antisense)
thick_start	Each genomic feature can contain a thickly drawn region. This value is the starting position of this region.
thick_end	The ending position of the thickly drawn region of the genomic feature.

Column	Description
item_rgb	The display color of the data for the interval. The color is defined in the RGB color space. Example: An RGB value of 255, 0, 255 produces a bright fuchsia color.
block_count	The number of separate blocks for the given region. Typically, blocks refer to exons.
block_size	The size of each block. Multiple values are separated by commas.
block_starts	The starting position of each block relative to the location defined in Start . Multiple values are separated by commas.
exp_count	These columns use experiment scores and IDs to set the colors of a track. They are not relevant to SureSelect Quality Analyzer.
exp_ids	
exp_scores	

Close Closes the dialog box.

User Preferences – Miscellaneous tab

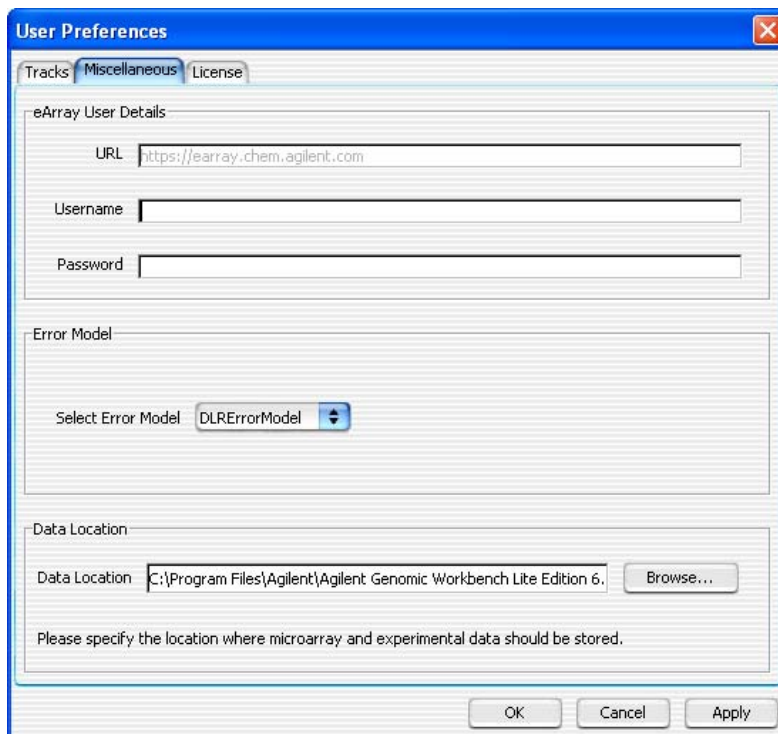


Figure 49 User Preferences dialog box – Miscellaneous tab

Purpose: For SureSelect Quality Analyzer, this dialog box lets you select a default location for analysis results and imported files. See [“To select where data and results are stored”](#) on page 19. This dialog box also lets you enter certain user preferences for other programs within Agilent Genomic Workbench.

To open: In the **Home** tab, click **User Preferences**, then click the **Miscellaneous** tab.

Alternatively, within the Genomic Viewer, right-click anywhere in Genome or Chromosome Views, or within the gene display area in Gene View, then click **User Preferences**. In the User Preferences dialog box, click the **Miscellaneous** tab.

- Data Location** Shows the default location for imported sequence read data and related files. The program also stores microarray sample data and analysis results for the CGH, ChIP-on-chip, and CH3 applications in this location.
- Browse** – Opens an Open dialog box, where you can select a default location.
- Other Options** All other options in this dialog box let you enter installation and configuration options for Agilent Genomic Workbench as a whole, or options that are specific for parts of the program other than the Target Enrichment Quality Analyzer.
- OK** Accepts any changes, and closes the dialog box.
- Cancel** Closes the dialog box, but does not save any changes.
- Apply** (Available after you make change(s) to any tab of the Preferences dialog box) Applies the changes that you made in any tab of the dialog box, but keeps the dialog box open.

User Preferences – Tracks tab

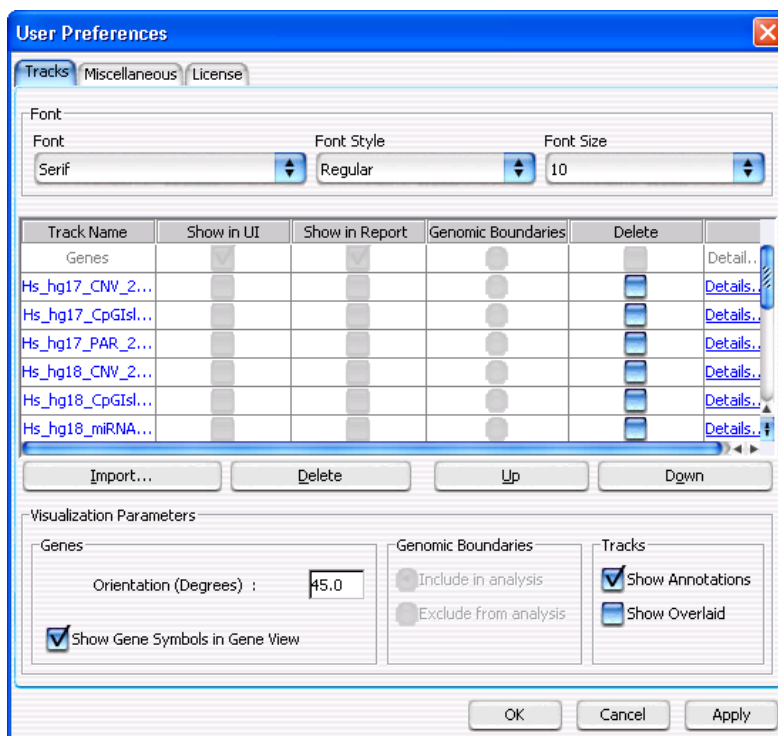


Figure 50 User Preferences dialog box – Tracks tab

Purpose: Lets you import and delete tracks from the program. Also lets you view the details of tracks and customize how genes and tracks appear in Gene View.

To open: In the **Home** tab, click **User Preferences**.

Alternatively, right-click anywhere within Genome or Chromosome Views, or within the gene display area in Gene View, then click **User Preferences**.

Font

Sets the font, style, and size of gene names and interval names in Gene View. These settings are available:

Setting	Description
Font	Select the desired font family.
Style	Select Regular , Bold , Italic , or Bold Italic .
Size	The size (in points) of gene and interval names. This is an absolute measure—the size of names does not change as you zoom in or out in Gene View.

List of Tracks – Columns

Shows a list of the tracks that are available in the program.

Track Name	The name of each track.
Show in UI	For the SureSelect Target Enrichment application type, the check boxes in this column are unavailable. The program automatically marks only Genes and the target interval track that has been defined for the selected experiment.
Show in Report	This column is not available for the SureSelect Target Enrichment application type.
Genomic Boundaries	The program automatically selects the target interval track that has been defined for the selected experiment.
Delete	Mark a check box to select a given track for deletion. To delete the selected track(s), click Delete .
Details	Opens the Track dialog box for the given track, where you can view a list of intervals in the track, as well as the header and other data that is associated with the track. See “Track” on page 139.

List of Tracks – Buttons

Import	Opens the Import Track dialog box, where you can select a BED format track file, and configure the track before you import it. See “Import Track” on page 131.
---------------	--

Delete Permanently removes selected tracks from the program. To select a track for deletion, mark its check box in the Delete column of the list of tracks.

Up, Down Changes the display order of tracks in Gene View. These commands are not relevant to the SureSelect Target Enrichment application type.

Visualization Parameters

Genes Orientation (Degrees) – The angle at which gene and interval names appear in Gene View. A value of 0° orients the names horizontally, and higher values rotate the name in a counterclockwise direction, up to a maximum of 360°.

Show Gene Symbols in Gene View – If you mark this option, the names of genes, as well as blue bars that show their genomic locations, appear in Gene View. If you clear this option, neither of these appears.

Genomic Boundaries These options are not available for the SureSelect Target Enrichment application type.

Tracks Mark any of these options:

- **Show Annotations** – In the target intervals track, displays the names of target intervals along with bars that represent their genomic locations.
- **Show Overlaid** – This option is only relevant to other programs in Agilent Genomic Workbench that can display more than one track at a time.

Commands that appear in all tabs

OK Accepts the change(s) that you made in any tab of the dialog box, and closes the dialog box.

Cancel Closes the dialog box, but does not make any changes.

Apply (Available if you have made changes to one or more preferences) Applies any change(s) that you have made in any tab of the dialog box, and leaves the dialog box open.

View Preferences

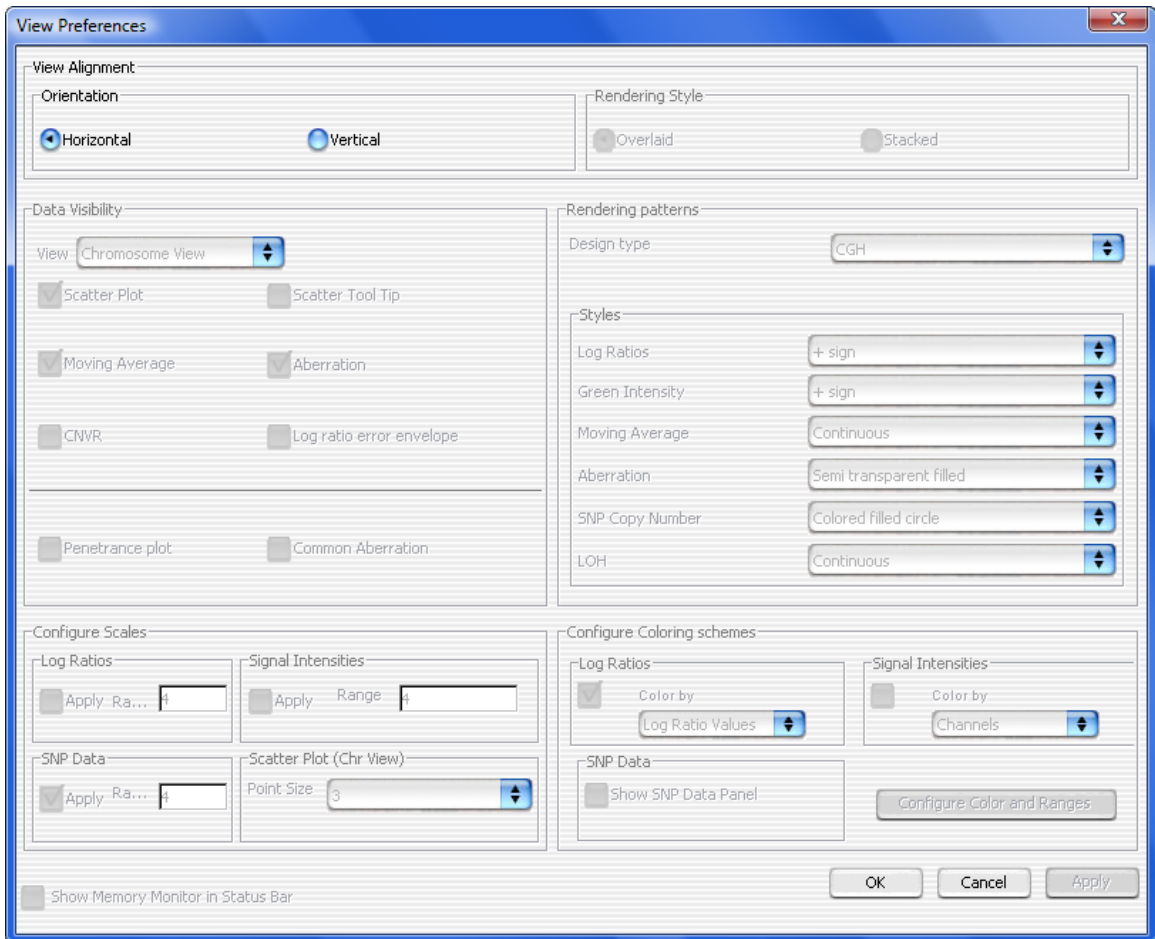


Figure 51 View Preferences dialog box

Purpose: For the SureSelect Target Enrichment application type, this dialog box lets you switch the Genomic Viewer between horizontal and vertical orientation.

To open: Click View > View Preferences.

View Alignment

Orientation The layout of Genome, Chromosome, and Gene Views. Select one of these options:

- **Horizontal** – Stacks the three views horizontally, from top to bottom. Because of the oblong shape of a typical monitor, Horizontal orientation can be somewhat more useful when you use the program to view sequence reads and the target interval track.
- **Vertical** – Stacks the three views vertically, from left to right.

In both cases, Tab view appears horizontally across the bottom of the Genomic Viewer.

Other commands

- OK** Accepts any changes you made, and closes the dialog box.
- Cancel** Closes the dialog box, but does not accept any changes.
- Apply** (Available only if you changed the orientation) Changes the orientation of the Genomic Viewer, and keeps the dialog box open.

No other options in this dialog box are applicable to the SureSelect Target Enrichment application type.

Target Enrichment Reports

This section describes the analysis results that are produced by SureSelect Quality Analyzer. These results/reports are available:

Result/Report	Description/Comments
Summary – PDF	<p>For this report, the program creates a PDF format report that contains tables of QC metrics. It also contains the results of read depth analysis shown both as tables and as a graph. See these topics:</p> <ul style="list-style-type: none"> • “To view and save reports” on page 56 • “QC metrics” on page 150 • “Read depth distribution” on page 152
Summary – Text	<p>For this report, the program creates a *.txt file that contains a table of QC metrics and the results of read depth analysis. You can open this file with a text editor or word processor. See these topics:</p> <ul style="list-style-type: none"> • “To view and save reports” on page 56 • “QC metrics” on page 150
Enrichment Analysis– UCSC	<p>For this report, the program creates a track file that contains enrichment analysis results for the chromosome that is currently selected in Genome View, and uploads it to the UCSC Genome Browser. See these topics:</p> <ul style="list-style-type: none"> • “To view and save reports” on page 56 • “Enrichment analysis results – UCSC” on page 153
Enrichment Analysis – XLS	<p>For this report, the program creates a *.xls file that contains enrichment analysis results for the chromosome that is currently selected in Genome View. You can view this file in Microsoft Excel. See these topics:</p> <ul style="list-style-type: none"> • “To view and save reports” on page 56 • “Enrichment analysis results – XLS report” on page 154

QC metrics

The table below describes the QC metrics that appear in both the PDF and *.txt format reports that are produced by Target Enrichment Quality Analyzer.

Metric	Description/comments
Total reads (good and bad)	Total number of reads contained in the read file in the selected experiment.
Total HQ uniquely mapped reads	The total number of high quality (HQ) input reads that map to a single genomic location. HQ reads are those that meet the quality criteria in SureSelect Quality Analyzer.
Read length	The number of bases in each of the input reads. SureSelect Quality Analyzer assumes all reads in the input file have the same length.
Total number of bases mapped	The total number of bases to which the reads in the selected experiment map. Typically, not all bases map to genomic regions.
Number of reads in target regions	The number of reads that overlap the regions that are defined in the target intervals track in the selected experiment. The program considers a read to be in a targeted region if the midpoint of the read falls within the targeted region.
Percent of reads in target regions	The fraction of the total reads that overlap one of the regions that are defined in the input target intervals track. The program considers a read to be in a targeted region if the midpoint of the read falls within the targeted region.
Percent of genome targeted	The fraction of the genome that is represented by the regions that are defined in the target intervals track in the selected experiment.
Enrichment in target regions	The overall fold-increase in representation of the target intervals over that expected for a completely random process.
Average read depth	The average number of reads that cover each targeted region. The program considers a read to be in a targeted region if the midpoint of the targeted region falls within the read.

Metric	Description/comments
Uniformity (3/4 mean with upper tail)	<p>A measure of how evenly reads are distributed. It is the percentage of all reads that fall in regions where the read depth is within a window defined by $\pm 75\%$ of the mean, around the mean, plus the entire upper tail of the distribution. Effectively, this is the percentage of all reads that fall in regions where the read depth is at least 25% of the mean read depth.</p> <p>Example: If the mean read depth is 38.42 reads, this metric gives the percentage of reads that fall in regions where the read depth is at least 9.61 reads.</p> <p>As the distribution of reads becomes broader, the Uniformity metric decreases. This metric can be useful in applications where you need a certain minimum level of coverage—for example, to make SNP calls. It lets you know the fraction of reads that fall in regions where the coverage is close to the mean, or above.</p>
Number of target regions	The number of intervals that were defined in the target intervals track in the selected experiment.
Number of target regions covered by at least n read(s)	<p>The number of regions in the target intervals track in the selected experiment that are covered by at least the indicated number of reads.</p> <p>Totals appear for n = 1, 5, 10, and 20.</p>
Percentage of target regions covered by at least n read(s)	<p>The percentage of regions in the target intervals track in the selected experiment that are covered by at least the indicated number of reads.</p> <p>Percentages appear for n = 1, 5, 10, and 20.</p>
Number of target bases covered by at least n read(s)	<p>The number of bases in targeted regions that are covered by at least the indicated number of reads.</p> <p>Totals appear for n = 1, 5, 10, 20, and 1000.</p>
Percentage of target bases covered by at least n read(s)	<p>The percentage of bases in targeted regions that are covered by at least the indicated number of reads.</p> <p>Totals appear for n = 1, 5, 10, 20, and 1000.</p>
Number of Bases in Target Regions	The total number of bases defined in the target intervals *.bed file for the selected experiment.
Percent target regions with zero coverage	The percentage of target regions that have no reads that overlap them by at least 1 bp.

Read depth distribution

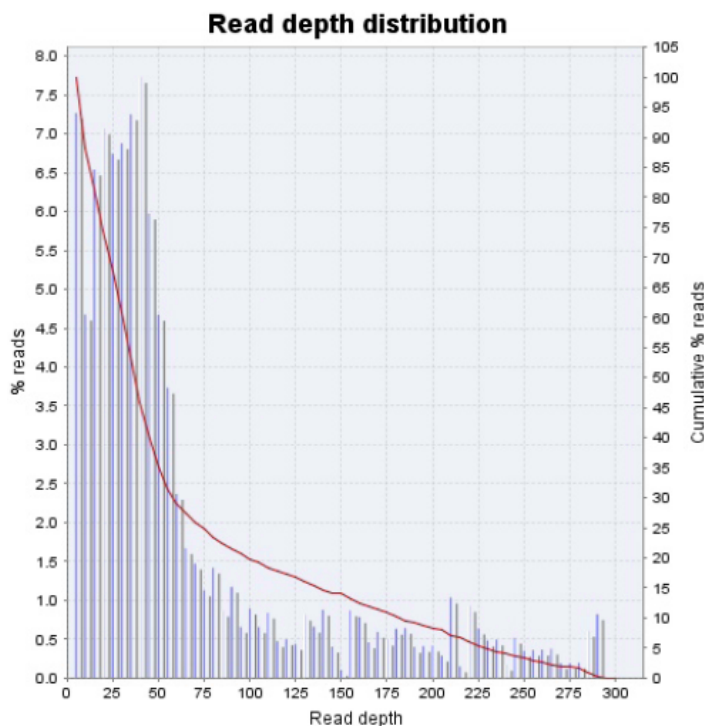


Figure 52 Distribution of read depths

The PDF format SureSelect QC report contains a graph that shows the distribution of read depths for the selected experiment. The program partitions the range of read depths into a fixed number of bins.

Bars (% Reads) – Shows the percentage of reads in each read depth bin. In general, a desirable distribution is gaussian, skewed toward higher values, with the peak at the desired level of coverage.

Red line (Cumulative % reads) – Shows the cumulative distribution function for read depths for the selected experiment. This function shows the probability that the read depth will be greater than or equal to a given depth.

Enrichment analysis results – UCSC

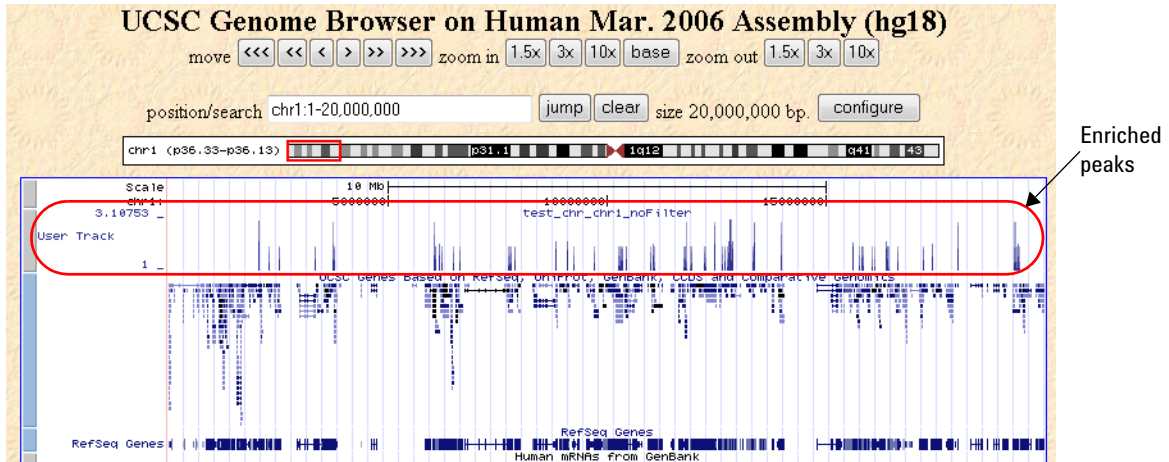


Figure 53 UCSC Genome Browser – Enriched peaks appear in the User Track area

The program displays enriched peaks for the selected chromosome as a histogram in the User Track area of the UCSC Genome Browser. It lets you see the fold-increase in the representation of each interval over that expected for a random process. See [“To view and save reports”](#) on page 56.

When you ask the program to display enrichment analysis results, it creates a WIG format file for the currently selected chromosome and transfers it to the UCSC Genome Browser. The UCSC Genome Browser opens in your Web browser.

You can also create an *.xls format file that contains a list of all of the enriched peaks for the selected chromosome. See [“Enrichment analysis results – XLS report”](#) on page 154.

Enrichment analysis results – XLS report

	A	B	C	D	E	F
1	Read File: s_1_1_export_top					
2	Target Regions: Example Targets					
3	Chromosome Definition: chrMapping.txt					
4	Chromosome: chr1					
5						
6	Interval Number	Chromosome	Start	Stop	Mid Point	Enrichment
7	1	chr1	879221	879520	879371	1
8	2	chr1	881200	881499	881350	1
9	3	chr1	947492	947791	947642	1
10	4	chr1	970915	971445	971411	2
11	5	chr1	1129627	1129926	1129777	1
12	6	chr1	1142911	1143210	1143061	1
13	7	chr1	1169203	1169502	1169353	1
14	8	chr1	1182149	1182448	1182299	1
15	9	chr1	1221278	1221577	1221428	1
16	10	chr1	1223817	1224116	1223967	1
17	11	chr1	1224568	1224867	1224718	1

Figure 54 Partial view of an enrichment analysis report (XLS format)

This report contains the results of enrichment analysis for the selected chromosome, organized by interval. You view this report in Microsoft Excel.

To create this report, see [“To view and save reports”](#) on page 56.

You can also view enriched intervals graphically in the UCSC Genome Browser. See [“Enrichment analysis results – UCSC”](#) on page 153.

Header information

Item	Description
Read File	The name of the sequence read file that was analyzed.
Target Regions	The name of the target intervals track that was used to analyze the read file. You supply this name when you import the track. See “To import target genomic regions as a track” on page 41.
Chromosome Definition	The name of the chromosome definitions file. This is a *.txt file that contains the names of the chromosomes that are referenced in the read file, correlated with the names of the chromosomes in the desired genome build of the species of interest. For details, see “To import sequence read files” on page 38.
Chromosome	The chromosome that was selected for enrichment analysis. Before you request an enrichment analysis report, you select the desired chromosome in Genome View of the Genomic Viewer.
Report generated by	The name and login ID of the person who requested the enrichment analysis report.

Data columns

Below the header information, these columns of data appear:

Column	Description
Interval Number	Intervals are numbered sequentially.
Chromosome	The name of the chromosome that contains the enriched interval
Start	The location of the first base of the enriched interval in the given chromosome
Stop	The location of the last base of the enriched interval in the given chromosome
Mid Point	The location of the midpoint of the enriched interval in the given chromosome
Enrichment	The fold-increase in representation of each interval over that expected for a purely random process.

www.agilent.com

In this book

This User Guide describes how to use the SureSelect Quality Analyzer application of Agilent Genomic Workbench to assess the quality of fragment pull-down for target enrichment experiments, both interactively and through workflows.

© Agilent Technologies, Inc. 2010

Revision A0, December 2010



G3800-90041



Agilent Technologies